


## 使用说明

### 2 系列充电触点 - 2 系列固定器


中文 (ZH)

## 1 安装

### 1.1 将充电触点或固定器固定在墙上

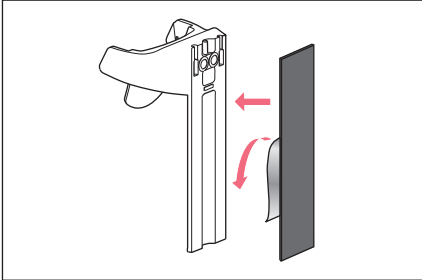

1. 清洁平滑的玻璃或不锈钢平面（用酒精等），并待其干燥。
2. 将胶带一面上的保护膜移除。
3. 将胶带牢固按压到充电触点或固定器背面。

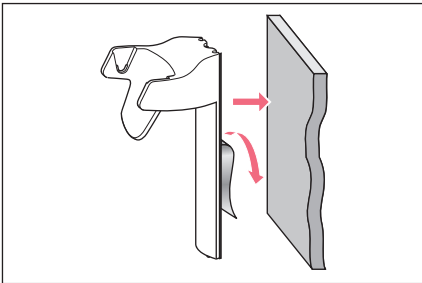

4. 将胶带第二面上的保护膜移除。
5. 将充电触点或固定器牢固按压到已清洁的表面上。  
须过 24 小时之后，才可对充电触点或固定器施加负载。

### 1.2 将充电触点或固定器从墙上取下

- ▶ 旋转充电触点或固定器并揭下胶带。

### 1.3 将充电触点或固定器装入旋转支架

充电触点或固定器无需胶带即可装入旋转支架。

## 2 操作

### 2.1 为电子定量器充电

在壁挂式安装的情况下无需通过固定器充电。定量器须以相应电源组件充电。

## 3 维护

### 3.1 清洁充电触点或固定器

- ▶ 用湿布和温和清洁剂拭去污物。

## 使用说明

2 系列充电触点 - 2 系列固定器

中文 (ZH)

### 3.2 材质

壁挂式支架	材料
开放平面	丙烯腈 - 苯乙烯 - 丙烯酸酯共聚物及聚碳酸酯 (ASA/PC)
充电适配器	镍镀层

## 4 订购信息

### 4.1 充电触点

订购信息 (International)	订购信息 (North America)	描述
3116 602.007	3116602007	<b>Charger Shell 2</b> for one Eppendorf Xplorer/Xplorer plus for Charger Carousel 2 (with charging functionality)
3116 603.003	3116603003	<b>Charger Shell 2</b> for one electronic Eppendorf Multipette (Repeater) for Charger Carousel 2 (with charging functionality)

### 4.2 固定器

订购信息 (International)	订购信息 (North America)	描述
3116 000.112	3116000112	<b>Holder 2</b> for one Eppendorf Research/plus, Reference/2 or Biomaster for Carousel 2, Charger Carousel 2 or wall mounting
3116 000.120	3116000120	<b>Holder 2</b> for one Eppendorf Xplorer/Xplorer plus for Carousel 2, Charger Carousel 2 or wall mounting (without charging functionality)
3116 000.139	3116000139	<b>Holder 2</b> for one electronic Eppendorf Multipette (Repeater) for Carousel 2, Charger Carousel 2 or wall mounting (without charging functionality)
3116 000.147	3116000147	<b>Holder 2</b> for one Eppendorf Multipette (Repeater) M4 for Carousel 2, Charger Carousel 2 or wall mounting

Your local distributor: [www.eppendorf.com/contact](http://www.eppendorf.com/contact)

Eppendorf AG · Barkhausenweg 1 · 22339 Hamburg · Germany

[eppendorf@eppendorf.com](mailto:eppendorf@eppendorf.com) · [www.eppendorf.com](http://www.eppendorf.com)