










Adjustable dispensing volumes

Applies to following dispensers:

- Multipette M4/Repeater M4
- Multipette plus/Repeater plus

Selection dial	Dispensing steps	 white	 light blue	 purple	 yellow	 green	 blue	 orange	 red	 light gray
•	100	1.0 µL	2.0 µL	5.0 µL	10 µL	25 µL	50 µL	0.1 mL	0.25 mL	0.5 mL
1	50	2.0 µL	4.0 µL	10 µL	20 µL	50 µL	100 µL	0.2 mL	0.50 mL	1.0 mL
•	33	3.0 µL	6.0 µL	15 µL	30 µL	75 µL	150 µL	0.3 mL	0.75 mL	1.5 mL
2	25	4.0 µL	8.0 µL	20 µL	40 µL	100 µL	200 µL	0.4 mL	1.00 mL	2.0 mL
•	20	5.0 µL	10 µL	25 µL	50 µL	125 µL	250 µL	0.5 mL	1.25 mL	2.5 mL
3	16	6.0 µL	12 µL	30 µL	60 µL	150 µL	300 µL	0.6 mL	1.50 mL	3.0 mL
•	14	7.0 µL	14 µL	35 µL	70 µL	175 µL	350 µL	0.7 mL	1.75 mL	3.5 mL
4	12	8.0 µL	16 µL	40 µL	80 µL	200 µL	400 µL	0.8 mL	2.00 mL	4.0 mL
•	11	9.0 µL	18 µL	45 µL	90 µL	225 µL	450 µL	0.9 mL	2.25 mL	4.5 mL
5	10	10 µL	20 µL	50 µL	100 µL	250 µL	500 µL	1.0 mL	2.50 mL	5.0 mL
•	9	11 µL	22 µL	55 µL	110 µL	275 µL	550 µL	1.1 mL	2.75 mL	5.5 mL
6	8	12 µL	24 µL	60 µL	120 µL	300 µL	600 µL	1.2 mL	3.00 mL	6.0 mL
•	7	13 µL	26 µL	65 µL	130 µL	325 µL	650 µL	1.3 mL	3.25 mL	6.5 mL
7	7	14 µL	28 µL	70 µL	140 µL	350 µL	700 µL	1.4 mL	3.50 mL	7.0 mL
•	6	15 µL	30 µL	75 µL	150 µL	375 µL	750 µL	1.5 mL	3.75 mL	7.5 mL
8	6	16 µL	32 µL	80 µL	160 µL	400 µL	800 µL	1.6 mL	4.00 mL	8.0 mL
•	5	17 µL	34 µL	85 µL	170 µL	425 µL	850 µL	1.7 mL	4.25 mL	8.5 mL
9	5	18 µL	36 µL	90 µL	180 µL	450 µL	900 µL	1.8 mL	4.50 mL	9.0 mL
•	5	19 µL	38 µL	95 µL	190 µL	475 µL	950 µL	1.9 mL	4.75 mL	9.5 mL
10	5	20 µL	40 µL	100 µL	200 µL	500 µL	1000 µL	2.0 mL	5.00 mL	10.0 mL