

Safety notice

Rotor S-4xUniversal (Centrifuge 5910 R/Centrifuge 5910 Ri)

English (EN)


This document is a supplement to the operating manual for the Centrifuge 5910 R and Centrifuge 5910 Ri and does not replace it. Therefore also read the operating manual of the centrifuge before using the rotors and accessories for the first time. The current version of the manual for all available languages can be found on our webpage www.eppendorf.com/manuals.



NOTICE! Material damage due to incorrect equipping of the swing-bucket rotor.

Incomplete equipping of the swing-bucket rotor or an uneven load will reduce the life span of the rotor and the corresponding buckets considerably.

- ▶ Always load all positions of a swing-bucket rotor with buckets.
- ▶ Load opposite buckets with the same weight (adapter, tubes, or plates and content).

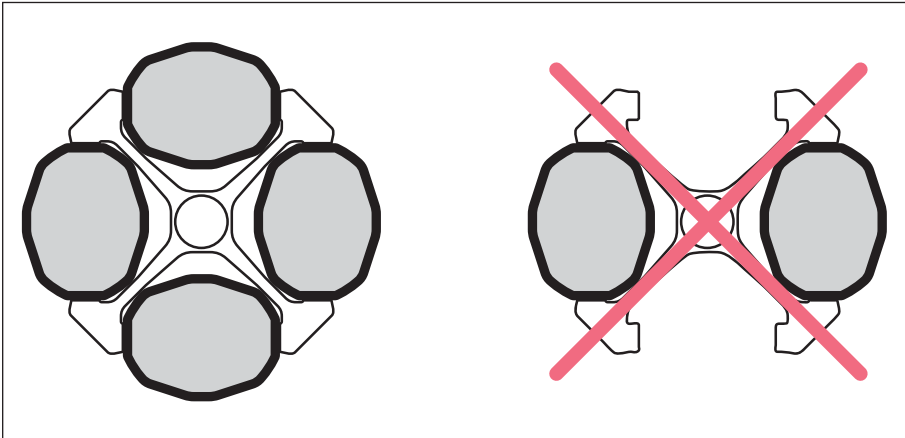

load difference
max. 1100 g



- ▶ Load adjoining buckets with a maximum of 1100 g difference in weight.

Your local distributor: www.eppendorf.com/contact

Eppendorf AG · Barkhausenweg 1 · 22339 Hamburg · Germany

eppendorf@eppendorf.com · www.eppendorf.com