## 1 Installation

### 1.1 System requirements

- Windows Vista SP2 (or higher)
- 32-bit processor (or higher), clock frequency 1 GHz (2 GHz recommended)
- USB 2.0 interface (or higher)
- 1 GB RAM
- 500 MB of free hard disk space

### 1.2 Installing the software

1. Insert the USB storage medium into a USB interface.
2. Open the directory and start the **setup.exe** program.  
The installation wizard starts.
3. Select the installation language and confirm with *OK*.  
The installation wizard is prepared.
4. Click on *Next* to continue.
5. Enter the user information and confirm with *Next*.
6. Select *Complete Installation* and confirm with *Next*.
7. Select *Install*.  
The software is installed.
8. Wait until the installation is complete and confirm with *Finish*.  
TrackIT has been installed and can be started.

### 1.3 Starting the software and setting up the reader

1. Start the software.
2. Select *Admin* as the login name.
3. Enter *admin* as the password.
4. Confirm with *Log in*.
5. Connect the reader to a USB interface.
6. In the *Configuration* menu, select the *Adjust reader* submenu.
7. In the *Configuration* window, select the serial port (ComPort) and confirm with *Close*.
8. In the *Configuration* menu, select the *Connect reader* submenu.  
The status display on the main screen shows: *Reader active*.
9. Select another ComPort if the reader remains inactive.

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