

1 Installation

1.1 System requirements

- Windows 7 SP1 (or higher)
- Windows 8
- Windows Vista SP2 (or higher)
- Windows XP SP3 (not recommended)
- Windows 10
- 32 bit processor (or higher), clock frequency 1 GHz (2 GHz recommended)
- USB 2.0 interface (or higher)
- 1 GB RAM
- 500 GB of free hard disk space

1.2 Installing the software or a software update

- ▶ Insert the USB storage medium in a USB interface.
- ▶ Open the directory and start the **setup.exe** program.
The installation wizard is launched.
- ▶ Select the installation language and confirm with *OK*.
The installation wizard is prepared.
- ▶ Click on *Next* to continue.
- ▶ Enter the user information and confirm with *Next*.
- ▶ Select *Complete installation* and confirm with *Next*.
- ▶ Select *Install*.
The software is installed.

- ▶ Wait for the installation to complete and confirm with *Finish*.

Eppendorf TrackIT is now installed and can be started.

1.3 Starting the software and setting up the reader

- ▶ Start the software.
- ▶ Select *Admin* as the login name.
- ▶ Enter *admin* as the password.
- ▶ Confirm with *Log in*.
- ▶ Connect the reader to a USB interface.
- ▶ In the *Configuration* menu, select the *Adjust reader* submenu.
- ▶ In the *Configuration* window, select the serial port (ComPort) and confirm with *Close*.
- ▶ In the *Configuration* menu, select the *Connect reader* submenu.
The status display on the main screen shows: *Reader active*.
- ▶ Select another ComPort if the reader remains inactive.

Your local distributor: www.eppendorf.com/contact

Eppendorf AG · Barkhausenweg 1 · 22339 Hamburg · Germany
eppendorf@eppendorf.com · www.eppendorf.com