

Concentrator plus/Vacufuge® plus – Instructions for Setting Up the Device

These Instructions supplement the Concentrator plus/Vacufuge® plus operating manual. They do not replace it. Be sure to read the operating manual before you begin to use the device.

Transport

The device weighs 31 kg (complete system) or 17 kg (basic device).


- ▶ Transport the device only when it is in its original packaging.
- ▶ Two persons are always required to transport it.
- ▶ Use a dolly or other mechanical help for longer distances.

Setting Up

- ▶ Follow these steps in order; two persons are required.

1. Open the box.

- a) Cut the straps.
- b) Cut open the tape.
- c) Open all 4 box flaps.
- d) Remove the paperboard covering.
- e) Take out the accessories.

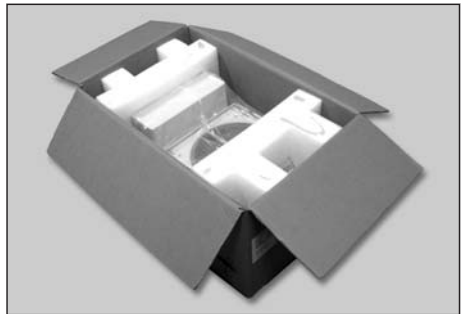

2. Take the device out of the box.

- a) With its transport protection pads still around it, lift the device out of the box.
- b) Carefully lay the device on its side, with transport protection pads.
- c) Gently slide back the rear transport protection pad until the second transport protection screw is visible.



1 2 2 3 4


- 1 Rear transport protection pad
- 2 Transport protection screws
- 3 Bottom transport protection pad
- 4 Front transport protection pad

eppendorf

Concentrator plus/Vacufuge® plus – Instructions for Setting Up the Device

3. Remove the transport protection pads.

- a) Pull out the bottom piece of transport protection pad.
- b) Unscrew and remove both transport protection screws.
- c) Set the device upright.
- d) Remove the front and rear transport protection pads.
- e) Remove the plastic wrapping from the device.
- f) Remove the plastic wrapping from the cover.
- g) Set the device on a stable, level, and resonance-free laboratory workbench. Let it rest 3 hours to acclimatize.

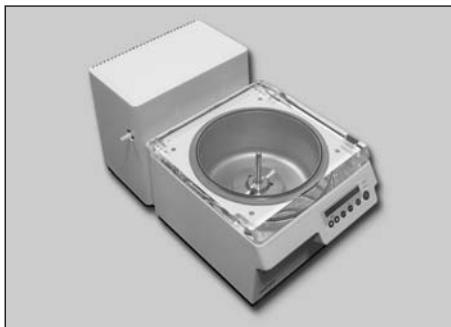

- ▶ Retain the box as well as the transport protection pads and screws to transport or store the device later.



Caution: It is dangerous to ignore the operating manual.

The device is not yet ready to operate. You must attach further components.

- ▶ Be sure to read the operating manual included here before you begin to use the device.

eppendorf
In touch with life

Your local distributor: www.eppendorf.com/worldwide

Eppendorf AG · 22331 Hamburg · Germany · Tel: +49 40 538 01-0 · Fax: +49 40 538 01-556
E-Mail: eppendorf@eppendorf.com

Eppendorf North America, Inc. · One Cantiague Road · P.O. Box 1019 · Westbury, N.Y. 11590-0207 · USA
Tel: +1 516 334 7500 · Toll free phone: +1 800 645 3050 · Fax: +1 516 334 7506 · E-Mail: info@eppendorf.com

Application Support

Europe, International: Tel: +49 1803 666 789 · E-Mail: support@eppendorf.com
North America: Tel: +1 800 645 3050 ext. 2258 · E-Mail: support_na@eppendorf.com
Asia, Pacific: Tel: +60 3 8023 6869 · E-Mail: support_asiapacific@eppendorf.com