



Take Full Advantage of Versatility

Missing a second incubation temperature for your Eppendorf ThermoMixer®?

The Eppendorf SmartExtender offers a comfortable incubation tool which can easily be used as an add-on to your existing SmartBlocks and the Eppendorf ThermoMixer C/Fx, as well as Eppendorf ThermoStat™ C.

- > Up to 12 vessels (1.5 mL) for higher throughput and higher flexibility
- > Independent heating function for second temperature
- > Parallel usage of Eppendorf SmartBlock™ and SmartExtender



Benefits

- > Temperature control is independent from Eppendorf SmartBlock (second temperature) for second application
- > SmartExtender can actively heat from 3 °C above room temperature up to 110 °C to cover all common applications
- > All SmartBlock sizes are usable in parallel to the SmartExtender by their full functionality (heating, cooling, mixing) for high flexibility

Conditions

- > Fits on Eppendorf ThermoMixer C, ThermoStat C, and ThermoMixer Fx for flexible usage
- > Instrument firmware update for installed unit needed, available on www.eppendorf.com/thermomixer
- > Take advantage of existing Eppendorf ThermoTop® power support connectors (SmartExtender or ThermoTop)



Ordering information

Description	Order no.
Eppendorf ThermoMixer® C, basic device without Eppendorf SmartBlock™, 220 V–240 V	5382 000.015
Eppendorf ThermoMixer® F1.5, with Eppendorf SmartBlock™ for 24 reaction vessels 1.5 mL, incl. Transfer Rack, 220 V–240 V	5384 000.012
Eppendorf ThermoMixer® FP, with Eppendorf SmartBlock™ for microplates and deepwell plates, including lid, 220 V–240 V	5385 000.016
Eppendorf ThermoMixer® F0.5, with Eppendorf SmartBlock™ for 24 reaction vessels 0.5 mL, incl. Transfer Rack, 220 V–240 V	5386 000.010
Eppendorf ThermoMixer® F2.0, with Eppendorf SmartBlock™ for 24 reaction vessels 2.0 mL, incl. Transfer Rack, 220 V–240 V	5387 000.013
Eppendorf ThermoStat™ C, basic device without Eppendorf SmartBlock™, 220 V–240 V	5383 000.019
Eppendorf ThermoTop®, with condens.protect® technology	5308 000.003
Eppendorf SmartBlock™ 0.5 mL, thermoblock for 24 × 0.5 mL tubes, incl. Transfer Rack	5361 000.031
Eppendorf SmartBlock™ 1.5 mL, thermoblock for 24 × 1.5 mL tubes, incl. Transfer Rack	5360 000.038
Eppendorf SmartBlock™ 2.0 mL, thermoblock for 24 × 2.0 mL tubes, incl. Transfer Rack	5362 000.035
Eppendorf SmartBlock™ 5.0 mL, thermoblock for 8 × Eppendorf Tubes® 5.0mL	5309 000.007
Eppendorf SmartBlock™ 15 mL, thermoblock for 8 × 15 mL conical tubes	5366 000.021
Eppendorf SmartBlock™ 50 mL, thermoblock for 4 × 50 mL conical tubes	5365 000.028
Eppendorf SmartBlock™ 12 mm, thermoblock for 24 reaction tubes, diameter up to 12 mm	5364 000.024
Eppendorf SmartBlock™ cryo, thermoblock for 24 cryo tubes, 1.5–2.0 mL, all base shapes	5367 000.025
Eppendorf SmartBlock™ plates, thermoblock for microplates and deepwell plates, incl. lid	5363 000.039
Eppendorf SmartBlock™ PCR 96, thermoblock for PCR plates 96, incl. lid	5306 000.006
Eppendorf SmartBlock™ PCR 384, thermoblock for PCR plates 384, incl. lid	5307 000.000
Eppendorf SmartBlock™ DWP 500, thermoblock for Eppendorf Deepwell Plates 96/500 µL, incl. lid	5316 000.004
Eppendorf SmartBlock™ DWP 1,000, thermoblock for Eppendorf Deepwell Plates 96/1,000 µL, incl. lid	5310 000.002
Eppendorf SmartExtender™, heating rack for 12x 1.5/ 2.0 mL	5322 000.008

Your local distributor: www.eppendorf.com/contact

Eppendorf AG · Barkhausenweg 1 · 22339 Hamburg · Germany
eppendorf@eppendorf.com · www.eppendorf.com

www.eppendorf.com/thermomixer