Transporting, Installing, Transferring and Disposing of the Centrifuge 5424 R

These instructions for use complement the operating manual for the Centrifuge 5424 R and do not replace it. You should therefore also read the operating manual before using the centrifuge for the first time. The current version of the operating manual can be found in the Internet at www.eppendorf.com.

1 Transport

- ▶ Only transport the device in the original packaging.
- ▶ The centrifuge must always be transported by two people.
- ▶ Use a transport aid for transporting over longer distances.

2 Installation

The weight of the centrifuge is 21 kg. The centrifuge must always be transported by two people.

2.1 Open the carton and lift the device

1. Cut the adhesive tape.
2. Open all 4 flaps of the carton.
3. Remove the accessories.
4. Reach with your hands under the device and lift the centrifuge from the carton with another person.

2.2 Place the device on a stable, horizontal and non-resonant lab bench.

1. Remove the front and back transport protection pads.
2. Remove the plastic sleeve
3. Observe a minimum distance of 30 cm to adjoining devices at the sides and from the rear side to the wall.
4. Install the device in a well-ventilated location which is protected from direct sunlight to prevent it from overheating.

2.3 Connect the device

1. After installation, wait for four hours before switching the centrifuge on in order to avoid damage to the compressor.
2. Check that the mains voltage and frequency match the requirements on the device name plate (see rear side of the device) and then connect the device to the power supply.

2.4 Remove the transport protection pad from the rotor chamber

1. Switch on the device at the mains power switch.
2. Open the centrifuge lid using the open key.
3. Only for the device version with rotor (left Fig.):
 - ▶ Turn the rotor nut counterclockwise using the rotor key supplied.
 - ▶ Remove the rotor by lifting vertically.
 - ▶ Remove the transport protection pad.
4. Only for the device version without rotor (right Fig.):
 - ▶ Remove the transport protection pad.
5. Place the rotor vertically onto the motor shaft.
6. Turn the rotor nut using the rotor key clockwise until the rotor nut is tightened.

2.5 The device is now ready to use.

- ▶ Retain the packaging and all transport protection pads for shipping the device at a later date.

3 Packing

- ▶ Pack the centrifuge in reverse order.

4 Passing On the Device

- ▶ When passing the device on to third parties, be sure to include the operating manual and these instructions for use.

5 Disposal

Observe the applicable legal regulations when disposing of the device. The disposal is regulated within the European Community by national regulations based on EU Directive 2002/96/EC on waste electrical and electronic equipment (WEEE). According to these regulations, any devices supplied after August 13, 2005 on a business-to-business basis, to which this product is assigned, may no longer be disposed of in municipal or household waste. For more information, please contact your supplier.

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