



ThermoStat plus – Short Instructions

These short instructions may only be used by trained personnel who are familiar with the operating manual and the device. Observe the safety precautions.


Control element overview

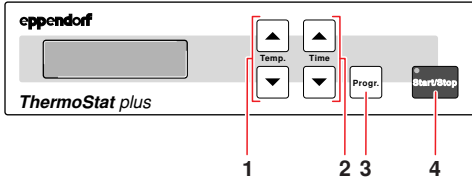

- 1 Adjust temperature
- 2 Adjust time
- 3 Open program function
- 4 Start or stop run

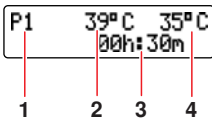



- 1 Phase number
- 2 Target temperature value
- 3 Tempering duration
- 4 Actual temperature value

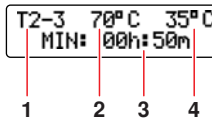


- 1 Temperature change phase number
- 2 Target temperature value
- 3 Temperature change phase duration
- 4 Current actual temperature value

Simple tempering


After the device has been switched on, it is in simple tempering mode.

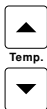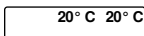



1. Adjust the temperature using the "Temp." arrow keys. The temperature between -5°C and 99°C is readjusted immediately. In addition you can switch on the timer.

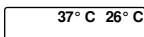

2. To start the timer, press the "Start/Stop" key. The control lamp illuminates in green. The tempering duration ("--h : --m") is displayed.

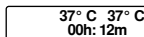


3. To stop the timer, press the "Start/Stop" key. The timer is no longer shown in the display and the control lamp goes out. The tempering continues with the target value set. The values remain stored even after the device has been switched off.

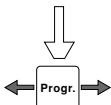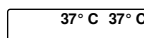

ThermoStat plus – Short Instructions

Programmable tempering

In addition to simple tempering you can also create a program with up to 4 different automatically sequenced phases. Here you can adjust the desired temperature and the desired tempering duration.

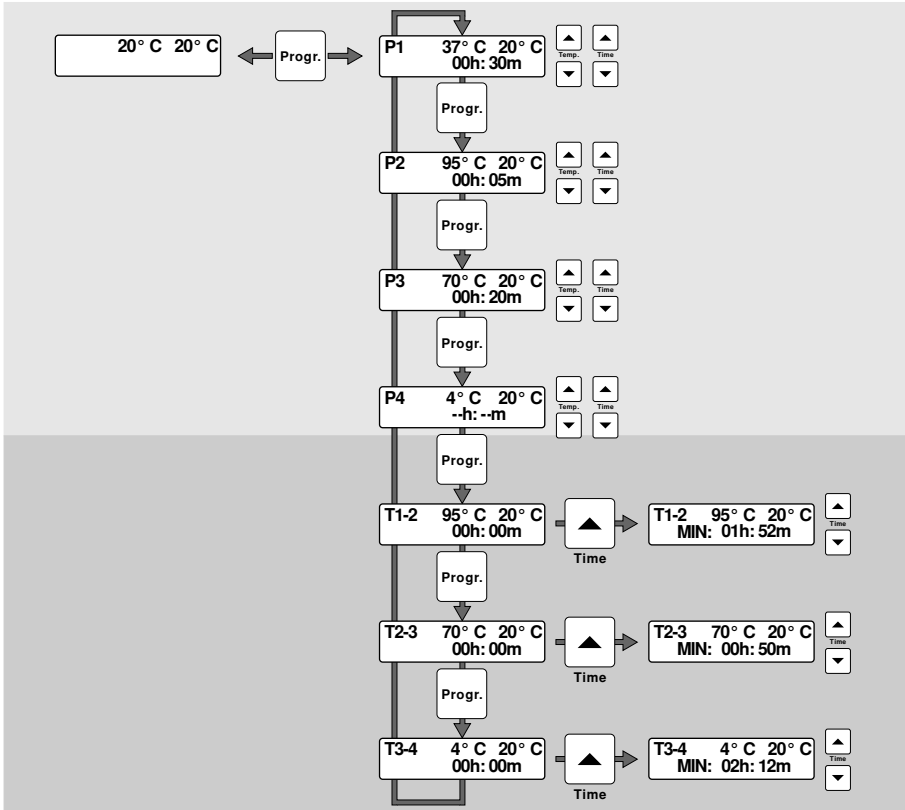

eppendorf
In touch with life

Your local distributor: www.eppendorf.com/worldwide

Eppendorf AG · Germany · Tel. +49 40 538 01-0 · E-Mail: eppendorf@eppendorf.com
Eppendorf North America, Inc. · USA · Toll free phone 800 645 3050 · E-Mail: info@eppendorf.com

Application Support

Europe, International: support@eppendorf.com · North America: support_NA@eppendorf.com
Asia, Pacific: support_AsiaPacific@eppendorf.com