

RS-232 OPC Server Update – Version B

Description of changes:

Update resolves communication errors with PC COM ports.

Instructions:

Before running the update:

- > End all runs on BioCommand.
- > Close BioCommand and any other software that is related to OPC server.

Perform the upgrade as listed below. You may want to print this page for easy access to the instructions while you work.

1. Go to the Saved file on your hard drive and double click the file
2. Pop up menu comes up – Select Run
3. From the pop up menu - under "Unzip to folder" – click "Browse" and select desired location on your hard-drive (suggest "desktop")
4. Select Unzip button and Okay
5. After download is complete select "Close" to close the dialog box
6. Open the folder 'M1295 RS232 Device Interface Kit 1252 Rev B.00 – Update' created on your hard drive.
7. Select setup.exe and follow instructions.
8. After the installation is complete, close all.

Congratulations! You have installed the latest performance enhancements. Check the Eppendorf Website regularly for new revisions.