

Analog Input/Output Module Update – Version Analog B.00

Description of changes:

Analog I/O software revision B.00 includes the ability to recognize and control up to 8 Analog I/O modules from a single installed version, and addresses previous issues with communication loss.

Instructions:

Determine which software version is currently installed in your controller:

- > Close BioCommand and any other client software that is related to Analog Input Output Module.
- > Select Help -> About on Analog Input Output Module.
- > View version information
- > Close the Analog I/O Module OPC Server Software.

If the installed software is lower than the latest version listed above, perform the upgrade as listed below. You may want to print this page for easy access to the instructions while you work.

1. Go to the Saved file on your hard drive and double click the file
2. Pop up menu comes up – Select Run
3. From the pop up menu - under "Unzip to folder" – click "Browse" and select desired location on your hard-drive (suggest "desktop")
4. Select Unzip button and Okay
5. After download is complete select "Close" to close the dialog box
6. Open the folder 'Analog IO B.00' created on your hard drive.
7. Select setup.exe and follow instructions.
8. After the installation is complete, start Analog Input Output Module.
9. Verify that the new version is installed by viewing Version information in Help -> About.
10. Analog Input Output Module will now be running with the new version that was installed on your PC.

Congratulations! You have installed the latest performance enhancements. Check the Eppendorf Website regularly for new revisions.