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eppendorf is a registered trademark.  
4860 900.179-01/0208 · Printed in Germany · Printed on chlorine-free bleached paper

**eppendorf**

## Charging stand / Support-chargeur

### Charging stand

The charging stand is recommended for storing and recharging the Research pro. Two different versions of the charging stand are available, for one or four Research pro instruments. The delivery package of the charging stand contains a power unit. Before using the power unit for the first time, please check that the voltage requirements in your lab correspond to the specifications on the I.D. plate. Do not use any other power unit for the charging stand. The hollow plug of the power unit is inserted into the charging stand. The power unit is then plugged into a mains socket. The charging stand is now ready to use.

 Make sure that the battery is inserted before recharging the Research pro.

For charging purposes, the Research pro is placed in the charging stand. Before recharging, please compare your voltage requirements with the voltage specifications on the plug-in unit. For charging purposes, the charging adapter is placed upon the Research pro. Alternatively, the Research pro can be placed in the charging stand.

During the start and termination of the charging process, "CON" appears briefly in the display. The display is switched on during the charging process. The battery symbol appears in a rolling form. During charging, the Dispensing function is frozen. When the charging process has finished, the battery symbol appears in the display without moving. When the pipette has been charged, it is stored in the charging stand.

#### Technical data for power unit

Input voltage (country-specific): 240 V  $\pm$ 10 %, 230 V  $\pm$ 10 %, 120 V  $\pm$ 10 %, 100 V  $\pm$ 10 %; 50/60 Hz

Output voltage: 7.5 V DC; 800 mA/6 VA

Technical specifications subject to change!

#### Ordering information

	International	North America
Charging stand for one pipette	4860 000.267	022461541
Charging stand for four pipettes	4860 000.860	022461559

If the delivery address is not in the country in which the device shall be used:

When placing your order, please specify the country in which the device is to be used and the voltage used in that country. Only then is it possible for us to deliver the power unit of the charging unit/stand with the necessary input voltage.

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