Keep this rotor instruction manual and the centrifuge manual in the file.

Angle Rotors for Ultracentrifuge

P50AT2

Before using these rotors, please carefully read this instruction manual, "ROTORS, TUBES, BOTTLES AND CAPS", and the centrifuge instruction manual for its efficient operation and for your safety.

Keep this instruction manual for your reference and refer to it as required.

Contents

| | | |
|-----------------------------|-------|---|
| 1. Specifications | _____ | 1 |
| 2. Applicable Tubes/Bottles | _____ | 2 |
| 3. How to Use | _____ | 3 |
| 4. Maintenance | _____ | 5 |
| 5. Using Cesium Chloride | _____ | 7 |
| 6. Packing List | _____ | 8 |
| 7. Decontamination | _____ | 9 |
| 8. Rotor Retirement | _____ | 9 |

Order no. 5720211107 (P50AT2 with RLM adapter)
Order no. 5720211117 (P50AT2 with Optical adapter)
S999181-18/2025-05

Original instruction manual

Safety Reminder

- Centrifuge rotors rotating at high speed have considerable potential for damage to personal properties if used improperly. For safe and proper use of this rotor, carefully read the centrifuge instruction manual and this rotor instruction manual before use and observe the instructions.

 **WARNING :** and  **CAUTION :** notes are used to call your attention in this manual to prevent personal injury or damage to the rotor and the centrifuge. These notes are defined as follows.

 **WARNING:** indicates a potentially hazardous situation which, if not avoided, could result in personal severe injury or possible death.

 **CAUTION:** indicates a hazardous situation which, if not avoided, could result in personal injury or severe damage to the instrument.

WARNING

- Never use any material capable of producing flammable or explosive vapors. Your centrifuge and rotor have no explosion-proof construction.
- Your centrifuge and rotor are not designed to confine any sample particles dispersed due to leakage. Therefore, when using toxic or radioactive samples or pathogenic or infectious blood samples, make sure to prepare necessary safety measures at your own responsibility.
- Never exceed the maximum speed of the rotor (mentioned on the rotor). Always reduce rotor speed as instructed in this manual when rotor speed is limited due to sample density or kinds of tubes.
- Check the chemical resistance chart attached the rotor, and do not use any sample inapplicable to the rotor. Using such a sample could corrode the rotor.
- Do not remove the RLM adapter or optical adapter from rotor, or replace it with the adapter for another rotor. The adapter is critical component that detects the over-speed of rotor. If an adapter that is not compatible with the rotor is attached, the rotor could break, resulting in damage to the centrifuge.
- Do not allow the rotor temperature to rise 100°C.: This would cause the material brittle.
- Never sterilize the rotor by autoclaving or boiling. Otherwise the strength of rotor may be significantly decreased.
- If the centrifuge, rotor, or an accessory is contaminated by samples that toxic or radioactive, or blood samples that pathogenic or infectious, be sure to decontaminate the item according to good laboratory procedures and methods.
- If there is a possibility that the centrifuge, rotor, or an accessory is contaminated by samples that might impair human health (for example, samples that are toxic or radioactive, or blood samples that are pathogenic or infectious), it is your responsibility to sterilize or decontaminate the centrifuge, rotor, or the accessory properly before requesting repairs from an authorized sales or service representative.
- It is your responsibility to sterilize and/or decontaminate the centrifuge, rotor, or parts properly before returning them to an authorized sales or service representative.

CAUTION

- Check the chemical resistance chart attached to the rotor, and do not use any sample inapplicable to the tubes/bottles, the tube adapter, or the cap for the tubes/bottles. Using such a sample could deteriorate them.
- The allowable speed may be lower depending on the density of sample, use of salt-such as cesium chloride or the combination of tubes or bottles and tube adapters and caps. Refer to this instruction manual.
- Always balance the tubes/bottles and samples within the allowable imbalance of rotor, including the caps and the tube adapters. Do not exceed the allowable imbalance. Otherwise the rotor and the centrifuge may be damaged.
- Do not load only one tube or load tubes asymmetrically : Asymmetrically loading may cause imbalance operation and damage the centrifuge and rotor.
- Clean the inside of the drive hole (crown hole) of the rotor and the surface of the drive shaft (crown) of the centrifuge once a month. If the drive hole or the drive shaft is stained or any foreign matter is adhered, the rotor may be improperly installed and come off during operation.
- Use the rotor tubes and bottles within their actual capacities.
- Do not use tubes/bottles that have exceeded their life expectancy. Failure to do so could result in damage of tubes/bottles and the rotor and the centrifuge. The life expectancy of tubes/bottles depends on factors such as the characteristics of samples, speed of the rotor used, and temperature. Always check for deterioration and damage (cracks, deformation, and so on) on tubes/bottles before using them. Do not use the tubes/bottles if you find such a problem.
- Inspect and maintain the rotor after use. If abnormality is observed, do not use it. Contact an authorized sales or service representative.

1. Specifications

1.1 Construction

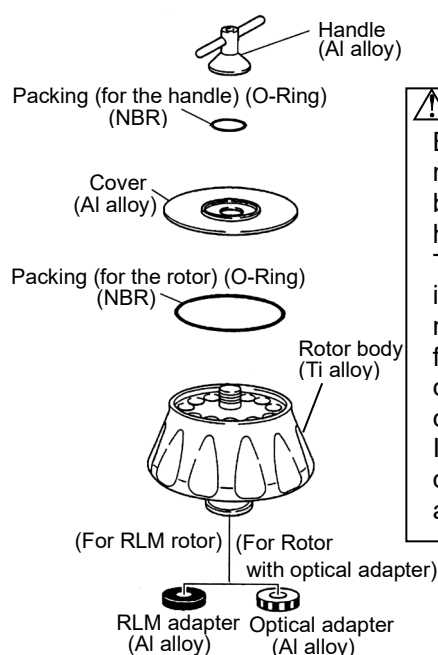
Fig.1-1. Construction of the P50AT2 Rotor

1.2 Specifications

- Max. speed: 50,000 rpm
- Max. RCF: 303,000 xg
- Allowable density at maximum speed of the rotor: 1.2 g/mL
- K Factor at maximum speed of the rotor: 70
- Nominal rotor capacity: 40 mL × 12 tubes = 480 mL
- Material of rotor: Ti alloy
- Material of cover: Al alloy
- Tube size: $\phi 2.6 \times 9.0$ cm
- Weight: 13.0 kg
- Max. diameter: 22.9 cm
- Max. height: 22.2 cm

⚠ CAUTION:

- Do not use a RLM rotor with the centrifuge other than a series of a CP-NX/WX/MX/ α ultracentrifuge.
- Do not use rotor with optical adapter with model 55P and 40P (old type centrifuge).

1.3 Cross-Sectional View of the Rotor

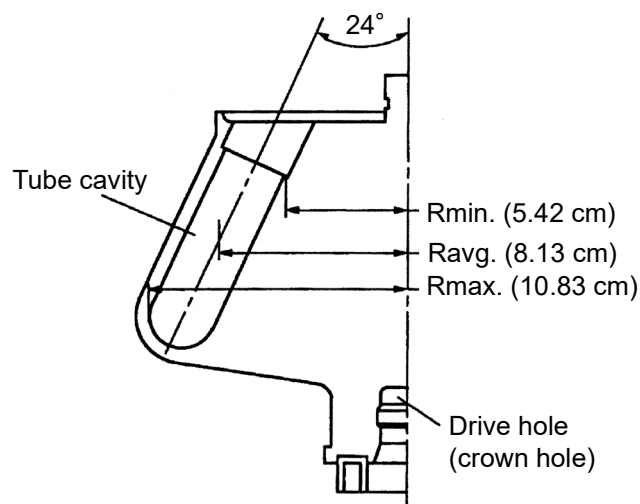
Fig.1-2. Cross-Sectional View of the P50AT2 Rotor

⚠ CAUTION:

Clean the inside of the drive hole (crown hole) of the rotor and the surface of the drive shaft (crown) of the centrifuge once a month. If the drive hole or the drive shaft is stained or any foreign matter is adhered, the rotor may be improperly installed and come off during operation.

1.4 Characteristic of the Rotor

Table 1-1. Characteristic of the P50AT2

| Rotor Speed (rpm) | RCF (xg) | | | K factor |
|-------------------|----------|---------|---------|----------|
| | Rmax. | Ravg. | Rmin. | |
| 10,000 | 12,100 | 9,090 | 6,060 | 1,753 |
| 20,000 | 48,400 | 36,400 | 24,200 | 438 |
| 30,000 | 109,000 | 81,800 | 54,500 | 195 |
| 40,000 | 194,000 | 145,000 | 97,000 | 110 |
| 50,000 | 303,000 | 227,000 | 151,000 | 70 |

※ For detail about K factor, refer to "ROTORS, TUBES, BOTTLES AND CAPS (Part No.S999204)" attached to the centrifuge.

Calculation formula

$$RCF = 1.118 \times 10^{-5} \times R \times N^2 \text{ (xg)}$$

R: Rotating radius (cm)

N: Speed (rpm)

2. Applicable Tubes/Bottles

2.1 Applicable Tubes/Bottles

Table 2-1. Applicable Tubes/Bottles

| Tubes/Bottles | | | | | Tube adapter Float spacer | | Cap | | Max. speed (rpm) | Max. RCF (xg) |
|-----------------------------|------------|-------------------------------------|--------------------------|---------------------------------|------------------------------|---------------------|------------|--------------------------------|------------------------|---------------------|
| Nominal capacity (mL) | Order no. | Name | Size (ϕ X L cm) | Actual capacity (mL/tube) | Order no. | Name | Order no. | Name | | |
| 40 | 5720411139 | 40PP tube | 2.6 X 9.0 | 35.7 | --- | --- | 5720411212 | E3-AL cap ass'y | 50,000 | 303,000 |
| | | | | 32.5 | --- | --- | 5720410102 | S-40AL cap ass'y ^{*1} | 50,000 | 303,000 |
| | 5720411141 | 40PP seal tube | 2.6 X 9.4 | 39.0 | --- | --- | 5720411256 | E space cap | 50,000 | 303,000 |
| | | | | 35.7 | --- | --- | 5720411212 | E3-AL cap ass'y | 40,000 | 194,000 |
| | 5720411146 | 40PE tube | 2.6 X 8.9 | 35.7 | --- | --- | 5720411223 | E4-AL cap ass'y | 40,000 | 194,000 |
| | | | | 31.8 | --- | --- | 5720410102 | S-40AL cap ass'y ^{*1} | 40,000 | 194,000 |
| | | | | 37.0 | --- | --- | 5720411212 | E3-AL cap ass'y | 40,000 | 194,000 |
| | 5720411148 | 40PET tube | 2.6 X 9.0 | 35.7 | --- | --- | 5720411212 | E3-AL cap ass'y | 45,000 | 245,000 |
| | | | | 24.4 | --- | --- | 5720411219 | E-SST Lid (B) | 38,000 | 175,000 |
| | 5720411149 | 40TI tube | 2.6 X 9.0 | 24.4 | --- | --- | --- | --- | 38,000 | 175,000 |
| | | | | 24.4 | --- | --- | 5720411219 | E-SST Lid (B) | 32,000 | 124,000 |
| | | | | 24.4 | --- | --- | --- | --- | 32,000 | 124,000 |
| | 5720411147 | 40SST tube (B) | 2.6 X 9.0 | 24.4 | --- | --- | --- | --- | 32,000 | 124,000 |
| | | | | 24.4 | --- | --- | --- | --- | 32,000 | 124,000 |
| 33 | 5720411137 | 33PC thick-walled tube | 2.6 X 8.9 | 20.4 | --- | --- | --- | --- | 40,000 | 194,000 |
| | 5720411138 | 33PP thick-walled tube | 2.6 X 8.8 | 20.4 | --- | --- | --- | --- | 25,000 | 75,700 |
| 30 | 5720411126 | 30PC bottle (C3) ass'y | 2.6 X 9.7 | 26.0 | --- | --- | --- | --- | 50,000 | 303,000 |
| | 5720411127 | 30PC bottle (B3) ass'y | 2.6 X 9.7 | 26.0 | --- | --- | --- | --- | 30,000 | 109,000 |
| | 5720411124 | 30PP bottle (C) ass'y ^{*2} | 2.6 X 9.7 | 26.0 | --- | --- | --- | --- | 41,000 | 204,000 |
| | 5720411123 | 30PP bottle ass'y | 2.6 X 9.7 | 26.0 | --- | --- | --- | --- | 28,000 | 95,000 |
| | 5720411121 | 30PP seal tube | 2.6 X 7.5 | 31.7 | 5720410250 | E40-30 float spacer | 5720411256 | E space cap | 50,000 | 303,000 |
| 20 | 5720411116 | 20PP seal tube | 2.6 X 5.5 | 22.2 | 5720411341 | E40-20 float spacer | 5720411256 | E space cap | 50,000 | 303,000 |
| | 5720411020 | 12PP tube | 1.6 X 7.6 | 10.5 | 5720411185 | 12A40 adapter | 5720411208 | C-AL cap ass'y | 40,000 | 183,000 |
| 12 | 5720411021 | 12PE tube | 1.6 X 7.5 | 10.5 | 5720411185 | 12A40 adapter | 5720411208 | C-AL cap ass'y | 40,000 | 183,000 |
| | 5720411022 | 12PC tube | 1.6 X 7.6 | 10.5 | 5720411185 | 12A40 adapter | 5720411208 | C-AL cap ass'y | 40,000 | 183,000 |
| | 5720411106 | 12PET tube | 1.6 X 7.6 | 11.4 | 5720411185 | 12A40 adapter | 5720411209 | C4-AL cap ass'y | 40,000 | 183,000 |
| | | | | 10.5 | 5720411185 | 12A40 adapter | 5720411208 | C-AL cap ass'y | 34,000 | 132,000 |
| | 5720411023 | 12SST tube | 1.6 X 7.6 | 10.5 | 5720411185 | 12A40 adapter | 5720411210 | C-SST cap ass'y | 30,000 | 103,000 |
| | | | | 8.3 | 5720411185 | 12A40 adapter | --- | --- | 40,000 | 183,000 |
| 10 | --- | 10PC bottle (B) ass'y | 1.6 X 8.2 | 8.4 | 5720411185 | 12A40 adapter | --- | --- | 40,000 | 183,000 |
| | 5720411024 | 10PC thick-walled tube | 1.6 X 7.7 | 6.3 | 5720411185 | 12A40 adapter | --- | --- | 40,000 | 183,000 |
| | 5720411027 | 10PC bottle ass'y | 1.6 X 8.2 | 8.0 | 5720411185 | 12A40 adapter | --- | --- | 40,000 | 183,000 |
| | 5720411028 | 10PP bottle ass'y | 1.6 X 8.0 | 8.0 | 5720411185 | 12A40 adapter | --- | --- | 35,000 | 140,000 |
| | 5720411025 | 10PP thick-walled tube | 1.6 X 7.6 | 6.5 | 5720411185 | 12A40 adapter | --- | --- | 30,000 | 103,000 |
| | 5720411029 | 6.5PP tube | 1.3 X 6.4 | 5.8 | 5720411184 | 6.5A40 adapter | 5720411201 | B-AL cap ass'y | 40,000 | 173,000 |
| 6.5 | 5720411030 | 6.5PC tube | 1.3 X 6.4 | 5.8 | 5720411184 | 6.5A40 adapter | 5720411201 | B-AL cap ass'y | 40,000 | 173,000 |
| | 5720411080 | 6.5PET tube | 1.3 X 6.4 | 6.2 | 5720411184 | 6.5A40 adapter | 5720411207 | B2-AL cap ass'y | 40,000 | 173,000 |
| 5 | 5720411031 | 5PP tube | 1.3 X 5.2 | 5 | 5720411183 | 5A40 adapter | 5720411201 | B-AL cap ass'y | 40,000 | 181,000 |
| 4.7 | 5720411032 | 4.7PC thick-walled tube | 1.3 X 6.4 | 3.4 | 5720411184 | 6.5A40 adapter | --- | --- | 40,000 | 173,000 |
| 4 | 5720411033 | 4PP tube | 1.1 X 6.0 | 4 | 5720411182 | 4A40 adapter | 5720411199 | A3-AL cap ass'y | 40,000 | 180,000 |
| 3 | 5720411012 | 3PC tube | 1.3 X 5.1 | 2.5 | 5720411183 | 5A40 adapter | --- | --- | 40,000 | 179,000 |

※The material of tube adapters: Polyacetal

※The material of float spacers: Modified polyphenylene ether

* 1: When using tubes with S-40AL cap ass'y, tubes can not be reused.

* 2: Use a 30PP bottle (C) at a temperature between 4°C and 20°C.

CAUTION :

Chemical resistance and the strength of tubes/bottles varies with speed, temperature, and so on. Before using sample, fill tubes/bottles with water, buffer solution, etc. instead of sample and run them at the intended speed, temperature, and so on to ensure that there is no abnormality.


When using 40PP tubes, 40PP seal tubes, 30PP seal tubes, and 20PP seal tubes, we recommend the purchase of the tube kit, what is the set of tubes and necessary tools. Refer to Chapter 6 for the tube kit.

Refer to "ROTORS, TUBES, BOTTLES AND CAPS (Part No.S999204)" for the cleaning and sterilization of the tubes/bottles.

3. How to Use

How to use tubes/bottles varies in the kinds of tubes/bottles. The brief procedure of tubes/bottles handling is described as follows.

For the detail about handling tubes/bottles, refer to “ROTORS, TUBES, BOTTLES AND CAPS (Part No.S999204)” attached to the centrifuge.

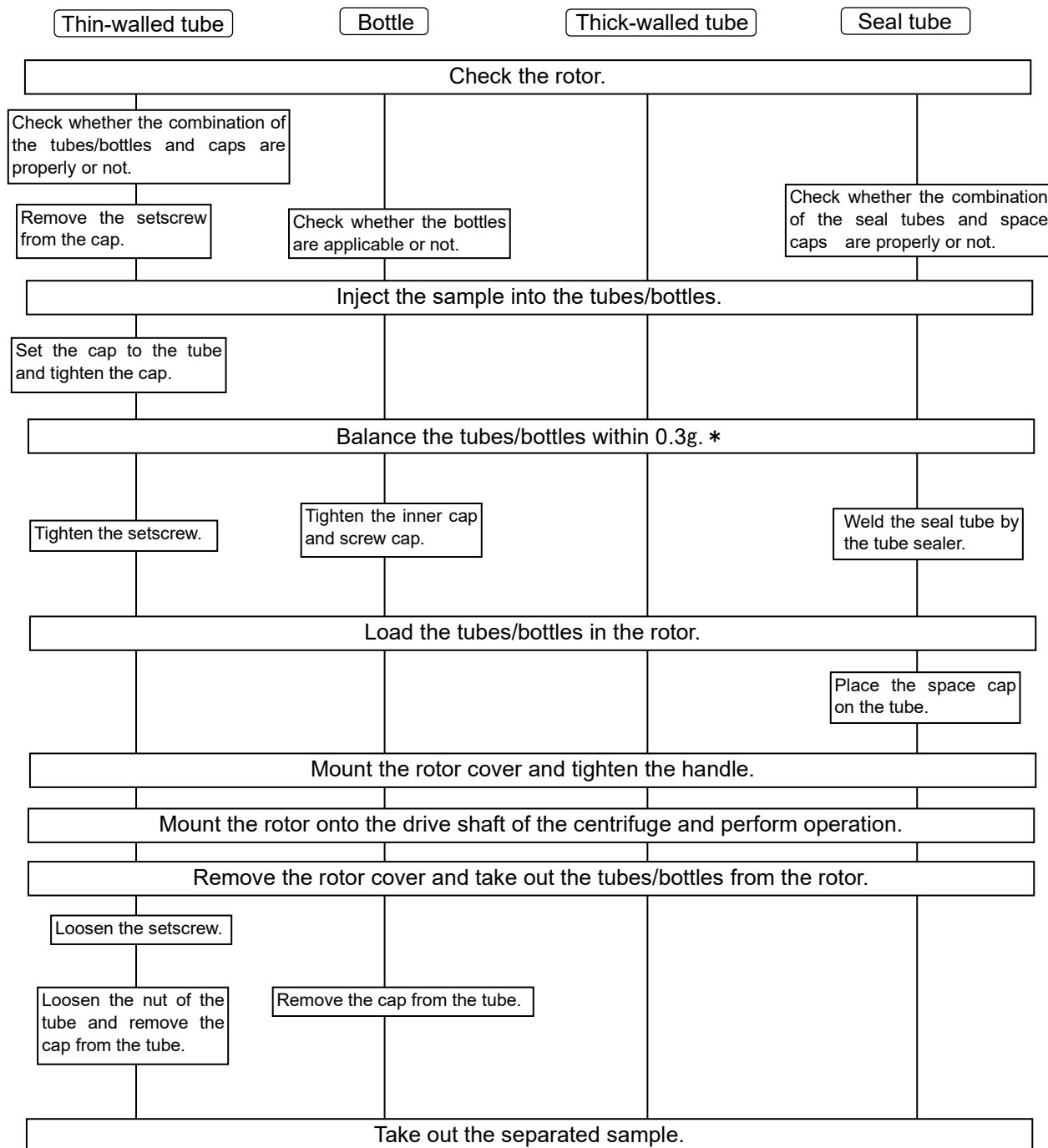(*) When using CP-NX series, CP-WX series, CP-MX series, or CP- α/β series, approximately even quantities of sample in the tubes are sufficient for balancing, and extremely different sample quantities must be avoided.

(The levels of samples in the tubes should be approximately equal. Their difference should be within 5 mm.)

For some kinds of tubes, if a small quantity of a sample is injected to them, the rotor can not be used at the allowable maximum speed. Tube deformation may occur due to a small quantity of a sample during operation.

=For handling=

● Allowable speed of the rotor

1. To centrifuge a sample of an average density more than 1.2 g/mL.
Determine the allowable speed from the following equation.

$$\text{Allowable speed (rpm)} = 50,000 \text{ (rpm)} \times \sqrt{\frac{1.2 \text{ (g / mL)}}{\text{Mean density of sample (g/mL)}}}$$

2. The allowable speed may be lower depending on the density of a sample when using of salt-such as cesium chloride. Refer to “5. Using Cesium Chloride” for the detail.
3. The allowable speed may be lower depending on the combination of tubes/bottles and tube adapter, etc. Refer to “2. Applicable Tubes/Bottles”

● Tubes/Bottles

Before operation, check whether the combination of tubes/bottles, tube adapter, etc. is properly or not. When using seal tubes and thin-walled tubes, fill them with a sample. When using bottles and thick-walled tubes, use them with optional volume less than net capacity. However, when using the 30PP bottle (C) and the 30PC bottle (C3) at 100,000xg or more, fill it with the sample until the sample level is more than the shoulder line of them.

Tighten the cap securely. Weld the seal tube completely according to the instruction manual of the STF-1, the STF2, or the STF3 tube sealer.

Refer to “2. Applicable Tubes/Bottles” in this manual and “ROTORS, TUBES, BOTTLES AND CAPS (Part No. S999204)”.

● Handling the rotor

WARNING :

- Never exceed the maximum speed of the rotor (mentioned on the rotor). Otherwise, rotor may be damaged and result in damage to the centrifuge.
- Check the chemical resistance chart attached the rotor, and do not use any sample inapplicable to the rotor. Using such a sample could corrode the rotor.

CAUTION :

- Do not load only one tube or load tubes asymmetrically : Asymmetrically loading may cause imbalance operation and damage the centrifuge and rotor.
- Balance the tubes/bottles and the samples including the caps and the tube adapters within the allowable imbalance of the rotor.
Do not exceed the allowable imbalance. Otherwise, the centrifuge or the rotor may be damaged due to excessive imbalance operation.
- If foreign matter adheres to the rotor, wipe it off promptly. Otherwise, the rotor may be corroded.
- Be sure to attach the rotor cover to the rotor body and tighten the handle securely.
Tightening the handle insufficiently or removing the handle from the rotor can cause the rotor cover to come off. And it may result in the damage to the centrifuge and the rotor.
- Inspect and maintain the rotor after use. If abnormality is observed, do not use it.
Contact an authorized sales or service representative.

● Sample limitation

WARNING :

Never use any material capable of producing flammable or explosive vapors.
Your centrifuge and rotor are not designed to confine any sample particles dispersed due to leakage. Therefore, when using toxic or radioactive samples or pathogenic or infectious blood samples, make sure to prepare necessary safety measures at your own responsibility.

● Rotor life management of the rotor with optical adapter

CAUTION :

Each time the rotor with optical adapter is used, be sure to enter the results in the “rotor log book”. Failure to do so will disable management of the life of rotor.
No entry is need for RLM rotor.

● RLM adapter and Optical adapter

⚠ WARNING :

- Do not remove the RLM adapter or optical adapter from rotor, or replace it with the adapter for another rotor. The adapter is a critical component that detects the over-speed of rotor : If an adapter that is compatible with the rotor is attached, the rotor could break, resulting in damage to the ultracentrifuge.

⚠ CAUTION :

- Do not bring the RLM adapter near a magnet, or scratch it: Doing so will erase the memory stored in adapter, and make the rotor unusable. To prevent the adapter from being scratched, store the rotor with RLM adapter, using the stand provided with the rotor (rotor stand for protecting adapter). If the adapter is scratched, immediately stop using rotor and contact an authorized sales or service representative.
- Take great care not to scratch the optical adapter: Any scratch on adapter will make the rotor unusable. To prevent the adapter from being scratched, store the rotor with optical adapter, using the stand provided with the rotor (rotor stand for protecting adapter). If the adapter is scratched, immediately stop using rotor and contact an authorized sales or service representative.
- If foreign matter adheres to the RLM adapter or optical adapter, wipe it off promptly. Otherwise, the RLM adapter or optical adapter may be corroded.

4. Maintenance


● Maintenance of the rotor

⚠ WARNING :

Do not allow the rotor temperature to rise over 100°C. Otherwise, the rotor can be brittle.

⚠ CAUTION :

Use a neutral detergent having a pH between 5 and 9. Otherwise, the rotor can be corroded.

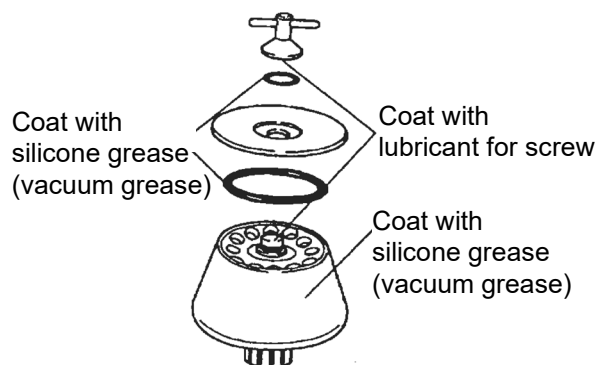Wash the rotor with tap water or a dilute solution of neutral detergent and rinse it out with distilled water. Turn the rotor upside down with the rotor removed to dry it well. Check the rotor is completely dried and then put the light coat of silicone grease (vacuum grease) on the surface of the rotor, the light coat of lubricant for screw on the thread portion, the light coat of silicone grease (vacuum grease) on the packing, Store them in a dry place. If foreign matter adheres to the rotor, soak the rotor in warm water in one or two hours and remove foreign matter with a soft brush etc. Use a neutral detergent having a pH between 5 and 9. If you cannot remove it, contact an authorized sales or service representative.

● Sterilizing the rotor

⚠ WARNING :

Never sterilize the rotor by autoclaving or boiling. Otherwise the strength of rotor may be significantly decreased.

⚠ CAUTION :

Do not sterilize the tube adapters made of plastic by autoclaving or boiling. Otherwise the tube adapters may be deformed. Such the tube adapters can not be used.

Sterilize the rotor according to either gas sterilization method (ethylene oxide or formaldehyde) or chemical sterilization method (70% ethanol, 3% hydrogen peroxide, 3% formalin).

Sterilize the tube adapters made of plastic according to either gas sterilization method (ethylene oxide or formaldehyde) or chemical sterilization method (70% ethanol, 3% hydrogen peroxide, 3% formalin).

* Do not dip the rotor and the tube adapters in the formalin (3%) solution more than 2 hours.

● Inspecting the rotor

Check the rotor after every run as follows.

- The thread portion is worn. Contact an authorized sales or service representative.
- The tube cavity is corroded. Contact an authorized sales or service representative.
- The packing is deteriorated. Replace it with the new one.
- The adapter is stained. Wipe it out with a soft cloth.

Check the rotor once a month as follows.

- The inside of drive hole (crown hole) is stained. . . . Wipe it out with a soft cloth dampened with water.

CAUTION :

Clean the inside of the drive hole (crown hole) of the rotor and the surface of the drive shaft (crown) of the centrifuge once a month. If the drive hole or the drive shaft is stained or any foreign matter is adhered, the rotor may be improperly installed and come off during operation.

● Life of the rotor

WARNING :

Do not use rotors that have exceeded their lifetime; to do so might cause the rotor failure and the centrifuge might be damaged.

While using rotor repeatedly, its strength decreases gradually due to fatigue and creep of material, by which the rotor life is determined.

This rotor must not be used when they reached 5,000 operation times or 10,000 accumulated hours. If you use the RLM rotor at a speed lower than its maximum permissible speed, the life of the RLM rotor will be prolonged automatically by calculating from the strength of the rotor material.

5. Using Cesium Chloride

The cesium chloride (CsCl) solution is frequently used as a density gradient medium, but if the CsCl solution with a high density is used, it may saturate during the rotation of the rotor depending on the rotation condition which may cause CsCl crystals to be educed. Educated crystals have a high density (approx. 4g/mL) and apply an excessive load to the rotor, which is very dangerous.

The crystallization will also greatly affect the density gradient of the CsCl solution and separation status of the sample. Therefore, always use a solution which will not crystallize the allowable speed.

Fig.5-1 shows the relationship between the maximum allowable density and the speed at which cesium chloride does not crystallize in the condition that standard tube is filled with the cesium chloride solution. In this diagram, crystallization will occur above the curves.

Be sure to select the density of the cesium chloride solution and the rotor speed at which cesium chloride does not crystallize. The curve in this diagram varies in temperature etc. Before actual operation, centrifuge the cesium chloride solution with the same density to ensure that crystallization will not occur.

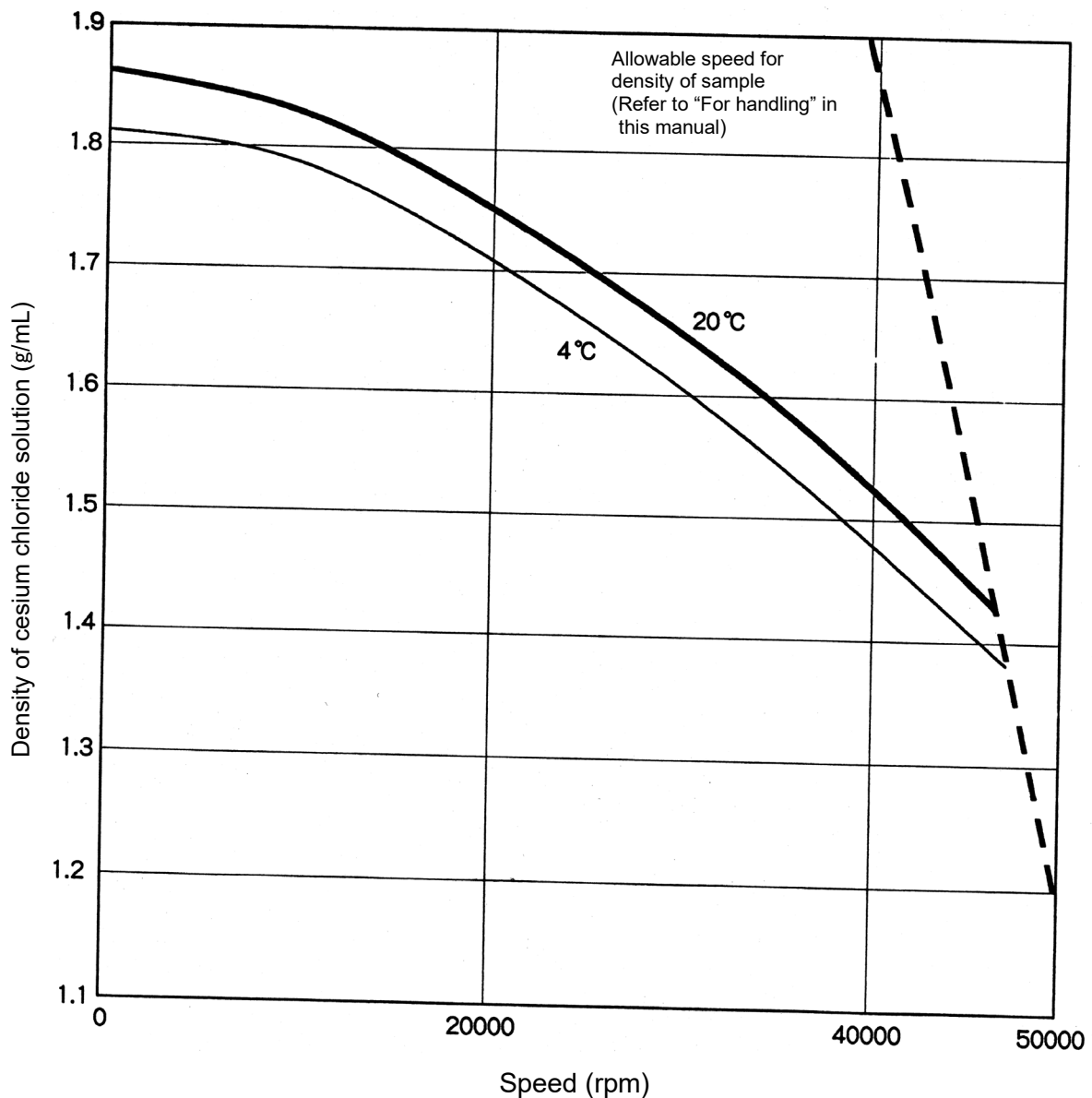
Fig.5-1 Relationship Between Density of Cesium Chloride and Speed

6. Packing List

Check the parts and accessories of rotor when receiving the rotor.
Contact an authorized sales or service representative when the need arises.

| No. | Name | Qty | Remark | Order no. (Qty) | No. | Name | Qty | Order no. (Qty) |
|-----|--------------------------------------|-----|-----------|--------------------|-----|------------------------------------|-----|--------------------|
| ① | Rotor | 1 | | | ⑩ | Rotor stand | 1 | 5720410232 (1) |
| ② | Handle | 1*1 | | | ⑪ | Lubricant for screw | 1 | 5720411230 (1) |
| ③ | Packing (for the handle) (O-Ring) | 1*1 | | 5720410154 (5) | ⑫ | Silicone grease (vacuum grease) | 1 | 5720850317 (1) |
| ④ | Cover | 1*1 | | | ⑬ | Tool box (C) | 1 | |
| ⑤ | Packing (for the rotor) (O-Ring) | 1*1 | | 5720410160 (3) | ⑭ | | | |
| ⑥ | | | | | ⑮ | | | |
| ⑦ | | | | | ⑯ | | | |
| ⑧ | Packing (for the handle) (O-Ring) | 2 | For spare | | ⑰ | Rotor log book *2 | 1 | 5720411287 (1) |
| ⑨ | Packing (for the rotor) (O-Ring) | 2 | For spare | | ⑱ | Instruction manual | 1 | |

*1: the quantities of the parts that are mounted to the rotor

*2: The "rotor log book" is not included in RLM rotor.

Tubes are not included in this rotor. Please purchase the applicable tubes.

When using 40PP tubes, 40PP seal tubes, 30PP seal tubes, and 20PP seal tubes, we recommend the purchase of under-mentioned tube kit, what is the set of tubes and necessary tools.

●40PP tube kit (B) (Order no.5720411144)

(component)

| | | | |
|------------|-----------------------------------|----|----------------|
| 5720411139 | 40PP tube | 2 | (50 pcs./pack) |
| 5720411212 | E3-AL Cap ass'y | 12 | |
| --- | Packing (for the cap) (O-Ring) | 2 | (10 pcs./pack) |
| 5720411365 | Tube vise(E) | 1 | |
| 5720411282 | Torque wrench | 1 | |
| 5720411283 | Socket adaptor | 1 | |
| 5720411410 | 17mm socket | 1 | |
| 5720411267 | Tube setter | 1 | |

●40PP seal tube kit (A) (Order no.5720411143)

(component)

| | | | |
|------------|----------------|----|----------------|
| 5720411141 | 40PP seal tube | 2 | (50 pcs./pack) |
| 5720411256 | E-space cap | 12 | |
| 5720411232 | Tube setter | 1 | |

※When using 40PP seal tubes, Tube sealer (STF3,STF2,STF-1) and Tube rack (E) (Order no.5720410008) are needed.

●30PP seal tube kit (Order no.5720411122)

(component)

| | | | |
|------------|---------------------|---|----------------|
| 5720411121 | 30PP seal tube | 2 | (50 pcs./pack) |
| 5720411256 | E- space cap | 4 | |
| 5720410250 | E40-30 float spacer | 4 | |
| 5720411232 | Tube setter | 1 | |
| 5720411267 | Tube setter (B) | 1 | |


※ When using 30PP seal tubes, Tube sealer (STF3,STF2) and Tube rack (E3) (Order no. 5720410015) are needed.

●20PP seal tube kit (Order no.5720411117)

(component)

| | | | |
|------------|---------------------|---|----------------|
| 5720411116 | 20PP seal tube | 2 | (50 pcs./pack) |
| 5720411256 | E- space cap | 4 | |
| 5720411341 | E40-20 float spacer | 4 | |
| 5720411232 | Tube setter | 1 | |
| 5720411267 | Tube setter (B) | 1 | |

※ When using 20PP seal tubes, Tube sealer (STF3,STF2) and Tube rack (E4) (Order no. 5720410016) are needed.

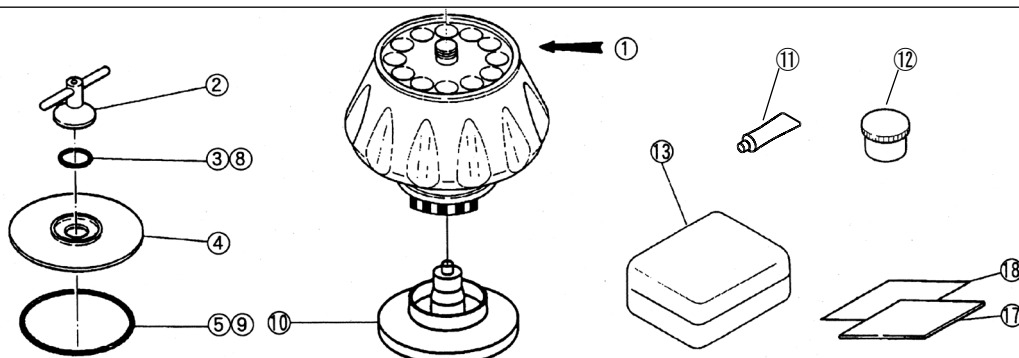7. Decontamination

⚠ WARNING:

- If the centrifuge, rotor or an accessory is contaminated by samples that are toxic or radioactive, or blood samples that are pathogenic or infectious, be sure to decontaminate the item according to good laboratory procedures and methods.
- If there is a possibility that the rotor or an accessory is contaminated by samples that might impair human health (for example, samples that are toxic or radioactive, or blood samples that are pathogenic or infectious), it is your responsibility to sterilize or decontaminate the rotor or the accessory properly before requesting repairs from an authorized sales or service representative. Note that we cannot repair the centrifuge, rotor or the accessory unless sterilization or decontamination is completed.
- It is your responsibility to sterilize and/or decontaminate the rotor or parts properly before returning them to an authorized sales or service representative. In such cases, copy the attached decontamination sheet and fill out the copied sheet, then attach it to the item to be returned. We may ask you about the treatment for the rotor or parts if the decontamination is checked and judged as insufficient by us. It is your responsibility to bear the cost of sterilization or decontamination. Note that we cannot repair or inspect the rotor or the accessory unless sterilization or decontamination is completed.

8. Rotor Retirement

After many years of use, there will be inevitably some corrosion or stress corrosion. At some points, the combination of such damage and metal fatigue could make the rotor vulnerable to a failure. Although a rotor may appear to be in a good condition, you should follow the rotor retirement recommendation shown below.

| Rotor | Material | Retire After Years |
|-----------------------|----------------------------------|--------------------|
| Angle rotor | Titanium alloy | 12 |
| Vertical rotor | Aluminum alloy | 10 |
| Swing rotor | Titanium alloy Aluminum alloy | 10 |
| Zonal rotor | Titanium alloy | 10 |
| Continuous flow rotor | Titanium alloy | 10 |

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Copyright © 2025 Eppendorf Himac Technologies Co., Ltd. All rights reserved.

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