Keep this rotor instruction manual and the centrifuge manual in the file.

Angle Rotors for Ultracentrifuge

P80AT

Before using these rotors, please carefully read this instruction manual, "ROTORS, TUBES, BOTTLES AND CAPS", and the centrifuge instruction manual for its efficient operation and for your safety.

Keep this instruction manual for your reference and refer to it as required.

Contents

1. Specifications	_____	1
2. Applicable tubes/bottles	_____	2
3. How to use	_____	3
4. Maintenance	_____	5
5. Using cesium chloride	_____	7
6. Packing list	_____	8
7. Decontamination	_____	9
8. Rotor retirement	_____	9

Part No. { RLM rotor
9123092M
Rotor with optical adapter
91230920
Rotor with optical disk
9123093C

2022.01
S99924016
Original instruction manual

Safety Reminder

- Centrifuge rotors rotating at high speed have considerable potential for damage to personal properties if used improperly.
For safe and proper use of this rotor, carefully read the centrifuge instruction manual and this rotor instruction manual before use and observe the instructions.

 **WARNING :** and  **CAUTION :** notes are used to call your attention in this manual to prevent personal injury or damage to the rotor and the centrifuge. These notes are defined as follows.

 **WARNING:** indicates a potentially hazardous situation which, if not avoided, could result in personal severe injury or possible death.

 **CAUTION:** indicates a hazardous situation which, if not avoided, could result in personal injury or severe damage to the instrument.

WARNING

- Never use any material capable of producing flammable or explosive vapors. Your centrifuge and rotor have no explosion-proof construction.
- Your centrifuge and rotor are not designed to confine any sample particles dispersed due to leakage. Therefore, when using toxic or radioactive samples or pathogenic or infectious blood samples, make sure to prepare necessary safety measures at your own responsibility.
- Never exceed the maximum speed of the rotor (mentioned on the rotor).
Always reduce rotor speed as instructed in this manual when rotor speed is limited due to sample density or kinds of tubes.
- Check the chemical resistance chart attached the rotor, and do not use any sample inapplicable to the rotor. Using such a sample could corrode the rotor.
- Do not remove the RLM adapter or optical adapter/disk from rotor, or replace it with the adapter/disk for another rotor. The adapter/disk is critical component that detects the over-speed of rotor. If an adapter/disk that is not compatible with the rotor is attached, the rotor could break, resulting in damage to the centrifuge.
- Do not allow the rotor temperature to rise 100°C.: This would cause the material brittle.
- Never sterilize the rotor by autoclaving or boiling. Otherwise the strength of rotor may be significantly decreased.
- If the centrifuge, rotor, or an accessory is contaminated by samples that toxic or radioactive, or blood samples that pathogenic or infectious, be sure to decontaminate the item according to good laboratory procedures and methods.
- If there is a possibility that the centrifuge, rotor, or an accessory is contaminated by samples that might impair human health (for example, samples that are toxic or radioactive, or blood samples that are pathogenic or infectious), it is your responsibility to sterilize or decontaminate the centrifuge, rotor, or the accessory properly before requesting repairs from an authorized sales or service representative.
- It is your responsibility to sterilize and/or decontaminate the centrifuge, rotor, or parts properly before returning them to an authorized sales or service representative.

CAUTION

- Check the chemical resistance chart attached to the rotor, and do not use any sample inapplicable to the tubes/bottles, the tube adapter, or the cap for the tubes/bottles. Using such a sample could deteriorate them.
- The allowable speed may be lower depending on the density of sample, use of salt-such as cesium chloride or the combination of tubes or bottles and tube adapters and caps. Refer to this instruction manual.
- Always balance the tubes/bottles and samples within the allowable imbalance of rotor, including the caps and the tube adapters. Do not exceed the allowable imbalance. Otherwise the rotor and the centrifuge may be damaged.
- Do not load only one tube or load tubes asymmetrically : Asymmetrically loading may cause imbalance operation and damage the centrifuge and rotor.
- Clean the inside of the drive hole (crown hole) of the rotor and the surface of the drive shaft (crown) of the centrifuge once a month. If the drive hole or the drive shaft is stained or any foreign matter is adhered, the rotor may be improperly installed and come off during operation.
- Use the rotor tubes and bottles within their actual capacities.
- Do not use tubes/bottles that have exceeded their life expectancy. Failure to do so could result in damage of tubes/bottles and the rotor and the centrifuge. The life expectancy of tubes/bottles depends on factors such as the characteristics of samples, speed of the rotor used, and temperature. Always check for deterioration and damage (cracks, deformation, and so on) on tubes/bottles before using them. Do not use the tubes/bottles if you find such a problem.
- Inspect and maintain the rotor after use. If abnormality is observed, do not use it.
Contact an authorized sales or service representative.

1. Specifications

1.1 Construction

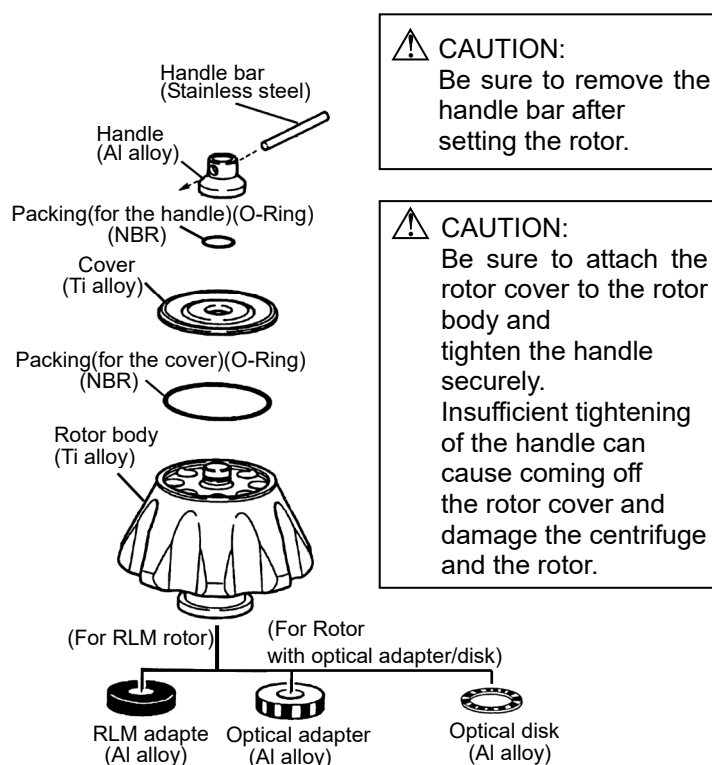
Fig.1-1. Construction of P80AT rotor

1.2 Specifications

- Max. speed: 80,000 rpm
- Max. RCF: 615,000 xg
- Allowable density at maximum speed of the rotor: 1.2 g/mL
- K Factor at maximum speed of the rotor: 27
- Nominal rotor capacity: 12 mL × 8 tubes = 96 mL
- Material of rotor: Ti alloy
- Tube size: ϕ 1.6 × 7.6 cm
- Weight: 6.0 kg
- Max. diameter: 18.1 cm
- Max. height: 16.4 cm

⚠ CAUTION:

- Do not use a RLM rotor with the centrifuge other than a series of a CP-NX/WX/MX/ α ultracentrifuge.
- Do not use rotor with optical adapter/disk with model 55P and 40P (old type centrifuge).

1.3 Cross-sectional view of rotor

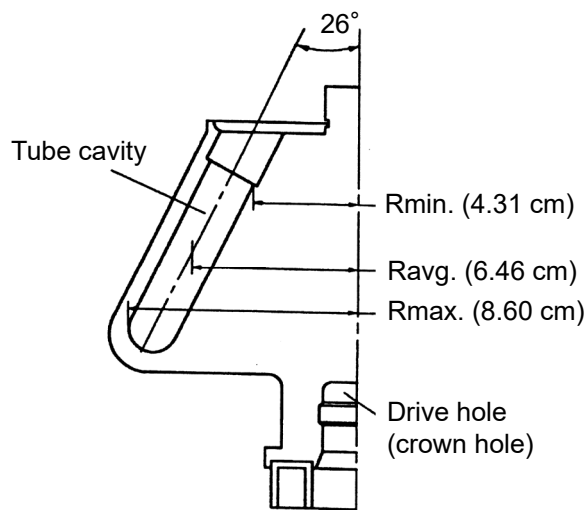
Fig.1-2. Cross-sectional view of P80AT rotor

⚠ CAUTION:

Clean the inside of the drive hole (crown hole) of the rotor and the surface of the drive shaft (crown) of the centrifuge once a month. If the drive hole or the drive shaft is stained or any foreign matter is adhered, the rotor may be improperly installed and come off during operation.

1.4 Characteristic of rotor

Table 1-1. Characteristic of P80AT

Rotor Speed (rpm)	RCF(xg)			K factor
	Rmax.	Ravg.	Rmin.	
10,000	9,620	7,220	4,820	1,750
20,000	38,500	28,900	19,300	437
30,000	86,500	65,000	43,400	194
40,000	154,000	116,000	77,100	109
50,000	240,000	181,000	120,000	70
60,000	346,000	260,000	173,000	49
70,000	471,000	354,000	236,000	36
80,000	615,000	462,000	308,000	27

※ For detail about K factor, refer to “ROTORS, TUBES, BOTTLES AND CAPS (Part No.S999204)” attached to the centrifuge.

Calculation formula

$$RCF = 1.118 \times 10^{-5} \times R \times N^2 (\times g)$$

R: Rotating radius (cm)

N: Speed (rpm)

2. Applicable tubes/bottles

2.1 Applicable tubes/bottles

Table 2-1. Applicable tubes/bottles

Tubes/Bottles					Tube adapter Float spacer		Cap		Max. speed (rpm)	Max. RCF (xg)
Nominal capacity (mL)	Part No.	Name	Size (ϕ X L cm)	Actual capacity (mL/tube)	Part No.	Name	Part No.	Name		
12	345320A	12PA Seal Tube	1.6 X 7.8	13.5	---	---	486582	C2 Space Cap	80,000	615,000
	329606A	12PA Tube	1.6 X 7.6	10.5	---	---	463577A	C-TI Cap Ass'y	80,000 ^{*1}	615,000
				10.2	---	---	S410542A	S-12AL Cap Ass'y ^{*2}	80,000	615,000
	325470A	12PE Tube	1.6 X 7.5	10.5	---	---	463577A	C-TI Cap Ass'y	60,000	346,000
				10.0	---	---	S410542A	S-12AL Cap Ass'y ^{*2}	60,000	346,000
	325751A	12PC Tube	1.6 X 7.6	10.5	---	---	463577A	C-TI Cap Ass'y	60,000	346,000
	S413604A	12TI Tube	1.6 X 7.6	10.5	---	---	463577A	C-TI Cap Ass'y	37,000	132,000
8.5				---	---	---	---	46,000	203,000	
326010A	12SST Tube	1.6 X 7.6	10.5	---	---	463577A	C-TI Cap Ass'y	32,000	98,500	
			8.5	---	---	---	---	38,000	139,000	
10	S303922A	10PC Bottle (B) Ass'y	1.6 X 8.2	8.4	---	---	---	---	62,000	370,000
	338455A	10PC Thick-Walled Tube	1.6 X 7.7	6.5	---	---	---	---	47,000	212,000
	325952A	10PC Bottle Ass'y	1.6 X 8.2	8.0	---	---	---	---	38,000	139,000
	334105A	10PA Bottle Ass'y	1.6 X 8.2	8.0	---	---	---	---	33,000	105,000
	338456A	10PA Thick-Walled Tube	1.6 X 7.6	6.7	---	---	---	---	28,000	75,400
8	S408831A	8PA Seal Tube	1.6 X 6.1	8.3	S413976	C12-8 Float Spacer	486582	C2 Space Cap	80,000	615,000
6.5	329445A	6.5PA Tube	1.3 X 6.4	5.8	336687A	6.5A12 Adapter	414429A	B-AL Cap Ass'y	38,000	124,000
	326415A	6.5PC Tube	1.3 X 6.4	5.8	336687A	6.5A12 Adapter	414429A	B-AL Cap Ass'y	38,000	124,000
6	S311555A	6PA Seal Tube	1.6 X 5.0	6.5	S413977	C12-6 Float Spacer	486582	C2 Space Cap	80,000	615,000
5	332245A	5PA Tube	1.3 X 5.2	5.0	336686A	5A12 Adapter	414429A	B-AL Cap Ass'y	38,000	115,000
4.7	S303141A	4.7PC Thick-Walled Tube	1.3 X 6.4	3.4	336687A	6.5A12 Adapter	---	---	38,000	124,000
4	329561A	4PA Tube	1.1 X 6.0	4.0	336685A	4A12 Adapter	336711A	A3-AL Cap Ass'y	38,000	132,000
3	336708A	3PA Tube	1.1 X 4.7	3.0	336684A	3A12 Adapter	336711A	A3-AL Cap Ass'y	38,000	123,000
	S301599A	3PC Tube	1.3 X 5.1	2.6	336686A	5A12 Adapter	---	---	38,000	115,000
0.9	S304296A	0.9PC Thick-Walled Tube	0.8 X 5.0	0.9	336683A	2A12 Adapter	---	---	42,000	151,000

*1: When using 12PA Tubes with C-TI Cap Ass'y at speed of 70,000 rpm and more than 70,000 rpm, 12PA Tubes can not be reused.

*2: When using tubes with S-12AL Cap Ass'y, tubes can not be reused.

1. When using 12PA Tubes, 12PA Seal Tubes, 8PA Seal Tubes, and 6PA Seal Tubes, we recommend the purchase of the tube kit, what is the set of tubes and necessary tools.
Refer to Chapter 6 for the tube kit.

2. Material of tube adapters: Polyacetal

3. Material of float spacers: Modified polyphenylene ether

CAUTION :


Chemical resistance and the strength of tubes/bottles varies with speed, temperature, and so on. Before using sample, fill tubes/bottles with water, buffer solution, etc. instead of sample and run them at the intended speed, temperature, and so on to ensure that there is no abnormality.

Refer to "ROTORS, TUBES, BOTTLES AND CAPS (Part No.S999204)" for the cleaning and sterilization of the tubes/bottles.

3. How to use

How to use tubes/bottles varies in the kinds of tubes/bottles. The brief procedure of tubes/bottles handling is described as follows.

For the detail about handling tubes/bottles, refer to "ROTORS, TUBES, BOTTLES AND CAPS (Part No.S999204)" attached to the centrifuge.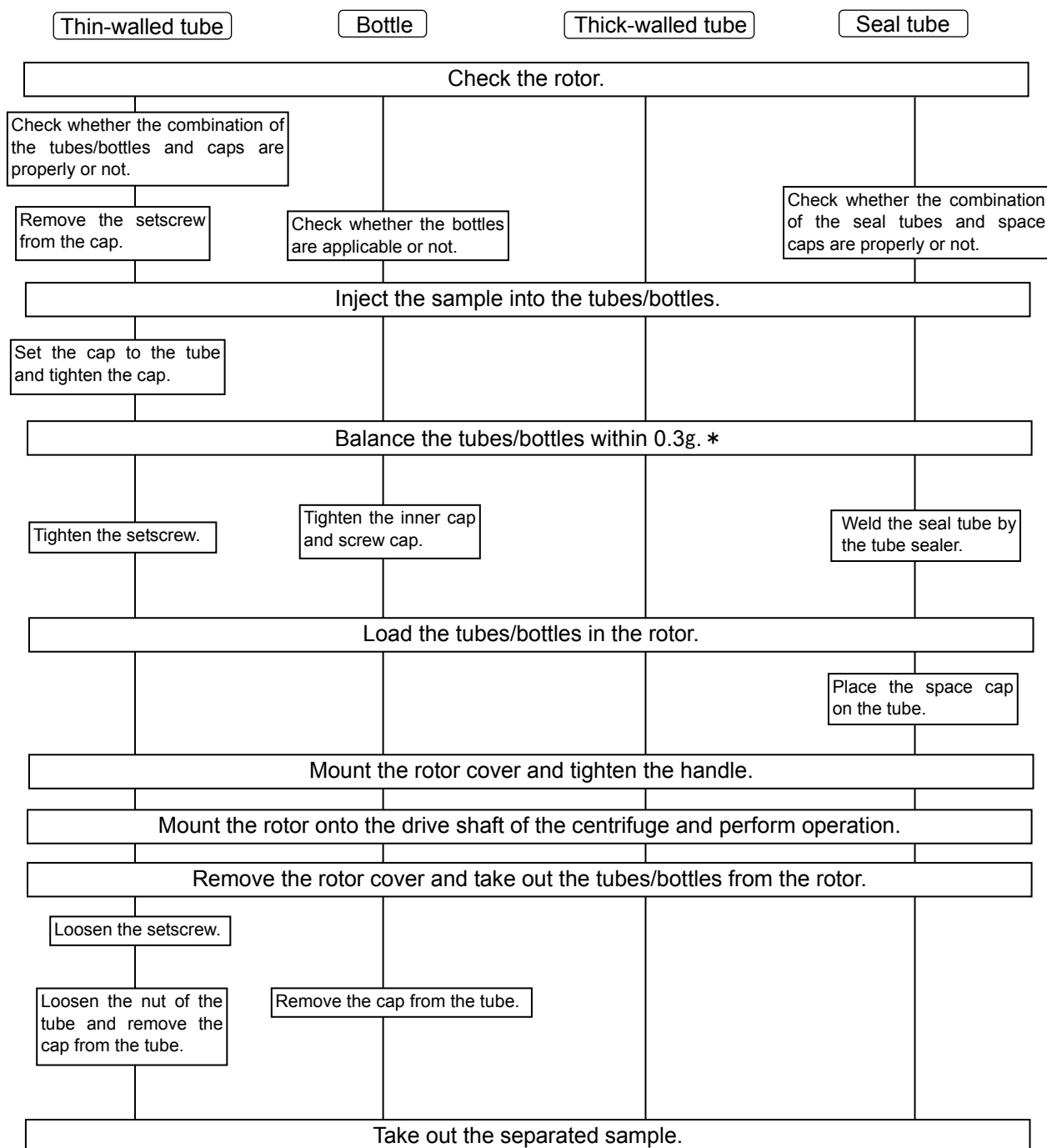

(*) When using CP-NX series, CP-WX series, CP-MX series, or CP- α/β series, approximately even quantities of sample in the tubes are sufficient for balancing, and extremely different sample quantities must be avoided.

(The levels of samples in the tubes should be approximately equal. Their difference should be within 5 mm.) For some kinds of tubes, if a small quantity of a sample is injected to them, the rotor can not be used at the allowable maximum speed. Tube deformation may occur due to a small quantity of a sample during operation.

=For handling=

● Allowable speed of rotors

1. To centrifuge a sample of an average density more than 1.2 g/mL.

Determine the allowable speed from the following equation.

$$\text{Allowable speed (rpm)} = 80,000 \text{ (rpm)} \times \sqrt{\frac{1.2 \text{ (g/mL)}}{\text{Mean density of sample (g/mL)}}}$$

2. The allowable speed may be lower depending on the density of a sample when using of salt-such as cesium chloride. Refer to “5. Using cesium chloride” for the detail.

3. The allowable speed may be lower depending on the combination of tubes/bottles and tube adapter, etc. Refer to “2. Applicable tubes/bottles”

● Tube/Bottle

Before operation, check whether the combination of tubes/bottles, tube adapter, etc. is properly or not. When using seal tubes and thin-walled tubes, fill them with a sample. When using bottles and thick-walled tubes, use them with optional volume less than net capacity. Tighten the cap securely. Weld the seal tube completely according to the instruction manual of the STF-1, the STF2, or the STF3 tube sealer.

Refer to “2. Applicable tubes/bottles” in this manual and “ROTORS, TUBES, BOTTLES AND CAPS (Part No. S999204)”.


● Handling the rotor

⚠ WARNING :

- Never exceed the maximum speed of the rotor (mentioned on the rotor). Otherwise, rotor may be damaged and result in damage to the centrifuge.
- Check the chemical resistance chart attached the rotor, and do not use any sample inapplicable to the rotor. Using such a sample could corrode the rotor.

⚠ CAUTION :

- Do not load only one tube or load tubes asymmetrically : Asymmetrically loading may cause imbalance operation and damage the centrifuge and rotor.
- Balance the tubes/bottles and the samples including the caps and the tube adapters within the allowable imbalance of the rotor.
Do not exceed the allowable imbalance. Otherwise, the centrifuge or the rotor may be damaged due to excessive imbalance operation.
- If foreign matter adheres to the rotor, wipe it off promptly. Otherwise, the rotor may be corroded.
- Be sure to attach the rotor cover to the rotor body and tighten the handle securely.
Insufficient tightening of the handle can cause coming off the rotor cover and damage the centrifuge and the rotor.
- Be sure to remove the handle bar after setting the rotor.
- Inspect and maintain the rotor after use. If abnormality is observed, do not use it.
Contact an authorized sales or service representative.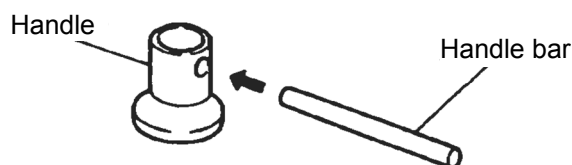

Insert the handle bar to the hole of the handle.

● Sample limitation

⚠ WARNING :

Never use any material capable of producing flammable or explosive vapors. Your centrifuge and rotor are not designed to confine any sample particles dispersed due to leakage. Therefore, when using toxic or radioactive samples or pathogenic or infectious blood samples, make sure to prepare necessary safety measures at your own responsibility.

● The rotor life management of the rotor with optical adapter/disk

⚠ CAUTION :

Each time the rotor with optical adapter/disk is used, be sure to enter the results in the “rotor log book”. Failure to do so will disable management of the life of rotor. No entry is need for RLM rotor.

● RLM adapter and Optical adapter/disk

⚠ WARNING :

- Do not remove the RLM adapter or optical adapter/disk from rotor, or replace it with the adapter/disk for another rotor. The adapter/disk is a critical component that detects the over-speed of rotor : If an adapter/disk that is compatible with the rotor is attached, the rotor could break, resulting in damage to the ultracentrifuge.

⚠ CAUTION :

- Do not bring the RLM adapter near a magnet, or scratch it: Doing so will erase the memory stored in adapter, and make the rotor unusable. To prevent the adapter from being scratched, store the rotor with RLM adapter, using the stand provided with the rotor (rotor stand for protecting adapter).
If the adapter is scratched, immediately stop using rotor and contact an authorized sales or service representative.
- Take great care not to scratch the optical adapter/disk: Any scratch on adapter/disk will make the rotor unusable. To prevent the adapter/disk from being scratched, store the rotor with optical adapter/disk, using the stand provided with the rotor (rotor stand for protecting adapter/disk). If the adapter/disk is scratched, immediately stop using rotor and contact an authorized sales or service representative.
- If foreign matter adheres to the RLM adapter or optical adapter/disk, wipe it off promptly. Otherwise, the RLM adapter or optical adapter/disk may be corroded.

4. Maintenance


● Maintenance of rotor

⚠ WARNING :

Do not allow the rotor temperature to rise over 100°C. Otherwise, the rotor can be brittle.

⚠ CAUTION :

Use a neutral detergent having a pH between 5 and 9. Otherwise, the rotor can be corroded.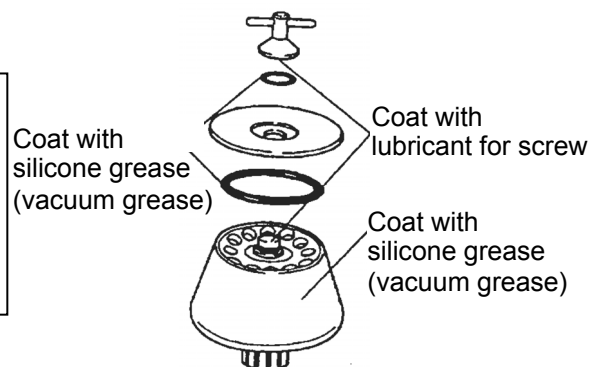

Wash the rotor with tap water or a dilute solution of neutral detergent and rinse it out with distilled water. Turn the rotor upside down with the rotor removed to dry it well. Check the rotor is completely dried and then put the light coat of silicone grease (vacuum grease) on the surface of the rotor, the light coat of lubricant for screw on the thread portion, the light coat of silicone grease (vacuum grease) on the packing (for the cover), Store them in a dry place. If foreign matter adheres to the rotor, soak the rotor in warm water in one or two hour and remove foreign matter with a soft brush etc. Use a neutral detergent having a pH between 5 and 9. If you cannot remove it, contact an authorized sales or service representative.

● Sterilizing rotor

⚠ WARNING :

Never sterilize the rotor by autoclaving or boiling. Otherwise the strength of rotor may be significantly decreased.

⚠ CAUTION :

Do not sterilize the tube adapters made of plastic by autoclaving or boiling. Otherwise the tube adapters may be deformed. Such the tube adapters can not be used.

Sterilize the rotor according to either gas sterilization method (ethylene oxide or formaldehyde) or chemical sterilization method (70% ethanol, 3% hydrogen peroxide, 3% formalin).

Sterilize the tube adapters made of plastic according to either gas sterilization method (ethylene oxide or formaldehyde) or chemical sterilization method (70% ethanol, 3% hydrogen peroxide, 3% formalin).

* Do not dip the rotor and the tube adapters in the formalin (3%) solution more than 2 hours.

● Inspecting rotor

Check the rotor after every run as follows.

- The thread portion is worn. Contact an authorized sales or service representative.
- The tube cavity is corroded. Contact an authorized sales or service representative.
- The packing (for the cover) is deteriorated. . . . Replace it with the new one.
- The adapter is stained. Wipe it out with a soft cloth.

Check the rotor once a month as follows.

- The inside of drive hole (crown hole) is stained. . . . Wipe it out with a soft cloth dampened with water.

⚠ CAUTION :
Clean the inside of the drive hole (crown hole) of the rotor and the surface of the drive shaft (crown) of the centrifuge once a month. If the drive hole or the drive shaft is stained or any foreign matter is adhered, the rotor may be improperly installed and come off during operation.

● Life of rotor

⚠ WARNING :
Do not use rotors that have exceeded their lifetime; to do so might cause the rotor failure and the centrifuge might be damaged.


While using rotor repeatedly, it's strength decreases gradually due to fatigue and creep of material, by which the rotor life is determined.

This rotor must not be used when they reached 5,000 operation times or 10,000 accumulated hours. If you use the RLM rotor at a speed lower than its maximum permissible speed, the life of the RLM rotor will be prolonged automatically by calculating from the strength of the rotor material.

● Replacing the optical disk

(1) Preparation

- Prepare a new optical disk and knife.
Make sure that the number of stripes of the new optical disk is right as follows.
- Wash and then dry the rotor well.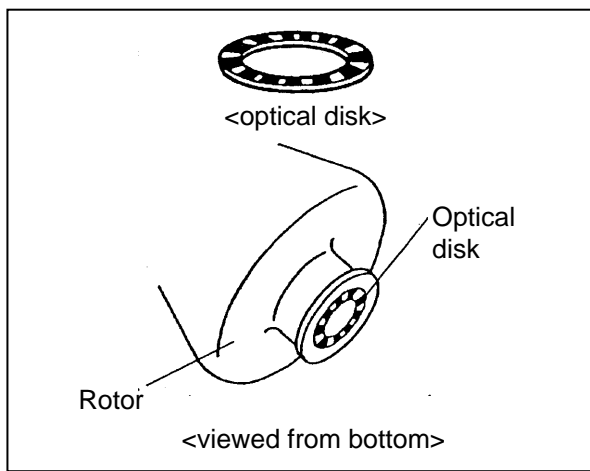

Rotor's maximum speed and optical disk's number of stripes

Rotor's maximum speed (rpm)	Optical disk's number of black bands	Rotor's maximum speed (rpm)	Optical disk's number of black bands	Rotor's maximum speed (rpm)	Optical disk's number of black bands
100,000	9	50,000	18	30,000	29
90,000	10	48,000	19	28,000	32
80,000~83,000	11	45,000	20	27,000	33
70,000	13	42,000	21	25,000	35
65,000	14	40,000	22	23,000	39
60,000	15	38,000	24	21,000	40
55,000~56,000	16	35,000	25	19,000	45
54,000	17	32,000	28	17,000	50

- (2) Hold the rotor upside down.
- (3) Pry the edge of the optical disk with the knife and remove the disk. Be very careful not to damage the rotor.
- (4) Clean the disk hole on the rotor by using alcohol.
- (5) Remove the backing paper from the new optical disk.
Place the optical disk so that it snugly fits into the groove of the disk hole on the rotor. Make sure the disk does not move.

5. Using cesium chloride

The cesium chloride (CsCl) solution is frequently used as a density gradient medium, but if the CsCl solution with a high density is used, it may saturate during the rotation of the rotor depending on the rotation condition which may cause CsCl crystals to be educed. Educated crystals have a high density (approx.4g/mL) and apply an excessive load to the rotor, which is very dangerous.

The crystallization will also greatly affect the density gradient of the CsCl solution and separation status of the sample. Therefore, always use a solution which will not crystallize the allowable speed.

Fig.5-1 shows the relationship between the maximum allowable density and the speed at which cesium chloride does not crystallize in the condition that standard tube is filled with the cesium chloride solution. In this diagram, crystallization will occur above the curves.

Be sure to select the density of the cesium chloride solution and the rotor speed at which cesium chloride does not crystallize. The curve in this diagram varies in temperature etc. Before actual operation, centrifuge the cesium chloride solution with the same density to ensure that crystallization will not occur.

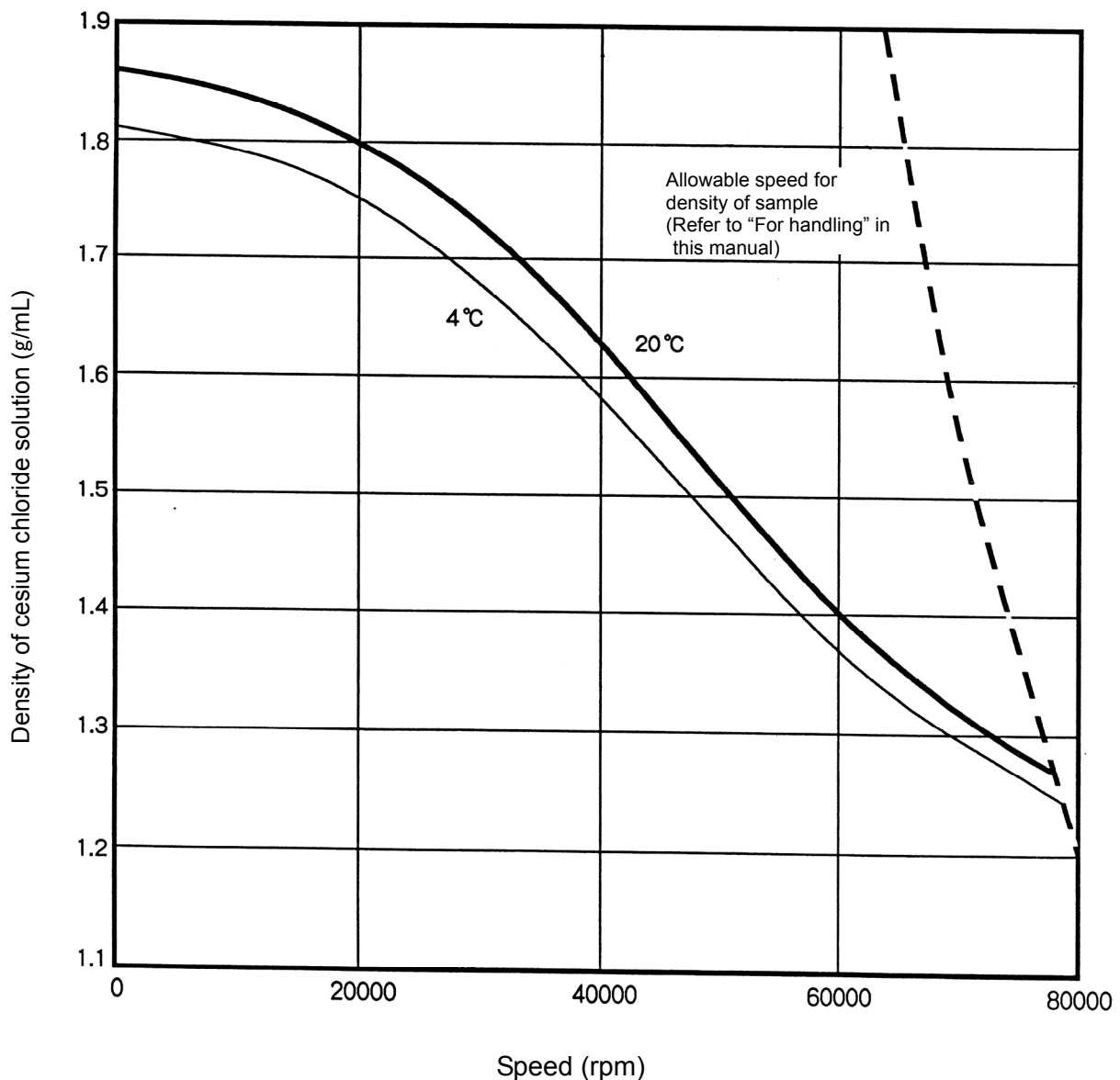
Fig.5-1. Relationship between density of cesium chloride and speed

6. Packing list

Check the parts and accessories of rotor when receiving the rotor.
Contact an authorized sales or service representative when the need arises.

No.	Name	Qty	Remark	Part No. (Qty)	No.	Name	Qty	Part No. (Qty)
①	Rotor	1			⑩			
②	Handle	1*1			⑪			
③	Packing (for the handle) (O-Ring)	1*1		8062026 (1)	⑫	Rotor stand	1	216715 (1)
④	Cover	1*1			⑬	Lubricant for screw	1	84810601 (1)
⑤	Packing (for the cover) (O-Ring)	1*1		S4017039 (1)	⑭	Silicone grease (vacuum grease)	1	483719 (1)
⑥	Handle bar	1		468395 (1)	⑮	Tool box(C)	1	
⑦	Packing (for the handle) (O-Ring)	2	For spare		⑯	Rotor log book *3	1	999494 (1)
⑧	Packing (for the cover) (O-Ring)	2	For spare		⑰	Instruction manual	1	S999240 (1)
⑨	Optical disk *2	2	For spare	346444	⑱			

*1: the quantities of the parts that are mounted to the rotor

*2: The "optical disk" is only provided with the rotor destined for Europe and U.S.A.

*3: The "rotor log book" is not included in RLM rotor.

Tubes are not included in this rotor. Please purchase the applicable tubes.

When using 12PA tubes, 12PA seal tubes, 8PA seal tubes, and 6PA seal tubes, we recommend the purchase of under-mentioned tube kit, what is the set of tubes and necessary tools.

●12PA Tube Kit (A) (Part No.S304986A)
(component)

329606A	12PA Tube :	2 (50 pcs./pack)
463577A	C-TI Cap Ass'y :	12
S401803A	Packing (for the cap): (O-ring)	2 (10 pcs./pack)
333885A	Tube vise (C2)	1
469283	Torque wrench	1
435823A	Tube setter	1

●12PA Seal Tube Kit (A) (Part No.S304991A)
(component)

345320A	12PA Seal Tube :	2 (50 pcs./pack)
486582	C2-Space Cap :	12
S407157	Tube setter :	1

※ When using 12PA seal tubes, tube sealer (STF3,STF2,STF-1) and tube rack (C) (Part No. S201778B) are needed.

●6PA seal tube kit (Part No.S311557A)
(component)


S311555A	6PA seal tube	2 (50 pcs./pack)
486582	C2 space cap	4
S413977	C12-6 float spacer	4
S407157	Tube setter	1
435823A	Tube setter (B)	1

※ When using 6PA seal tubes, Tube sealer (STF3,STF2) and Tube rack (C4) (Part No. S206017A) are needed.

●8PA seal tube kit (Part No.S311562B)
(component)

S408831A	8PA seal tube	2 (50 pcs./pack)
486582	C2 space cap	4
S413976	C12-8 float spacer	4
S407157	Tube setter	1
435823A	Tube setter (B)	1

※ When using 8PA seal tubes, Tube sealer (STF3,STF2) and Tube rack (C2) (Part No. S201778L) are needed.

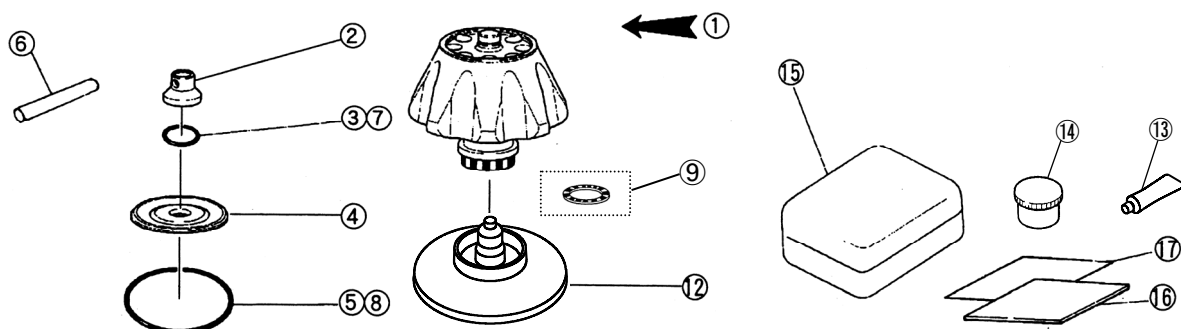7. Decontamination

⚠ WARNING:

- If the centrifuge, rotor or an accessory is contaminated by samples that are toxic or radioactive, or blood samples that are pathogenic or infectious, be sure to decontaminate the item according to good laboratory procedures and methods.
- If there is a possibility that the rotor or an accessory is contaminated by samples that might impair human health (for example, samples that are toxic or radioactive, or blood samples that are pathogenic or infectious), it is your responsibility to sterilize or decontaminate the rotor or the accessory properly before requesting repairs from an authorized sales or service representative. Note that we cannot repair the centrifuge, rotor or the accessory unless sterilization or decontamination is completed.
- It is your responsibility to sterilize and/or decontaminate the rotor or parts properly before returning them to an authorized sales or service representative. In such cases, copy the attached decontamination sheet and fill out the copied sheet, then attach it to the item to be returned. We may ask you about the treatment for the rotor or parts if the decontamination is checked and judged as insufficient by us. It is your responsibility to bear the cost of sterilization or decontamination. Note that we cannot repair or inspect the rotor or the accessory unless sterilization or decontamination is completed.

8. Rotor retirement

After many years of use, there will be inevitably some corrosion or stress corrosion. At some points, the combination of such damage and metal fatigue could make the rotor vulnerable to a failure. Although a rotor may appear to be in a good condition, you should follow the rotor retirement recommendation shown below.

Rotor	Material	Retire After Years
Angle rotor	Titanium alloy	12
Vertical rotor	Aluminum alloy	10
Swing rotor	Titanium alloy Aluminum alloy	10

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Copyright © 2022 Eppendorf Himac Technologies Co., Ltd. All rights reserved.

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