8. Rotor Retirement

(Appendix) K Factors

Keep this rotor instruction manual and the centrifuge manual in the file.

# Angle Rotor for Refrigerated Centrifuge

# R10A5

(Serial No. 1001 or later)

Before using this rotor, please carefully read this instruction manual and the centrifuge instruction manual for its efficient operation and for your safety. Keep this instruction manual for your reference and refer to it as required.

# Contents 1. Specifications — 1 2. Bottles — 3 3. How to Use — 4 4. Maintenance — 6 5. Packing List — 8 6. Spare Parts — 8 7. Decontamination — 9

2022.01 \$99983510 Original instruction manual

# 

• Centrifuge rotors rotating at high speed have considerable potential for damage to personal properties if used improperly.

For safe and proper use of this rotor, carefully read the centrifuge instruction manual and this rotor instruction manual before use and observe the instructions.

⚠ WARNING: and ⚠ CAUTION: notes are used to call your attention in this manual to prevent personal injury or damage to the rotor and the centrifuge.

These notes are defined as follows.

WARNING: indicates a potentially hazardous situation which, if not avoided, could result in personal severe injury or possible death.

CAUTION: indicates a hazardous situation which, if not avoided, could result in personal injury or severe damage to the instrument.

# **⚠** WARNING

- Never use any material capable of producing flammable or explosive vapors.
- Your centrifuge and rotor are not designed to confine any sample particles dispersed due
  to leakage. Therefore, when using toxic or radioactive samples or pathogenic or infectious
  blood samples, make sure to prepare necessary safety measures at your own
  responsibility.
- Never exceed the maximum speed of the rotor(mentioned on the rotor).
   Always reduce rotor speed as instructed in this manual when rotor speed is limited due to sample density or kinds of tubes.
- Check the attached chemical resistance chart, and do not use any sample inapplicable to the rotor.
- If the centrifuge, rotor, or an accessory is contaminated by samples that are toxic or radioactive, or blood samples that are pathogenic or infectious, be sure to decontaminate the item according to good laboratory procedures and methods.
- If there is a possibility that the centrifuge, rotor, or an accessory is contaminated by samples that might impair human health (for example, samples that are toxic or radioactive, or blood samples that are pathogenic or infectious), it is your responsibility to sterilize or decontaminate the centrifuge, rotor, or the accessory properly before requesting repairs from an authorized sales or service representative.
- It is your responsibility to sterilize and/or decontaminate the centrifuge, rotor, or parts properly before returning them to an authorized sales or service representative.

# **CAUTION**

- Do not run the rotor without the rotor cover in position.
   Tighten the cover-locking knob securely.
- Balance the tubes/bottles and the samples including the caps and the adapters within the allowable imbalance of the rotor. Do not exceed the allowable imbalance.
- Clean the inside of the drive hole (crown hole) of the rotor and the surface of the drive shaft (crown) of the centrifuge once a month. If the drive hole or the drive shaft is stained or any foreign matter is adhered, the rotor may be improperly installed and come off during operation.
- Use the rotor tubes and bottles within their actual capacities.
- Do not use tubes/bottles that have exceeded their life expectancy. Failure to do so could result in damage of tubes/bottles and the rotor and the centrifuge. The life expectancy of tubes/bottles depends on factors such as the characteristics of samples, speed of the rotor used, and temperature. Always check for deterioration and damage (cracks, deformation, and so on) on tubes/bottles before using them. Do not use the tubes/bottles if you find such a problem.
- Inspect and maintain the rotor after use. If abnormality is observed, do not use it.
   Contact an authorized sales or service representative.

# 1.Specifications

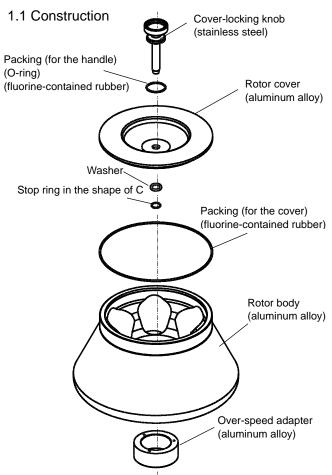
Figure 1-1 Construction of R10A5 rotor

### 1.2 Specifications

Applicable centrifuge\*1

(Some centrifuges are on sale in only Japan.

Visit our web site.)

Max. speed\*2: 10,000 rpm Max. RCF:  $18,900 \times g$ 

Nominal rotor capacity: 500 mL × 6 bottles = 3000 mL

Rotor Size:  $\phi$  348 × 219 mm

Weight: 15.5 kg
Material: Aluminum alloy

Rotor No.: 65

- \*1: For applicable centrifuges, refer to the "Applicable centrifuges (Rotors for high-speed refrigerated centrifuges) (Part No. S998611)".
- \*2: The maximum allowable speed varies with centrifuge models and so on. For further details, refer to the "Applicable centrifuges (Rotors for high-speed refrigerated centrifuges) (Part No. S998611)".

### 1.3 Cross-sectional view of rotor

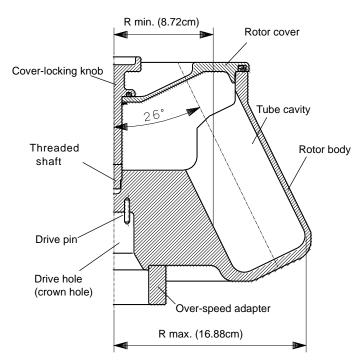
Figure 1-2 Cross-sectional view of R10A5 rotor

### 1.4 Characteristic of rotor

table1-1 Characteristics

Rotor	RCF		
speed (rpm)	R = R min. (8.72cm)	R = R max. (16.88cm)	K factor*
· · ·	` ,		44.00=
2000	390	756	41,827
4000	1,560	3,020	10,457
6000	3,510	6,800	4,647
8000	6,240	12,100	2,614
10000	9,760	18,900	1,673

<sup>\*</sup>For details about K factor, refer to "(Appendix) K Factors"

Calculation formula

 $RCF = 1.118 \times 10^{-5} \times R \times N^2 (\times g)$ 

R: Rotating radius (cm)

N: speed(rpm)

### 1.5 Allowable speed of the rotor

### WARNING:

Do not exceed the maximum speed of the rotor. The rotor speed should be limited depending on sample characteristic, kinds of bottles and centrifuge model.

Do not exceed the allowable speed of the rotor.

The maximum speed marked on the surface of the rotor cover is permitted under the conditions as follows.

- -Average sample density: less than 1.2g/mL
- -500mL bottles are used.

Depending on sample density, kinds of bottle and centrifuge model, the rotor speed may be limited as follows. Do not exceed the allowable speed of the rotor.

### (1) Allowable speed for sample density

### WARNING:

When sample density exceeds 1.2g/mL, calculate the allowable speed according to the following equation.

Allowable speed (rpm) = 
$$10,000$$
(rpm) ×  $\sqrt{\frac{1.2(g/mL)}{\text{Average density of sample}(g/mL)}}$ 

### (2) Allowable speed for kinds of bottles.

The rotor speed should be limited by kinds of bottles.

For details, refer to "2.1 Applicable bottles".

Perform operation under the allowable speed, otherwise bottles may be broken during operation.

### 2.Bottles

### 2.1 Applicable bottles

Table2-1 Applicable bottles

Name	Size (mm)	Actual capacity	Max. speed	Max. RCF
Nalgene <sup>™</sup> Polycarbonate Centrifuge Bottles (Nominal Capacity: 500 mL) Catalog# 3122-0500	φ 69.5 × 169	400 mL	10,000 rpm	18,900×g
Nalgene <sup>™</sup> PPCO Centrifuge Bottles (Nominal Capacity: 500 mL) Catalog# 3120-9500	φ 69.5 × 170.2	400 mL	10,000 rpm	18,900×g

If the bottle is filled to 90% of actual capacity at 4°C to 25°C, the allowable maximum speed and maximum RCF described in Table 2.1 are applicable. Bottles might be deformed and damaged depending on the characteristics of sample, speed of the rotor used, temperature, and so on (see the following.). In that case, perform the rotor at the lower speed.

- The sample volume that is less than 90% of actual capacity of the bottle
- The set temperature that is lower than 4°C
- The set temperature that is higher than 25°C

Refer to the attached instruction manual of the manufacturer for the details.

The allowable speed varies with centrifuge models and so on (see "1.5 Allowable Speed of the Rotor").

- (1) Bottles are not included in this rotor. Please purchase applicable ones selecting from the above list.
- (2) All applicable bottles for this rotor are on the market. 500 PA bottle (Part No. 330437A) can not be used in this rotor. For bottles other than the above, test them by filling with water instead of sample, and run them at the intended speed to ensure that there is no abnormality.
- Maximum speeds listed are guidelines only. Because of variances in user methodologies, no guarantee of performance is expressed or implied.

### 2.2 Cleaning and sterilization of bottles

Clean and sterilize bottles on the market according to the instructions of the manufacturer.

### 2.3 Life expectancy of plastic bottles

Plastic bottles are consumables.

The life of plastic bottle depends upon the sample, speed of rotor, temperature, etc. The life expectancies of the PC bottles specifically depend on the pre-treatment conditions such as cleaning and sterilization.

When the crazing (slight cracks) or the cracks are found on the surface of the bottle, stop using it immediately. Replace it with the new one.

### 3. How to use

### 3.1 Preparation of bottles

Check that the bottles are free from crack or deformation. Do not use remarkably deformed or cracked bottles.

### **WARNING:**

- · These rotors and the centrifuge are not explosion-proof. Never use explosive or flammable samples.
- For safety, there are limits on the use of bio-samples which require bio-isolation, such as pathogenic germs and DNA recombination, as well as RI substance in centrifuges. Perform strict safety controls when separating samples containing these substances.

### CAUTION:

- Do not use any sample inapplicable to the bottles and the caps in use referring to the attached "Chemical resistance chart". Otherwise, the bottles and the caps may be degraded.
- Chemical resistance and the strength of bottles varies with speed, temperature, and so on. Before using sample, fill bottles with water, buffer solution, etc. instead of sample and run them at the intended speed, temperature, and so on to ensure that there is no abnormality.
- Do not exceed the specified sample capacity of the bottles, otherwise sample leakage or deformation of bottles may occur.
- · Balance the sample within the allowable imbalance including the weight of the bottles, caps. Excessive imbalance can cause damage to the centrifuge and the rotor.
- Use the samples that are same in compositions and same in density when placing them symmetrically. Otherwise the precipitation levels may be different by centrifugation and such operation may increase the level of imbalance. And it may result in damage to the centrifuge and the rotor.
- · Balance between symmetrically arranged the bottles/tubes filled with the same-density sample. Variations in sample density cause imbalance operation.
- · Before using bottles on the market, test them by filling with water instead of sample, and run them at the intended speed to ensure that there is no abnormality.
- · Tighten the cap of the bottle securely, otherwise the sample leakage during operation can cause the excessive imbalance, and it may result in damage to the centrifuge and the rotor.

Inject the sample into the bottles and cap the bottles properly. Do not exceed the specified sample capacity of the bottles, otherwise sample leakage or deformation of bottles may occur.

Balance the bottles that are symmetrically arranged. The allowable imbalance of this rotor is within 6g/bottle.

### 3.2 Setting of bottles into rotor

### CAUTION:

Do not load only one bottle nor arrange the bottle asymmetrically with respect to the drive shaft, otherwise the centrifuge or the rotor may be damaged due to excessive imbalance operation.

Check that the tube cavities of the rotor are free from foreign substances. Set the balanced bottles symmetrically with respect to the drive shaft as shown in the figure below. When using bottles of even numbers, set bottles in pairs symmetrically with respect to the drive shaft. When using bottles of odd numbers, set the bottles so as to keep the same space between bottles.

Pre-cool the rotor when the operating temperature of the rotor is lower than the room temperature or pre-heat the rotor when the operating temperature of the rotor is higher than the room temperature before setting the bottles in the rotor.

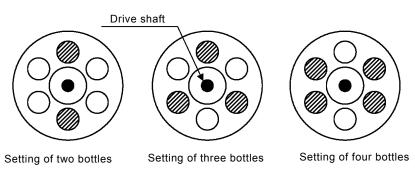
Fig.3-1 Setting of bottles

### 3.3Mounting the rotor cover

### ↑ CAUTION:

Be sure to mount the rotor cover to the rotor body and tighten the cover locking knob securely before operation, otherwise the rotor or the rotor cover may be removed during operation and result in damage to the centrifuge and the rotor.

- (1) Apply silicone grease (vacuum grease) (standard accessory of the centrifuge) to the packing (for the cover) and then fit the packing (for the cover) in the groove of the rotor body or the cover. Replace the packing (for the cover) with new one if damaged or deteriorated.
- (2) Apply lubricant for screw (Part No. 84810601) to the threaded portion of the cover. Put the rotor cover on the rotor body. Turn the cover locking knob to screw in the rotor body securely.

### 3.4 Operation

Mount the rotor onto the drive shaft of the centrifuge gently and securely. Check that the rotor cover is mounted and the cover-locking knob is securely tightened. Perform operation according to the instructions of the centrifuge instruction manual.

### 3.5 Taking out samples

After the operation, gently take out the rotor from the centrifuge and remove the rotor cover by turning the cover locking knob. Take out the bottles from the rotor and then take out the sample in a proper manner.

### 4. Maintenance

### 4.1 Corrosion Resistance of Rotor

### WARNING:

Check the attached chemical resistance chart, and do not use any sample inapplicable to the rotor.

This rotor is made of aluminum alloy. The rotors made of aluminum alloy have high corrosion resistance and they are covered with an anodic oxidation coating. However, use of inapplicable chemicals can corrode these rotors and decrease the strength. Use samples that will not affect the rotor referring to the attached chemical resistance chart.

### 4.2 Maintenance of Rotor

After the operation, properly maintain the rotor to prevent corrosion that can cause the rotor breakage.

• Normal maintenance

### CAUTION:

- Clean the inside of the drive hole (crown hole) of the rotor and the surface of the drive shaft (crown) of the centrifuge once a month. If the drive hole or the drive shaft is stained or any foreign matter is adhered, the rotor may be improperly installed and come off during operation.
- Do not soak the rotor in a dilute solution of neutral detergent or tap water. Failure to do so can cause the rotor to be damaged due to the remaining liquid inside of the rotor.

After operation, wipe the rotor and the rotor cover with a soft cloth which is dampened with a dilute solution of neutral detergent or tap water and wipe the rotor and the rotor cover with a dry cloth. Dry the rotor and the rotor cover separately.

Check that the rotor and the rotor cover are completely dried and then put a light coat of silicone grease (vacuum grease)(standard accessory of the centrifuge) on the rotor and the rotor cover. Store the rotor and the rotor cover in a dry place.

Maintenance after use of a corrosive sample

Wipe the rotor and the rotor cover with a soft cloth which is dampened with tap water soon after operation. Then perform the normal maintenance.

• Maintenance when foreign substances are adhered to the rotor

### CAUTION:

Use a neutral detergent having a pH between 5 and 9, otherwise the rotor can be discolored or corroded.

Wipe the inside of the tube cavities and the crown hole at the rotor bottom with a soft cloth which is dampened with tap water to remove foreign substances.

Be sure to remove foreign substances from the drive hole of the rotor bottom, otherwise the rotor and the drive shaft may be damaged.

### 4.3 Sterilizing Rotor

Sterilize this rotor and the rotor cover in any following methods; autoclaving (121°C, 20minutes), gas sterilization method (ethylene oxide or formaldehyde), chemical sterilization method (70% ethanol, 3% hydrogen peroxide, 3% formalin), or ultraviolet rays (200-300nm) sterilization.

### 

- Never sterilize the rotor whose serial number is 1000 or earlier by autoclaving.
- · Be sure to remove the rotor cover from the rotor before autoclaving, otherwise the rotor and the rotor cover may be deformed.
- Never sterilize the rotor and the rotor cover by boiling, otherwise they will be brittle.

### CAUTION:

- · After autoclaving, wait until the temperature in the autoclaving chamber is the same as the room temperature and then take out the rotor and the rotor cover.
- Take out the rotor and the rotor cover carefully from the autoclaving chamber as they are wet and slippery.
- · Inspect the packing (for the cover) after sterilization. If you observe that the packing (for the cover) is damaged and is deteriorated, replace it with the new one.
- Do not soak the rotor and the rotor cover in chemicals. Failure to do so can cause the rotor to be damaged due to the remaining liquid inside of the rotor.

### 4.4 Inspecting Rotor

Periodically check that the rotor is free from corrosion every 100 hours use. Especially be careful of the tube cavities of the rotor and the crown hole at the rotor bottom because the rotor can be brittle if these portions are corroded. If the rotor surface is discolored, dented or cracked, the rotor is Do not use such a corroded rotor and contact an authorized sales or service corroded. representative immediately for further inspection.

# 5. Packing list


Check the parts and accessories of the rotor when the rotor is delivered. Contact an authorized sales or service representative if there are a missing item and inferior goods.

No.	Name	Qty	Part No. of spare parts (Qty)
1	Rotor	1	_
2	/ Packing (for the handle)	1*	84521017 (3)
3	│ Packing (for the cover)	1*	S411565A (2)
4	Rotor set handle	1	455650B
5	Rotor instruction manual	1	S999835

<sup>\*:</sup> The quantities of the parts that are mounted to the rotor.

《Note》

Bottles are not included in this rotor. Please purchase applicable ones on the market.

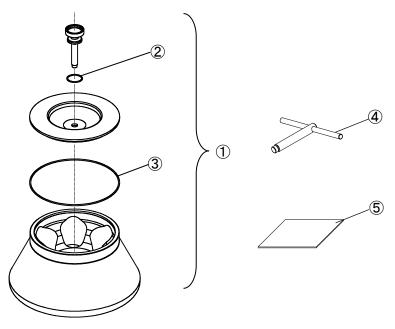# 6. Spare parts

PART No.	Name	Qty
84521017	Packing (for the handle)	3
S411565A	Packing (for the cover)	2
84810601	Lubricant for screw	1
483719	Silicone grease (vacuum grease)	1
455650B	Rotor set handle	1
S999835	Rotor Instruction manual	1

### 7. Decontamination

# MARNING:

- ●If the centrifuge, rotor or an accessory is contaminated by samples that are toxic or radioactive, or blood samples that are pathogenic or infectious, be sure to decontaminate the item according to good laboratory procedures and methods.
- If there is a possibility that the rotor or an accessory is contaminated by samples that might impair human health (for example, samples that are toxic or radioactive, or blood samples that are pathogenic or infectious), it is your responsibility to sterilize or decontaminate the rotor or the accessory properly before requesting repairs from an authorized sales or service representative. Note that we cannot repair the centrifuge, rotor or the accessory unless sterilization or decontamination is completed.
- It is your responsibility to sterilize and/or decontaminate the rotor or parts properly before returning them to an authorized sales or service representative. In such cases, copy the attached decontamination sheet and fill out the copied sheet, then attach it to the item to be returned.
  We may ask you about the treatment for the rotor or parts if the decontamination is checked and judged as insufficient by us. It is your responsibility to bear the cost of sterilization or decontamination. Note that we cannot repair or inspect the rotor or the accessory unless sterilization or decontamination is completed.

## 8. Rotor retirement

After many years of use, there will be inevitably some corrosion or stress corrosion. At some points, the combination of such damage and metal fatigue could make the rotor vulnerable to a failure. Although a rotor may appear to be in a good condition, you should follow the rotor retirement recommendation shown below.

Rotor	Material	Retire After Years
Defrigerated centrifuge retere	Aluminum alloy	15
Refrigerated centrifuge rotors	Stainless steel	15

# (Appendix) K Factors

### Calculation of separation time

The K factor can conveniently be used which makes it possible to estimate the settling time directly from Svedberg unit(S).

Namely, the time taken for particles of 1S to settle from Rmin. to Rmax. is K factor.

The K factor can be obtained from the formula show below.

$$K = \frac{\ln(R \max .) - \ln(R \min .)}{\omega^2} \cdot \frac{10^{13}}{3600}$$

ω: Angular velocity ω = 2π/60 × N

N : Speed (rpm)

Rmin. : Distance from axis of rotation to top of solution (cm) Rmax. : Distance from axis of rotation to bottom of tube (cm)

Using the K factor, the settling time t(in hours) is expressed as follows:

$$t = \frac{K}{S}$$

For example, the K factor at the maximum speed of 10,000rpm of Model R10A5 rotor is 1673. The time taken for particles of 1000 S to be settled from R min. to R max. using this rotor is as follows:

$$t = \frac{1673}{1000} = 1.7 \text{ hours}$$

The time taken for settling of particles of 1000 S by the same rotor at a speed of 8,000 rpm is:

$$K = \left(\frac{10,000}{8,000}\right)^2 \times 1673 = 2614$$
  $t = \frac{2614}{1000} = 2.6 \text{ hours}$ 

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このマニュアルの内容はすべて著作権により保護されています。このマニュアルの内容の一部または全部を、無断で転載すること は禁じられています。  $\label{lem:copyright} \verb§@ 2022 Eppendorf Himac Technologies Co., Ltd. All rights reserved.$ この取扱説明書に記載されている会社名および製品名は、一般に各社の商標または登録商標です。