Keep this rotor instruction manual and the centrifuge manual in the file.

Angle Rotor for Refrigerated Centrifuge

R13A

Before using this rotor, please carefully read this instruction manual and the centrifuge instruction manual for its efficient operation and for your safety.

Keep this instruction manual for your reference and refer to it as required.

Contents

1. Specifications	_____	1
2. Bottles/Tubes	_____	4
3. How to Use	_____	8
4. Maintenance	_____	10
5. Packing List	_____	12
6. Spare Parts	_____	12
7. Decontamination	_____	13
8. Rotor Retirement	_____	13
(Appendix) K Factors	_____	14

Safety Reminder

- Centrifuge rotors rotating at high speed have considerable potential for damage to personal properties if used improperly.
For safe and proper use of this rotor, carefully read the centrifuge instruction manual and this rotor instruction manual before use and observe the instructions.

 **WARNING :** and  **CAUTION :** notes are used to call your attention in this manual to prevent personal injury or damage to the rotor and the centrifuge.

These notes are defined as follows.

 **WARNING:** indicates a potentially hazardous situation which, if not avoided, could result in personal severe injury or possible death.

 **CAUTION:** indicates a hazardous situation which, if not avoided, could result in personal injury or severe damage to the instrument.

WARNING

- Never use any material capable of producing flammable or explosive vapors.
- Your centrifuge and rotor are not designed to confine any sample particles dispersed due to leakage. Therefore, when using toxic or radioactive samples or pathogenic or infectious blood samples, make sure to prepare necessary safety measures at your own responsibility.
- Never exceed the maximum speed of the rotor (mentioned on the rotor).
Always reduce rotor speed as instructed in this manual when rotor speed is limited due to sample density or kinds of tubes.
- Check the attached chemical resistance chart, and do not use any sample inapplicable to the rotor.
- If the centrifuge, rotor, or an accessory is contaminated by samples that are toxic or radioactive, or blood samples that are pathogenic or infectious, be sure to decontaminate the item according to good laboratory procedures and methods.
- If there is a possibility that the centrifuge, rotor, or an accessory is contaminated by samples that might impair human health (for example, samples that are toxic or radioactive, or blood samples that are pathogenic or infectious), it is your responsibility to sterilize or decontaminate the centrifuge, rotor, or the accessory properly before requesting repairs from an authorized sales or service representative.
- It is your responsibility to sterilize and/or decontaminate the centrifuge, rotor, or parts properly before returning them to an authorized sales or service representative.

CAUTION

- Do not run the rotor without the rotor cover in position.
Tighten the cover-locking knob securely.
- Balance the tubes/bottles and the samples including the caps and the adapters within the allowable imbalance of the rotor. Do not exceed the allowable imbalance.
- Clean the inside of the drive hole (crown hole) of the rotor and the surface of the drive shaft (crown) of the centrifuge once a month. If the drive hole or the drive shaft is stained or any foreign matter is adhered, the rotor may be improperly installed and come off during operation.
- Use the rotor tubes and bottles within their actual capacities.
- Do not use tubes/bottles that have exceeded their life expectancy. Failure to do so could result in damage of tubes/bottles and the rotor and the centrifuge. The life expectancy of tubes/bottles depends on factors such as the characteristics of samples, speed of the rotor used, and temperature.
Always check for deterioration and damage (cracks, deformation, and so on) on tubes/bottles before using them. Do not use the tubes/ bottles if you find such a problem.
- Inspect and maintain the rotor after use. If abnormality is observed, do not use it.
Contact an authorized sales or service representative.

1.Specifications

1.1 Construction

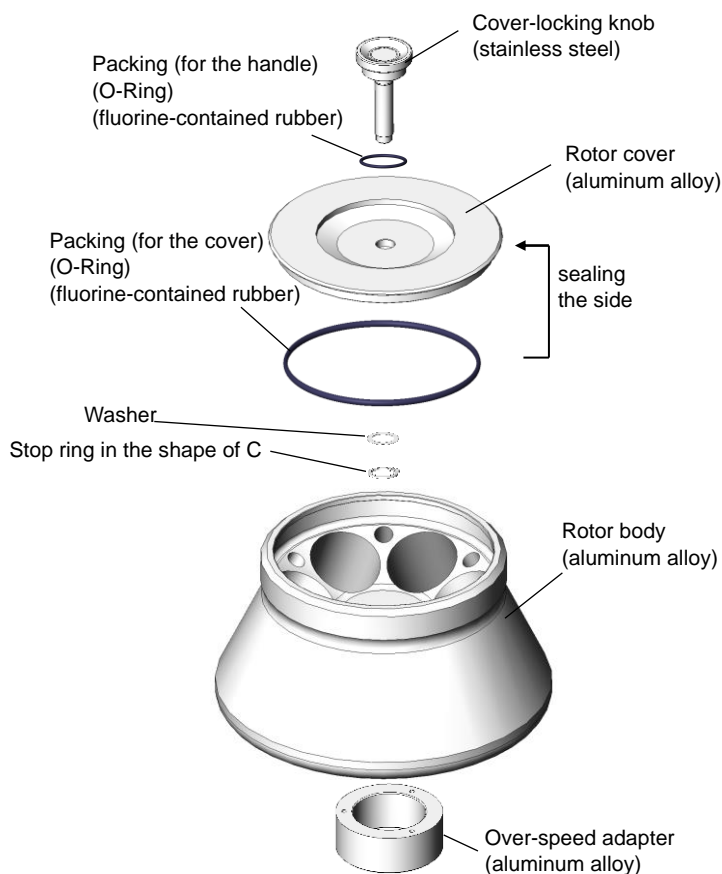
Figure1-1 Construction of R13A rotor

1.2 Specifications

Applicable centrifuge*1

(Some centrifuges are on sale in only Japan.
Visit to our web site.)

Max. speed*2: 13,000 rpm

Max. RCF: 26,400 × g (250PP Bottles)
23,400 × g (11PA Thick-Walled Tubes)

Nominal rotor capacity:

250 mL × 6 bottles = 1500 mL

Rotor Size: φ 290 × 193 mm

Weight: 11.0 kg

Material: Aluminum alloy

Rotor No.: 77

*1: For applicable centrifuges, refer to the “Applicable centrifuges (Rotors for high-speed refrigerated centrifuges (Part No. S998611))”.

*2: The maximum allowable speed varies with centrifuge models and so on. For further details, refer to the “Applicable centrifuges (Rotors for high-speed refrigerated centrifuges (Part No. S998611))”.

1.3 Cross-sectional view of rotor

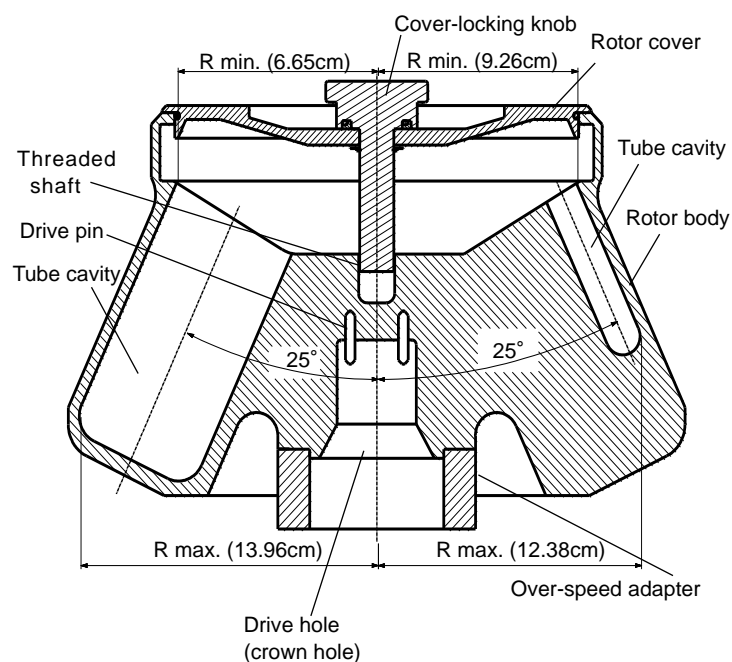
Figure1-2 Cross-sectional view of R13A rotor

1.4 Characteristic of rotor

Table1-1 Characteristics

Rotor speed (rpm)	250PP Bottle			11PA Tube		
	RCF (×g)		K factor ※	RCF (×g)		K factor ※
	R = Rmin. (6.65cm)	R = Rmax. (13.96cm)		R = Rmin. (9.26cm)	R = Rmax. (12.38cm)	
1,000	74	156	187,844	104	138	73,554
2,000	297	624	46,961	414	554	18,388
3,000	669	1,400	20,872	932	1,250	8,173
4,000	1,190	2,500	11,740	1,660	2,210	4,597
5,000	1,860	3,900	7,514	2,590	3,460	2,942
6,000	2,680	5,620	5,218	3,730	4,980	2,043
7,000	3,640	7,650	3,834	5,070	6,780	1,501
8,000	4,760	10,000	2,935	6,630	8,860	1,149
9,000	6,020	12,600	2,319	8,390	11,200	908
10,000	7,430	15,600	1,878	10,400	13,800	736
11,000	9,000	18,900	1,552	12,500	16,700	608
12,000	10,700	22,500	1,304	14,900	19,900	511
13,000	12,600	26,400	1,111	17,500	23,400	435

※For details about K factor, refer to “(Appendix) K Factors”

Calculation formula

$$RCF = 1.118 \times 10^{-5} \times R \times N^2 (\times g)$$

R: Rotating radius (cm)
N: speed(rpm)

1.5 Allowable speed of the rotor

WARNING :

Do not exceed the maximum speed of the rotor. The rotor speed should be limited depending on sample characteristic, kinds of bottles and centrifuge model.

Do not exceed the allowable speed of the rotor.

The maximum speed marked on the surface of the rotor cover is permitted under the conditions as follows.

-Average sample density : less than 1.2g/mL

-250mL bottles are used.

Depending on sample density, kinds of bottle and centrifuge model, the rotor speed may be limited as follows. Do not exceed the allowable speed of the rotor.

(1) Allowable speed for sample density

WARNING :

When sample density exceeds 1.2g/mL, calculate the allowable speed according to the following equation.

$$\text{Allowable speed (rpm)} = 13,000 \text{ (rpm)} \times \sqrt{\frac{1.2(\text{g/mL})}{\text{Average density of sample (g/mL)}}}$$

(2) Allowable speed for kinds of bottles/tubes

The rotor speed should be limited by kinds of bottles/tubes etc.

For details, refer to "2.2 Applicable bottles/tubes".

Perform operation under the allowable speed, otherwise bottles may be broken during operation.

1.6 Rotor registration data for the CR-G II series centrifuges




If this rotor is not registered in the CR-G II series centrifuges, register this rotor by entering the following data according to "3-5-5 Entry of New Rotor" on the instruction manual of the CR-G II series centrifuges.

Item/No.	DATA	Item/No.	DATA	Item/No.	DATA	Item/No.	DATA
ROTOR No.	77	6	2153	12	51065	18	0
1	4000	7	4106	13	32032	19	1154
2	50204	8	42	14	32032	20	0
3	96000	9	16000	15	50200	21	7006
4	5116	10	8001	16	4175	22	14255
5	5096	11	82049	17	0	23	42029

2. Bottles/Tubes

2.1 Kinds of Bottles and Tubes

Following table shows general types of bottles and tubes.

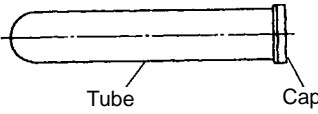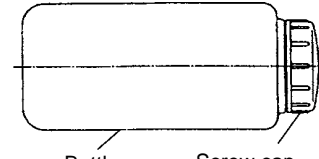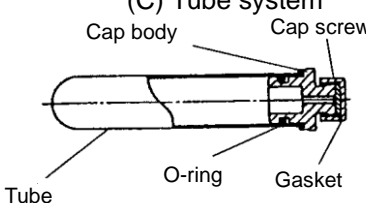Simplified types		Seal-capped type
(A) Tube system 	(B) Bottle system 	(C) Tube system 

Bottles and tubes are made of the following materials.

PC: Polycarbonate

Strong, transparent and can be autoclaved at 121°C or less. Use the tubes/bottles at a temperature between 4 °C and 25 °C. Weak in organic solution, alkali solution, and alkali detergent.

PP: Polypropylene

Strong but susceptible to low temperature (brittle point:0°C). It can be autoclaved at 121°C or less. Use the tubes/bottles at a temperature between 4 °C and 25 °C.

PA: Polypropylene copolymer

Strong and can be autoclaved at 115°C or less. Use the tubes/bottles at a temperature between 4 °C and 25 °C.

SST: Stainless steel

Excellent in chemical resistance and heat resistance. It can be autoclaved at 121°C or less.

Glass: It is transparent and highly heat-resistant.

Refer to the attached "Chemical resistance chart (Part No. S999313)" for chemical resistance of each material.

2.2 Applicable bottles/tubes

Table2-1 Applicable bottles/tubes

Tubes/Bottles				Tube Adapter				Cap		Max. Speed (rpm)	Max. RCF (xg)
Part No. (Qty)	Name	Size (φ X L cm)	Actual capacity (mL/tube)	Part No. (Qty)	Name	Size (φ X L cm)	Material *2	Part No. (Qty)	Name		
S306749A (10 pcs.)	250PP Bottle Ass'y	6.2X12.4	250	---	---	---	---	---	---	13,000	26,400
S306750A (10 pcs.)	250PC Bottle Ass'y	6.2X12.4	250	---	---	---	---	---	---	13,000	26,400
328343A (2 pcs.)	300SST Tube	6.2X12.8	130	---	---	---	---	---	---	13,000	26,400
464977A (50 pcs.)	10 Glass Tube	1.2X10.5	7.1	336717A (4 pcs.)	10G7A300 Adapter	6.1X11.7	POM	---	---	4,000	2,320
473866A (10 pcs.)	50 Glass Tube	3.5X10.0	30	336720A (4 pcs.)	50GA300 Adapter	6.1X11.7	NBR	---	---	4,000	2,180
S311187A (100 pcs.)	50 TC Tube Ass'y	3.0X11.5	43	S408203A (2 pcs.)	50TA300 Adapter	6.1X10.5	POM	---	---	11,000	16,200
Marketed	50mL PP CultureTube *1 (conical bottom)	---	43	S408203A (2 pcs.)	50TA300 Adapter	6.1X10.5	POM	---	---	11,000	16,200
328353A (10 pcs.)	50PA Tube	2.9X10.4	28	S407879A (2 pcs.)	50A300 Adapter (B)	6.1X11.5	PP	328479A (20 pcs.)	M-PP Lid	10,000	13,600
328354A (10 pcs.)	50PC Tube	2.9X10.3	28	S407879A (2 pcs.)	50A300 Adapter (B)	6.1X11.5	PP	328479A (20 pcs.)	M-PP Lid	10,000	13,600
S306352A (10 pcs.)	50PP Bottle Ass'y	2.9X10.3	30	S407879A (2 pcs.)	50A300 Adapter (B)	6.1X11.5	PP	---	---	10,000	13,600
328352A (10 pcs.)	50PC Bottle Ass'y	2.9X10.3	30	S407879A (2 pcs.)	50A300 Adapter (B)	6.1X11.5	PP	---	---	10,000	13,600
328342A (2 pcs.)	50SST Tube	2.9X10.2	34	S407879A (2 pcs.)	50A300 Adapter (B)	6.1X11.5	PP	---	---	10,000	13,600
332094A (10 pcs.)	80PA Bottle Ass'y	3.8X10.4	54	336722A (4 pcs.)	80A300 Adapter	6.1X11.7	NBR	---	---	10,000	13,600
S308432A (10 pcs.)	70PC Bottle Ass'y	3.8X10.4	49	336722A (4 pcs.)	80A300 Adapter	6.1X11.7	NBR	---	---	10,000	13,600
326012A (2 pcs.)	94SST Tube	3.8X10.1	40	336722A (4 pcs.)	80A300 Adapter	6.1X11.7	NBR	---	---	10,000	13,600
333959A (10 pcs.)	11PA Thick-Walled Tube	1.6X8.2	8	---	---	---	---	328478A (20 pcs.)	C2-PP Lid	13,000	23,400

*1: When using tubes on the market, perform operation under the allowable speed or the allowable RCF specified by the manufacturer. Otherwise the tubes may be broken during operation.

*2: POM: Polyacetal NBR: Nitrile rubber PP: Polypropylene

The maximum allowable speed varies with centrifuge models and so on. For further details, refer to the "Applicable centrifuges (Rotors for high-speed refrigerated centrifuges) (Part No. S998611)".

The rotor speed is limited by the sample density and so on (see Section1.5).

 **CAUTION:**
When using bottles/tubes outside 4-25°C, bottles/tubes may be broken or deformed during operation. Test them before actual operation and perform actual operation after checking that there is no inconvenience such as the above.

- Maximum speeds listed are guidelines only. Because of variances in user methodologies, no guarantee of performance is expressed or implied.

2.3 Cleaning and sterilization of bottles/tubes

Properly wash and sterilize the bottles/tubes according to Table 2-2.

Comply with instructions of each maker for washing and sterilization of commercially available tubes.

- Autoclaving PA, PP and PC bottles/tubes

Wash the bottles/tubes well. Put them on a heat-resistant tube stand facing their mouths upward. Remove the screw caps and the lids to prevent deformation or breakage during autoclaving. After autoclaving, wait until the temperature in the autoclaving chamber reduces to the room temperature, then take out the tubes/bottles.

CAUTION:

Do not place the screw cap or the lid on the mouth of the bottle. Otherwise the bottle may be deformed or damaged during autoclaving.

- Washing

Wash the bottles/tubes with tap water or a dilute solution of neutral detergent. If the bottles/tubes stain heavily, soak them in a dilute solution of neutral detergent or wash them with a soft brush being careful not to scratch them. Then rinse the bottles/tubes in tap water and finally rinse them with distilled water. Dry the bottles/tubes in the air. Use a detergent of pH 5 to 9 because the PC tubes are susceptible to alkaline solutions. Avoid soaking the tubes/bottles in a dilute solution of neutral detergent for a long time. Otherwise, the tubes/bottles may be degraded in strength. For the caps of the tubes/bottles, remove them from the tubes/bottles. After disassembling the caps, wash them, then dry and store the caps in the same manner as the tubes/bottles.

Table 2-2 Washing and sterilization of bottles/tubes

○ : Applicable x : Inapplicable

Conditions		Material	PE	PA	PC	PP	SST
Washing	Washing solution	pH 5 or lower	x	x	x	x	x
		pH 5 to 9	○	○	○	○	○
		pH 9 or higher	○	○	x	○	○
		Warm water (50°C or lower)	○	○	○	○	○
Ultrasonic washing	Neutral detergent (pH 7)	○	○	○	○	○	
Sterilization	High-pressure steam sterilization (Autoclaving)	115°C(0.7kg/cm ²), 30 minutes	x	○	○	○	○
		121°C (1.0kg/cm ²), 20 minutes	x	x	○	○	○
		126°C (1.4kg/cm ²), 15 minutes	x	x	x	x	x
	Boiling	For 15 to 30 minutes	x	○	○	○	○
	Ultraviolet ray	200-300nm	x	x	x	x	○
	Gas	Ethylene oxide	○	○	x	○	○
		Formaldehyde	○	○	○	○	x
Chemical	Ethanol (70%)	○	○	x	○	○	
	Hydrogen peroxide (3%)	○	○	○	○	○	
	Formalin (3%)	○	○	○	○	x	

PE: Polyethylene PA: Polypropylene copolymer PC: Polycarbonate
 PP: Polypropylene SST: Stainless steel

2.4 Life expectancy of plastic bottle and tube

The life of plastic bottle and tube depends upon the sample, speed of rotor, temperature, etc. The standard life of plastic bottle and tube when used with aqueous samples (of pH5 to pH9) is specified as follows:

- (1) Tubes (PA, PP, PC and PE) 5 times
- (2) Thick-walled tube and bottle (PA, PP, PC and PE) 50 times
- (3) 50TC tubes and Commercial tubes 1 time

However, the standard life of the PC thick-walled tube and PC bottle is specified as follows depending upon the pre-treating conditions such as washing, sterilization, etc.:

Table 2-3 Standard Life of tube and bottle

Sample \ Washing and sterilization	Sterilization in gas or washing in hot water	Autoclaving at 121°C x 20min
Neutral (pH7)	50 times	10 times
Weak alkali (pH7 to pH9)	30 times	5 times

Do not use crazed (cracked) tubes or bottles.

The standard life of plastic tubes and bottles as the above are an approximate guide when they are used repeatedly for 1 hour at the maximum speed of rotor. We do not warrant the life expectancies of bottles/tubes.

3. How to Use

3.1 Preparation of bottles

Check that the bottles are free from crack or deformation. Do not use remarkably deformed or cracked bottles.

⚠ WARNING :

- These rotors and the centrifuge are not explosion-proof. Never use explosive or flammable samples.
- For safety, there are limits on the use of bio-samples which require bio-isolation, such as pathogenic germs and DNA recombination, as well as RI substance in centrifuges. Perform strict safety controls when separating samples containing these substances.

⚠ CAUTION:

- Do not use any sample inapplicable to the bottles, the caps, and the tubes in use referring to the attached "Chemical resistance chart. Otherwise, the bottles/tubes and the caps may be degraded.
- Chemical resistance and the strength of bottles/tubes varies with speed, temperature, and so on. Before using sample, fill bottles/tubes with water, buffer solution, etc. instead of sample and run them at the intended speed, temperature, and so on to ensure that there is no abnormality.
- Do not exceed the specified sample capacity of the bottles/tubes, otherwise sample leakage or deformation of bottles/tubes may occur.
- Balance the sample within the allowable imbalance including the weight of the bottles/tubes, caps. Excessive imbalance can cause damage to the centrifuge and the rotor.
- Use the samples that are same in compositions and same in density when placing them symmetrically. Otherwise the precipitation levels may be different by centrifugation and such operation may increase the level of imbalance. And it may result in damage to the centrifuge and the rotor.
- Balance between symmetrically arranged the bottles/tubes filled with the same-density sample. Variations in sample density cause imbalance operation.
- Tighten the cap of the tube/bottle securely, otherwise the sample leakage during operation can cause the excessive imbalance, and it may result in damage to the centrifuge and the rotor.

Inject the sample into the bottles/tubes and cap the bottles/tubes properly.

If the bottle/tube lid without the O-ring is used, do not invert or shake the bottle/tube with the lid, otherwise the sample may leak.

Do not exceed the specified sample capacity of the bottles/tubes, otherwise sample leakage or deformation of bottles/tubes may occur.

Balance the bottles/tubes that are symmetrically arranged. The allowable imbalance of this rotor is within 4g/bottle.


3.2 Setting of bottles into rotor

⚠ CAUTION:

Do not load only one bottle nor arrange the bottle asymmetrically with respect to the drive shaft, otherwise the centrifuge or the rotor may be damaged due to excessive imbalance operation.

Check that the tube cavities of the rotor are free from foreign substances. Set the balanced bottles symmetrically with respect to the drive shaft as shown in the figure below. When using bottles of even numbers, set bottles in pairs symmetrically with respect to the drive shaft. When using bottles of odd numbers, set the bottles so as to keep the same space between bottles.

Pre-cool the rotor when the operating temperature of the rotor is lower than the room temperature or pre-heat the rotor when the operating temperature of the rotor is higher than the room temperature before setting the bottles in the rotor.

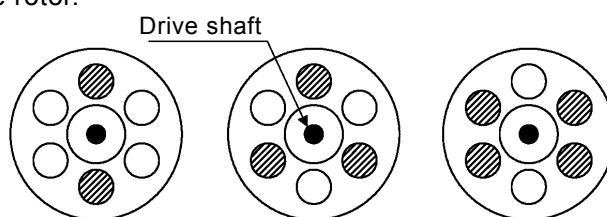Setting of two bottles Setting of three bottles Setting of four bottles

Fig.3-1 Setting of bottles

3.3 Mounting the rotor cover

CAUTION:

Be sure to mount the rotor cover to the rotor body and tighten the cover locking knob securely before operation, otherwise the rotor or the rotor cover may be removed during operation and result in damage to the centrifuge and the rotor.

- (1) Apply silicone grease (vacuum grease) (standard accessory of the centrifuge) to the packing (for the cover) and then fit the packing (for the cover) in the groove of the rotor body or the cover. Replace the packing (for the cover) with new one if damaged or deteriorated.
- (2) Apply lubricant for screw (Part No. 84810601) to the threaded portion of the cover. Put the rotor cover on the rotor body. Turn the cover locking knob to screw in the rotor body securely.

3.4 Operation

Mount the rotor onto the drive shaft of the centrifuge gently and securely. Check that the rotor cover is mounted and the cover-locking knob is securely tightened. Perform operation according to the instructions of the centrifuge instruction manual.

3.5 Taking out samples

After the operation, gently take out the rotor from the centrifuge and remove the rotor cover by turning the cover locking knob. Take out the bottles/tubes from the rotor and then take out the sample in a proper manner.

4. Maintenance

4.1 Corrosion resistance of rotor

WARNING:

Check the attached chemical resistance chart, and do not use any sample inapplicable to the rotor.

This rotor is made of aluminum alloy. The rotors made of aluminum alloy have high corrosion resistance and they are covered with an anodic oxidation coating. However, use of inapplicable chemicals can corrode these rotors and decrease the strength. Use samples that will not affect the rotor referring to the attached chemical resistance chart.

4.2 Maintenance of rotor

After the operation, properly maintain the rotor to prevent corrosion that can cause the rotor breakage.

- Normal maintenance

CAUTION:

Clean the inside of the drive hole (crown hole) of the rotor and the surface of the drive shaft (crown) of the centrifuge once a month. If the drive hole or the drive shaft is stained or any foreign matter is adhered, the rotor may be improperly installed and come off during operation.

Wash the rotor with tap water or a dilute solution of neutral detergent and rinse it out with distilled water. Wipe the rotor with a soft cloth. Turn the rotor upside down with the rotor cover removed to dry it well. Check that the rotor is completely dried and then put a light coat of silicone grease (vacuum grease) (standard accessory of the centrifuge) on the rotor. Store the rotor in a dry place.

- Maintenance after use of a corrosive sample

Wash the rotor with tap water immediately after the operation. Then perform the normal maintenance.

- Maintenance when foreign substances are adhered to the rotor

CAUTION:

Use a neutral detergent having a pH between 5 and 9, otherwise the rotor can be discolored or corroded.

Soak the rotor in warm water for one or two hours and wash the inside of the tube cavities and the crown hole at the rotor bottom with a soft brush to remove foreign substances. Do not forget to remove foreign substances adhered in the crown hole at the rotor bottom, otherwise the rotor and the drive shaft may be damaged.

4.3 Sterilizing rotor

- Sterilize this rotor according to either gas sterilization method (ethylene oxide or formaldehyde) or chemical sterilization method (70% ethanol, 3% hydrogen peroxide, or 3% formalin) or autoclaving (121°C, 20 minutes).
- Sterilize the plastic or rubber adapters according to either gas sterilization method (ethylene oxide or formaldehyde) or chemical sterilization method (70% ethanol, 3% hydrogen peroxide, or 3% formalin).

 **WARNING:** Be sure to remove the rotor cover from the rotor before autoclaving, otherwise the rotor or the rotor cover may be deformed.

 **WARNING:** Never sterilize the rotor by boiling, otherwise the rotor will be brittle.

 **CAUTION :** Never sterilize the adapters by autoclaving or boiling, otherwise the adapters may be deformed. Such adapters cannot be used.

 **CAUTION :** After autoclaving, wait until the temperature in the autoclaving chamber reduces to the room temperature, then take out the rotor and the rotor cover.

 **CAUTION :** Take out the rotor and the rotor cover carefully from the autoclaving chamber as they are wet and slippery.

 **CAUTION :** Check the packing (for the cover) after sterilizing. The packing (for the cover) may be damaged during operation due to deterioration. Replace the new one if a damage or deterioration is observed.

 **CAUTION :** Do not dip the rotor and the adapters in the formalin (3%) solution more than 2 hours.

4.5 Inspecting rotor


Periodically check that the rotor is free from corrosion every 100 hours use. Especially be careful of the tube cavities of the rotor and the crown hole at the rotor bottom because the rotor can be brittle if these portions are corroded. If the rotor surface is discolored, dented or cracked, the rotor is corroded. Do not use such a corroded rotor and contact an authorized sales or service representative immediately for further inspection.

5. Packing List

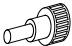
Check the parts and accessories of the rotor when the rotor is delivered. Contact an authorized sales or service representative if there are a missing item and inferior goods.

No.	Name	Qty	Part No. of spare parts (Qty)
①	Rotor	1	—
②	Packing (for the handle) (O-Ring)	1*	84521017 (3)
③	Packing (for the cover) (O-Ring)	1*	S411569A (3)
④	250PP Bottle Ass'y	6	S306749A (10)
⑤	Bottle setter	1	84550301 (1)
⑥	Rotor instruction manual	1	S999842

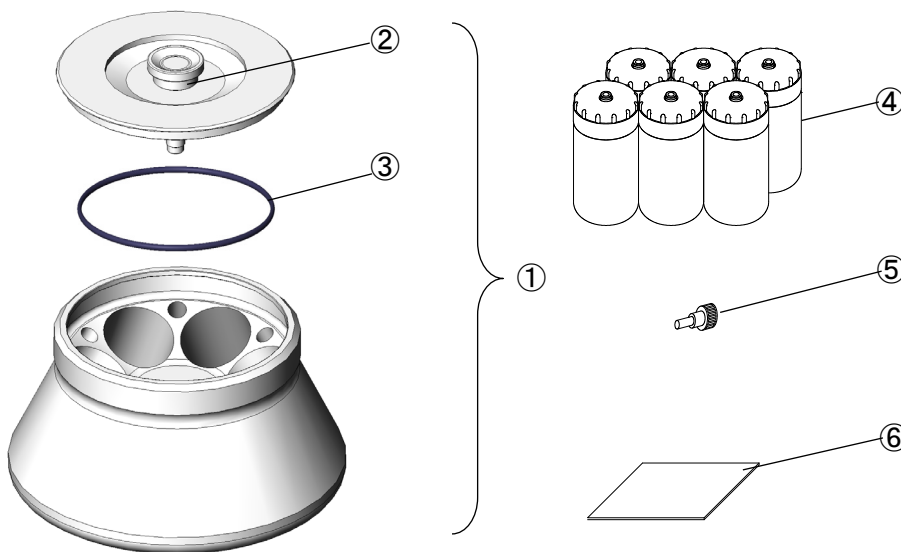*: The quantities of the parts that are mounted to the rotor.

6. Spare Parts

	Part No.	Name	Qty
	84521017	Packing (for the handle) (O-ring)	3
	S411569A	Packing (for the cover) (O-ring)	3
	84810601	Lubricant for screw	1
	483719	Silicone grease (vacuum grease)	1
	84550301	Bottle setter	1
	S999842	Rotor instruction manual	1

7. Decontamination

 **WARNING:**

- If the centrifuge, rotor or an accessory is contaminated by samples that are toxic or radioactive, or blood samples that are pathogenic or infectious, be sure to decontaminate the item according to good laboratory procedures and methods.
- If there is a possibility that the rotor or an accessory is contaminated by samples that might impair human health (for example, samples that are toxic or radioactive, or blood samples that are pathogenic or infectious), it is your responsibility to sterilize or decontaminate the rotor or the accessory properly before requesting repairs from an authorized sales or service representative. Note that we cannot repair the centrifuge, rotor or the accessory unless sterilization or decontamination is completed.
- It is your responsibility to sterilize and/or decontaminate the rotor or parts properly before returning them to an authorized sales or service representative. In such cases, copy the attached decontamination sheet and fill out the copied sheet, then attach it to the item to be returned. We may ask you about the treatment for the rotor or parts if the decontamination is checked and judged as insufficient by us. It is your responsibility to bear the cost of sterilization or decontamination. Note that we cannot repair or inspect the rotor or the accessory unless sterilization or decontamination is completed.

8. Rotor Retirement

After many years of use, there will be inevitably some corrosion or stress corrosion. At some points, the combination of such damage and metal fatigue could make the rotor vulnerable to a failure. Although a rotor may appear to be in a good condition, you should follow the rotor retirement recommendation shown below.

Rotor	Material	Retire After Years
Refrigerated centrifuge rotors	Aluminum alloy	15
	Stainless steel	

(Appendix) K Factors

- Calculation of separation time

The K factor can conveniently be used which makes it possible to estimate the settling time directly from Svedberg unit(S).

Namely, the time taken for particles of 1S to settle from Rmin. to Rmax. is K factor.
The K factor can be obtained from the formula show below.

$$K = \frac{\ln(R_{\max.}) - \ln(R_{\min.})}{\omega^2} \cdot \frac{10^{13}}{3600}$$

ω : Angular velocity $\omega = 2\pi/60 \times N$

N : Speed (rpm)

Rmin. : Distance from axis of rotation to top of solution (cm)

Rmax. : Distance from axis of rotation to bottom of tube (cm)

Using the K factor, the settling time t(in hours) is expressed as follows:

$$t = \frac{K}{S}$$

For example, when using 250PP bottle, the K factor at the maximum speed of 13,000rpm of Model R13A rotor is 1,111. The time taken for particles of 1000 S to be settled from R min. to R max. using this rotor is as follows:

$$t = \frac{1,111}{1,000} = 1.1 \text{ hours}$$

The time taken for settling of particles of 1000 S by the same rotor at a speed of 8,000 rpm is:

$$K = \left[\frac{13,000}{8,000} \right]^2 \times 1,111 = 2,934 \quad t = \frac{2,934}{1,000} = 2.9 \text{ hours}$$

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Copyright © 2022 Eppendorf Himac Technologies Co., Ltd. All rights reserved.

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