Keep this rotor instruction manual and the centrifuge manual in the file.

Angle Rotor for Ultracentrifuges

P32AT

Before using these rotors, please carefully read this instruction manual, "ROTORS, TUBES, BOTTLES AND CAPS", and the centrifuge instruction manual for its efficient operation and for your safety.

Keep this instruction manual for your reference and refer to it as required.

Contents

1. Specifications	_____	1
2. Applicable tubes/bottles	_____	2
3. How to use	_____	3
4. Maintenance	_____	5
5. Packing list	_____	7
6. Decontamination	_____	8
7. Rotor retirement	_____	8

Part No. { RLM rotor
9123150M
Rotor with optical adapter
91231500
Rotor with optical disk
9123151C

2022.01
S99801808
Original instruction manual

• The exterior or specification of this product is subject to change for improvement.

Safety Reminder

- Centrifuge rotors rotating at high speed have considerable potential for damage to personal properties if used improperly.
For safe and proper use of this rotor, carefully read the centrifuge instruction manual and this rotor instruction manual before use and observe the instructions.

 **WARNING :** and  **CAUTION :** notes are used to call your attention in this manual to prevent personal injury or damage to the rotor and the centrifuge. These notes are defined as follows.

 **WARNING:** indicates a potentially hazardous situation which, if not avoided, could result in personal severe injury or possible death.

 **CAUTION:** indicates a hazardous situation which, if not avoided, could result in personal injury or severe damage to the instrument.

WARNING

- Never use any material capable of producing flammable or explosive vapors. Your centrifuge and rotor have no explosion-proof construction.
- Your centrifuge and rotor are not designed to confine any sample particles dispersed due to leakage. Therefore, when using toxic or radioactive samples or pathogenic or infectious blood samples, make sure to prepare necessary safety measures at your own responsibility.
- Never exceed the maximum speed of the rotor (mentioned on the rotor).
Always reduce rotor speed as instructed in this manual when rotor speed is limited due to sample density or kinds of tubes.
- Check the chemical resistance chart attached the rotor, and do not use any sample inapplicable to the rotor. Using such a sample could corrode the rotor.
- Do not remove the RLM adapter or optical adapter/disk from rotor, or replace it with the adapter/disk for another rotor. The adapter/disk is critical component that detects the over-speed of rotor. If an adapter/disk that is not compatible with the rotor is attached, the rotor could break, resulting in damage to the centrifuge.
- Do not allow the rotor temperature to rise 100°C.: This would cause the material brittle.
- Never sterilize the rotor by autoclaving or boiling. Otherwise the strength of rotor may be significantly decreased.
- If the centrifuge, rotor, or an accessory is contaminated by samples that toxic or radioactive, or blood samples that pathogenic or infectious, be sure to decontaminate the item according to good laboratory procedures and methods.
- If there is a possibility that the centrifuge, rotor, or an accessory is contaminated by samples that might impair human health (for example, samples that are toxic or radioactive, or blood samples that are pathogenic or infectious), it is your responsibility to sterilize or decontaminate the centrifuge, rotor, or the accessory properly before requesting repairs from an authorized sales or service representative.
- It is your responsibility to sterilize and/or decontaminate the centrifuge, rotor, or parts properly before returning them to an authorized sales or service representative.

CAUTION

- Check the chemical resistance chart attached to the rotor, and do not use any sample inapplicable to the tubes/bottles, or the cap for the tubes/bottles. Using such a sample could deteriorate them.
- The allowable speed may be lower depending on the density of sample, or the combination of tubes or bottles and caps. Refer to this instruction manual.
- Always balance the tubes/bottles and samples within the allowable imbalance of rotor, including the caps. Do not exceed the allowable imbalance. Otherwise the rotor and the centrifuge may be damaged.
- Do not load only one tube or load tubes asymmetrically : Asymmetrically loading may cause imbalance operation and damage the centrifuge and rotor.
- Clean the inside of the drive hole (crown hole) of the rotor and the surface of the drive shaft (crown) of the centrifuge once a month. If the drive hole or the drive shaft is stained or any foreign matter is adhered, the rotor may be improperly installed and come off during operation.
- Use the rotor tubes and bottles within their actual capacities.
- Do not use tubes/bottles that have exceeded their life expectancy. Failure to do so could result in damage of tubes/bottles and the rotor and the centrifuge. The life expectancy of tubes/bottles depends on factors such as the characteristics of samples, speed of the rotor used, and temperature. Always check for deterioration and damage (cracks, deformation, and so on) on tubes/bottles before using them. Do not use the tubes/bottles if you find such a problem.
- Inspect and maintain the rotor after use. If abnormality is observed, do not use it.
Contact an authorized sales or service representative.

1. Specifications

1.1 Construction

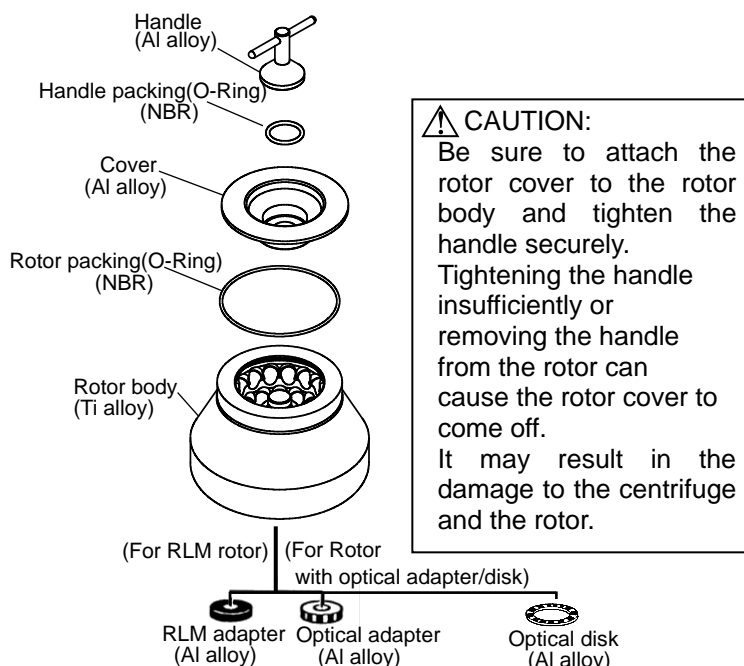
Fig.1-1. Construction of P32AT rotor

1.2 Specifications

- Max. speed: 32,000 rpm
- Max. RCF: 111,000 × g
- Allowable density at maximum speed of the rotor: 1.2g/mL
- K Factor at maximum speed of the rotor: 186
- Nominal rotor capacity: 12 mL × 32 tubes = 384mL
- Material of rotor: Ti alloy
- Material of cover: Al alloy
- Tube size: $\phi 1.6 \times 7.6$ cm
- Weight: 8.8 kg
- Max. diameter: 19.9cm
- Max. height: 19.1cm

CAUTION:

- Do not use a RLM rotor with the centrifuge other than a series of a CP-NX/WX/MX/ α ultracentrifuge.
- Do not use rotor with optical adapter/disk with model 55P and 40P (old type centrifuge).

NOTE:

When using a RLM rotor with a series of a CP-WX/MX/ α Ultracentrifuge, replacement of the ROM may be required (extra-cost option) depending on the serial number of your centrifuge. The serial number of your centrifuge is the SN or the MFG No. on the product label attached to the centrifuge.

1.3 Cross-sectional view of rotor

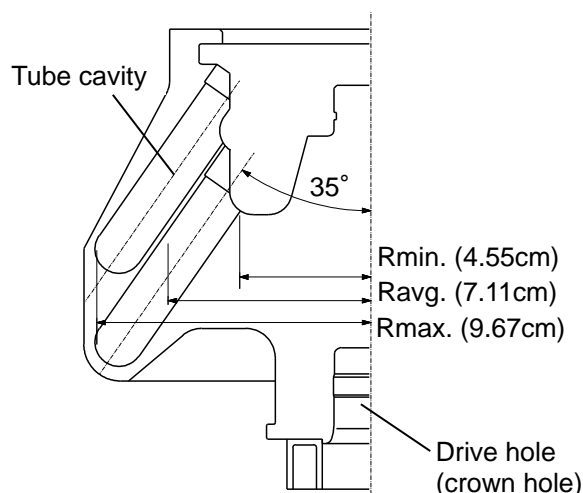
Fig.1-2. Cross-sectional view of P32AT rotor

CAUTION:

Clean the inside of the drive hole (crown hole) of the rotor and the surface of the drive shaft (crown) of the centrifuge once a month. If the drive hole or the drive shaft is stained or any foreign matter is adhered, the rotor may be improperly installed and come off during operation.

1.4 Characteristic of rotor

Table 1-1. Characteristic of P32AT

Rotor Speed (rpm)	RCF (×g)			K factor
	Rmax.	Ravg.	Rmin.	
5,000	2,700	1,990	1,270	7,639
10,000	10,800	7,950	5,090	1,910
15,000	24,300	17,900	11,400	849
20,000	43,200	31,800	20,300	477
25,000	67,600	49,700	31,800	306
30,000	97,300	71,500	45,800	212
32,000	111,000	81,400	52,100	186

※ For detail about K factor, refer to “ROTORS, TUBES, BOTTLES AND CAPS (Part No. S999204)” attached to the centrifuge.

Calculation formula

$$RCF = 1.118 \times 10^{-5} \times R \times N^2 (\times g)$$

R: Rotating radius (cm)

N: Speed (rpm)

2. Applicable tubes/bottles

2.1 Applicable tubes/bottles

Table 2-1. Applicable tubes/bottles

Nominal capacity (mL)	Tubes/Bottles				Cap		Max speed (rpm)	Max RCF (xg)
	Part No.	Name	Size (φ X L cm)	Actual capacity (mL/tube)	Part No.	Name		
12	329606A	12PA tube	1.6 X 7.6	10.2	S410542A	S-12AL cap ass'y	32,000	111,000
				10.5	441399A	C-AL cap ass'y	32,000	111,000
	325470A	12PE tube	1.6 X 7.5	10.0	S410542A	S-12AL cap ass'y	32,000	111,000
				10.5	441399A	C-AL cap ass'y	32,000	111,000
	325751A	12PC tube	1.6 X 7.6	10.5	441399A	C-AL cap ass'y	32,000	111,000
10	S303922A	10PC bottle(B)	1.6 X 8.2	8.4	---	---	32,000	111,000
	338455A	10PC thick-walled tube	1.6 X 7.7	6.3	---	---	32,000	111,000
	325952A	10PC bottle	1.6 X 8.2	8.0	---	---	32,000	111,000
	334105A	10PA bottle	1.6 X 8.2	8.0	---	---	32,000	111,000
	338456A	10PA thick-walled tube	1.6 X 7.6	6.3	---	---	32,000	111,000

1: When using 12PA tubes, we recommend the purchase of under-mentioned tube kit, what is the set of tubes and necessary tools.

Refer to Chapter 5 for the tube kit.

2: We will check the applicability of them upon request about tubes other than the above list. Contact an authorized sales or service representative.


CAUTION :

Chemical resistance and the strength of tubes/bottles varies with speed, temperature, and so on. Before using sample, fill tubes/bottles with water, buffer solution, etc. instead of sample and run them at the intended speed, temperature, and so on to ensure that there is no abnormality.

Refer to "ROTORS, TUBES, BOTTLES AND CAPS (Part No.S999204)" for the cleaning and sterilization of the tubes/bottles.

3. How to use

How to use tubes/bottles varies in the kinds of tubes/bottles. The brief procedure of tubes/bottles handling is described as follows.

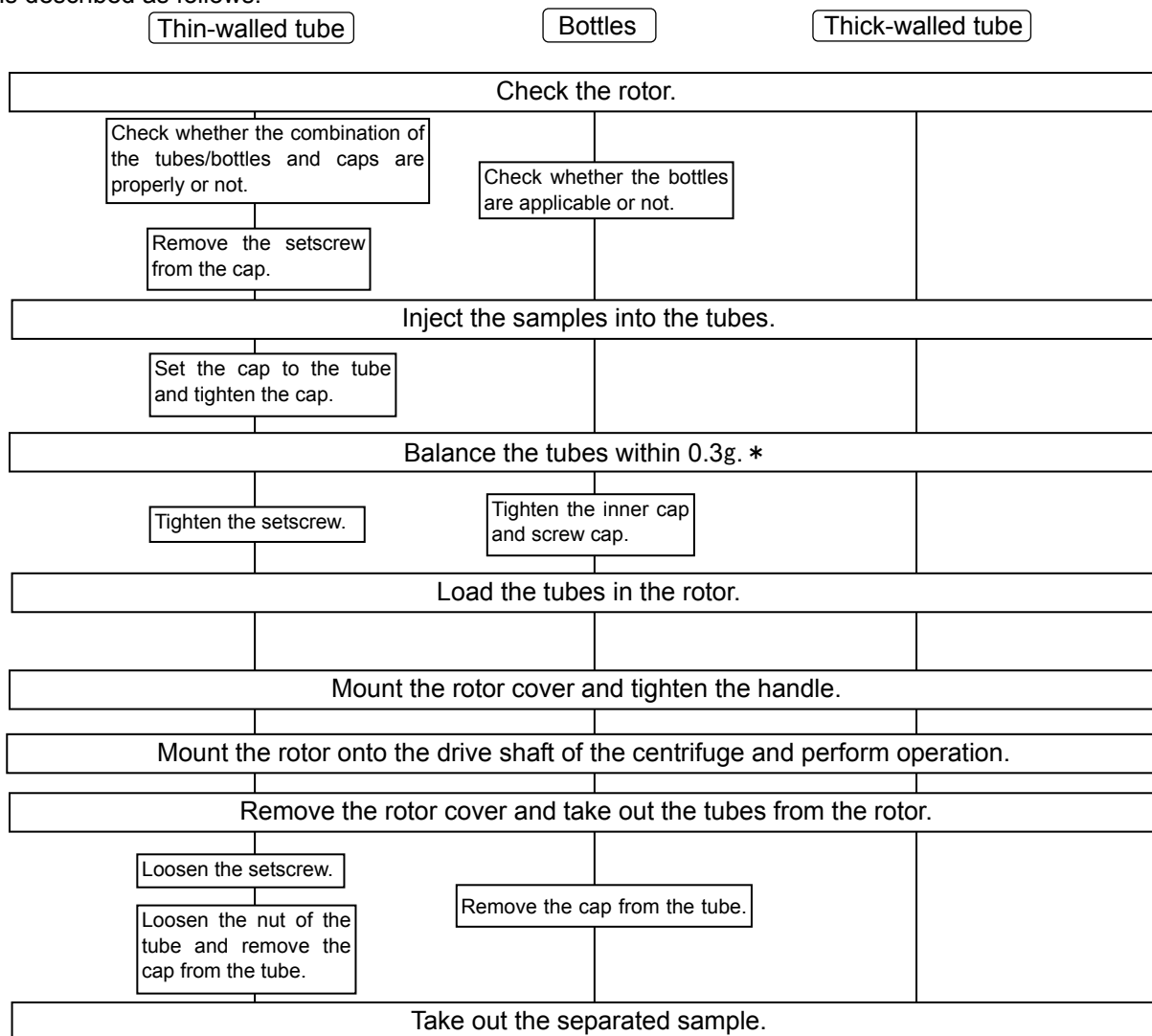(*) When using CP-NX series, CP-WX series, CP-MX series, or CP- α series, approximately even quantities of sample in the tubes are sufficient for balancing, and extremely different sample quantities must be avoided. (The levels of samples in the tubes should be approximately equal. Their difference should be within 5mm.)

For some kinds of tubes, if a small quantity of a sample is injected to them, the rotor can not be used at the allowable maximum speed. Tube deformation might occur due to a small quantity of a sample during operation.

=For handling=

● Allowable speed of rotors

1. To centrifuge a sample of an average density more than 1.2g/mL.

Determine the allowable speed from the following equation.

$$\text{Allowable speed(rpm)} = 32,000(\text{rpm}) \times \sqrt{\frac{1.2 (\text{g/mL})}{\text{Mean density of sample(g/mL)}}}$$

● Tube/Bottle

Before operation, check whether the combination of tubes/bottles, caps, etc. is properly or not.

When using thin-walled tubes, fill them with a sample. When using bottles and thick-walled tubes, use them with optional volume less than the actual capacity.

Tighten the cap securely. Refer to "2. Applicable tubes/bottles" in this manual and "ROTORS, TUBES, BOTTLES AND CAPS(Part No. S999204)".

=For handling=

● Handling the rotor


WARNING :

- Never exceed the maximum speed of the rotor(mentioned on the rotor). Otherwise, rotor may be damaged and result in damage to the centrifuge.
- Check the chemical resistance chart attached the rotor, and do not use any sample inapplicable to the rotor. Using such a sample could corrode the rotor.

CAUTION :

- Do not load only one tube or load tubes asymmetrically : Asymmetrically loading may cause imbalance operation and damage the centrifuge and rotor. This rotor has outside tube cavities and inside tube cavities. When you load the tube in the outside tube cavity, load the balanced other tube in the outside tube cavity symmetrically. When you load the tube in the inside tube cavity, load the balanced other tube in the inside tube cavity symmetrically.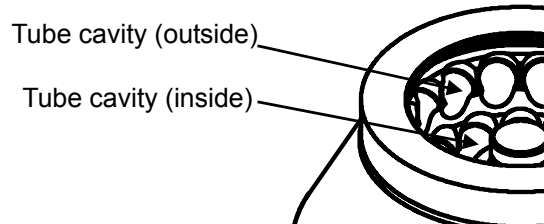

- Balance between symmetrically arranged tubes filled with the same-density sample.
- Balance the tubes and the samples within the allowable imbalance of the rotor.
Do not exceed the allowable imbalance. Otherwise, the centrifuge or the rotor may be damaged due to excessive imbalance operation.
- If foreign matter adheres to the rotor, wipe it off promptly. Otherwise, the rotor may be corroded.
- Be sure to attach the rotor cover to the rotor body and tighten the handle securely.
Tightening the handle insufficiently or removing the handle from the rotor can cause the rotor cover to come off and it might result in the damage to the centrifuge and the rotor.
- Check the packing (O-ring) every time. Replace it with new one if deterioration (crack, deformation, etc.) is observed. Otherwise, samples might be damaged due to poor rotor seal or the centrifuge or the rotor might be damaged during operation.
- Inspect and maintain the rotor after use. If abnormality is observed, do not use it.
Contact an authorized sales or service representative.

● Sample limitation

WARNING :

Never use any material capable of producing flammable or explosive vapors.
Your centrifuge and rotor are not designed to confine any sample particles dispersed due to leakage. Therefore, when using toxic or radioactive samples or pathogenic or infectious blood samples, make sure to prepare necessary safety measures at your own responsibility.

● The rotor life management of the rotor with optical adapter/disk

CAUTION :

Each time the rotor with optical adapter/disk is used, be sure to enter the results in the "rotor log book ". Failure to do so will disable management of the life of rotor. No entry is need for RLM rotor.

● RLM adapter and Optical adapter/disk

WARNING :

- Do not remove the RLM adapter or optical adapter/disk from rotor, or replace it with the adapter/disk for another rotor. The adapter/disk is a critical component that detects the over-speed of rotor : If an adapter/disk that is compatible with the rotor is attached, the rotor could break, resulting in damage to the ultracentrifuge.


CAUTION :

- Do not bring the RLM adapter near a magnet, or scratch it: Doing so will erase the memory stored in adapter, and make the rotor unusable. To prevent the adapter from being scratched, store the rotor with RLM adapter, using the stand provided with the rotor (rotor stand for protecting adapter). If the adapter is scratched, immediately stop using rotor and contact an authorized sales or service representative.
- Take great care not to scratch the optical adapter/disk: Any scratch on adapter/disk will make the rotor unusable. To prevent the adapter/disk from being scratched, store the rotor with optical adapter/disk, using the stand provided with the rotor (rotor stand for protecting adapter/disk). If the adapter/disk is scratched, immediately stop using rotor and contact an authorized sales or service representative.
- If foreign matter adheres to the RLM adapter and Optical adapter/disk, wipe it off promptly. Otherwise, the rotor can not be used.

4. Maintenance

● Maintenance of rotor

- ⚠ WARNING :**
Do not allow the rotor temperature to rise over 100°C.
Otherwise, the rotor can be brittle.
- ⚠ CAUTION :**
Use a neutral detergent having a pH between 5 and 9.
Otherwise, the rotor can be corroded.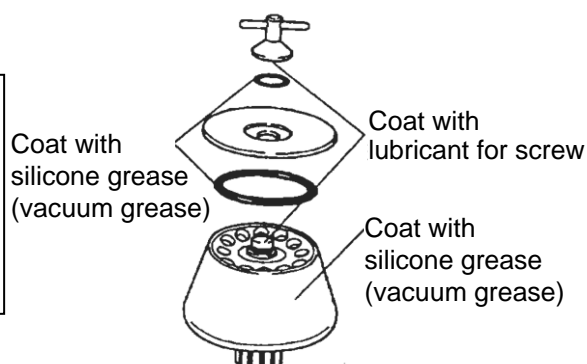

Wash the rotor with tap water or a dilute solution of neutral detergent and rinse it out with distilled water. Turn the rotor upside down with the rotor removed to dry it well. Check the rotor is completely dried and then put the light coat of silicone grease (vacuum grease) on the surface of the rotor, the light coat of lubricant for screw on the thread portion, the light coat of silicone grease (vacuum grease) on the rotor packing, Store them in a dry place. If foreign matter adheres to the rotor, soak the rotor in warm water in one or two hour and remove foreign matter with a soft brush etc. Use a neutral detergent having a pH between 5 and 9. If you cannot remove it, contact an authorized sales or service representative.

● Sterilizing rotor

- ⚠ WARNING :**
Never sterilize the rotor by autoclaving or boiling. Otherwise the strength of rotor may be significantly decreased.

Sterilize the rotor according to either gas sterilization method (ethylene oxide or formaldehyde), chemical sterilization method (70% ethanol, 3% hydrogen peroxide, 3% formalin), or ultraviolet rays sterilization.

* Do not dip the rotor and the tube adapters in the formalin (3%) solution more than 2 hours.

● Inspecting rotor

Check the rotor after every run as follows.

- The thread portion is worn. Contact an authorized sales or service representative.
- The tube cavity is corroded. Contact an authorized sales or service representative.
- The rotor packing is deteriorated. Replace it with the new one.
- The adapter is stained. Wipe it out with a soft cloth.

Check the rotor once a month as follows.

- The inside of drive hole (crown hole) is stained. Wipe it out with a soft cloth dampened with water.

- ⚠ CAUTION :**
Clean the inside of the drive hole (crown hole) of the rotor and the surface of the drive shaft (crown) of the centrifuge once a month. If the drive hole or the drive shaft is stained or any foreign matter is adhered, the rotor may be improperly installed and come off during operation.

● Life of rotor


While using rotor repeatedly, it's strength decreases gradually due to fatigue and creep of material, by which the rotor life is determined.

This rotor must not be used when they reached 5,000 operation times or 10,000 accumulated hours. If you use the RLM rotor at a speed lower than its maximum permissible speed, the life of the RLM rotor will be prolonged automatically by calculating from the strength of the rotor material.

● Replacing the optical disk

(1) Preparation

- Prepare a new optical disk and knife.
Make sure that the number of stripes of the new optical disk is right as follows.
- Wash and then dry the rotor well.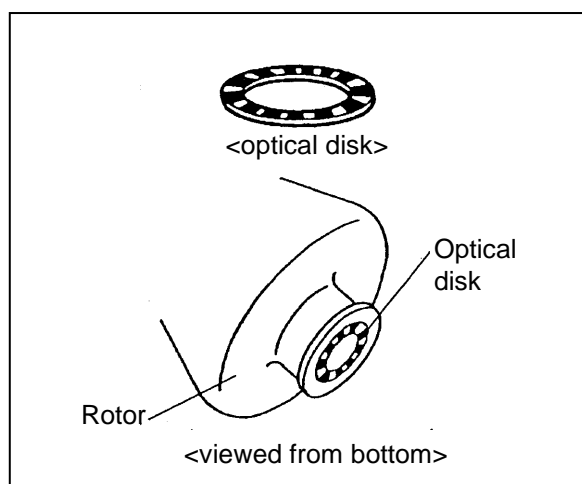

Rotor's maximum speed and optical disk's number of stripes

Rotor's maximum speed (rpm)	Optical disk's number of black bands	Rotor's maximum speed (rpm)	Optical disk's number of black bands	Rotor's maximum speed (rpm)	Optical disk's number of black bands
100,000	9	50,000	18	30,000	29
90,000	10	48,000	19	28,000	32
80,000~83,000	11	45,000	20	27,000	33
70,000	13	42,000	21	25,000	35
65,000	14	40,000	22	23,000	39
60,000	15	38,000	24	21,000	40
55,000~56,000	16	35,000	25	19,000	45
54,000	17	32,000	28	17,000	50

(2) Hold the rotor upside down.

(3) Pry the edge of the optical disk with the knife and remove the disk. Be very careful not to damage the rotor.

(4) Clean the disk hole on the rotor by using alcohol.

(5) Remove the backing paper from the new optical disk.

Place the optical disk so that it snugly fits into the groove of the disk hole on the rotor. Make sure the disk does not move.

5. Packing list

Check the parts and accessories of rotor when receiving the rotor.
Contact an authorized sales or service representative when the need arises.

No.	Name	Qty	Remark	Parts No. (Qty)	No.	Name	Qty	Parts No. (Qty)
①	Rotor	1			⑩	Rotor stand	1	216715 (1)
②	Handle	1*1			⑪	Lubricant for screw	1	84810601 (1)
③	Handle Packing (O-Ring)	1*1		433310 (1)	⑫	Silicone grease (vacuum grease)	1	483719 (1)
④	Cover	1*1			⑬	Tool box (C)	1	
⑤	Rotor Packing (O-Ring)	1*1		84520120 (1)	⑭			
⑥					⑮			
⑦	Optical disk *2	2	For spare	34644414	⑯			
⑧	Handle packing (O-Ring)	2	For spare		⑰	Rotor log book *3	1	999494 (1)
⑨	Rotor packing (O-Ring)	2	For spare		⑱	Instruction manual	1	S998018 (1)

*1: the quantities of the parts that are mounted to the rotor

*2: The "optical disk" is only provided with the rotor destined for Europe and U.S.A.

*3: The "rotor log book" is not included in RLM rotor.


Tubes are not included in this rotor. Please purchase the applicable tubes.

When using 12PA tubes, we recommend the purchase of under-mentioned tube kit, what is the set of tubes and necessary tools.

● 12PA tube kit (B) (Part No.S304986B)

(component)

329606A	12PA tube	2(50pcs/set)
441399A	C-AL Cap ass'y	12
S401796A	Packing for cap (gasket)	2(10pcs/set)
S300714A	Tube vise (C)	1
4361453	Box wrench	1
435823A	Tube setter (B)	1

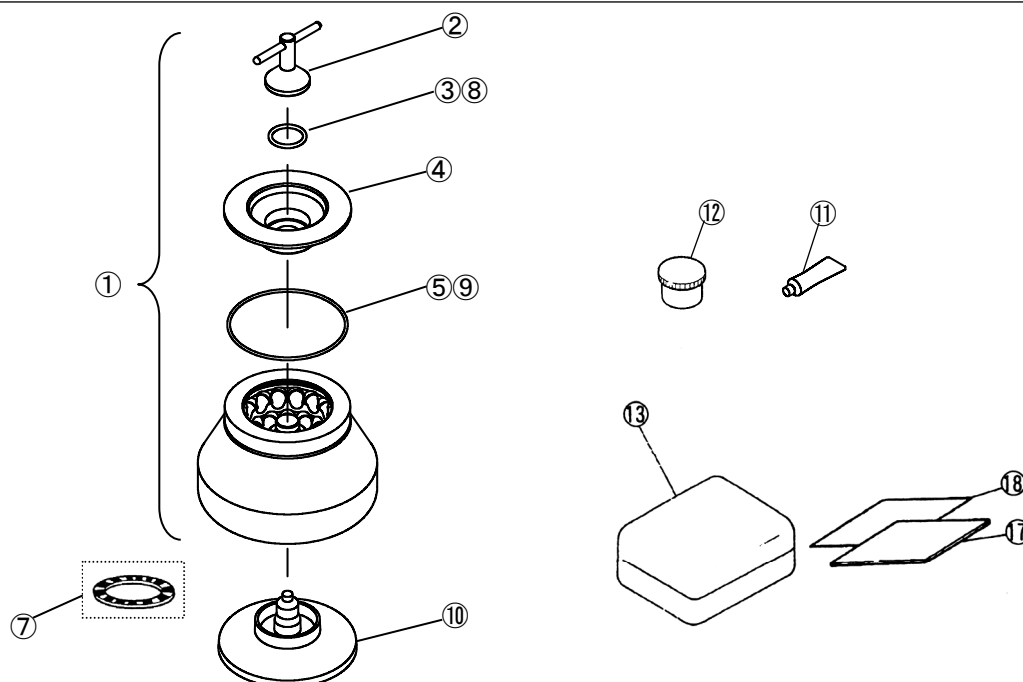6. Decontamination

⚠ WARNING:

- If the centrifuge, rotor or an accessory is contaminated by samples that are toxic or radioactive, or blood samples that are pathogenic or infectious, be sure to decontaminate the item according to good laboratory procedures and methods.
- If there is a possibility that the rotor or an accessory is contaminated by samples that might impair human health (for example, samples that are toxic or radioactive, or blood samples that are pathogenic or infectious), it is your responsibility to sterilize or decontaminate the rotor or the accessory properly before requesting repairs from an authorized sales or service representative. Note that we cannot repair the centrifuge, rotor or the accessory unless sterilization or decontamination is completed.
- It is your responsibility to sterilize and/or decontaminate the rotor or parts properly before returning them to an authorized sales or service representative. In such cases, copy the attached decontamination sheet and fill out the copied sheet, then attach it to the item to be returned. We may ask you about the treatment for the rotor or parts if the decontamination is checked and judged as insufficient by us. It is your responsibility to bear the cost of sterilization or decontamination. Note that we cannot repair or inspect the rotor or the accessory unless sterilization or decontamination is completed.

7. Rotor retirement

After many years of use, there will be inevitably some corrosion or stress corrosion. At some points, the combination of such damage and metal fatigue could make the rotor vulnerable to a failure. Although a rotor may appear to be in a good condition, you should follow the rotor retirement recommendation shown below.

Rotor	Material	Retire After Years
Angle rotor	Titanium alloy	12
Vertical rotor	Aluminum alloy	10
Swing rotor	Titanium alloy Aluminum alloy	10

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Copyright © 2022 Eppendorf Himac Technologies Co., Ltd. All rights reserved.

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