Keep this rotor instruction manual and the centrifuge manual in the file.

Vertical rotors for Ultracentrifuge

P100VT

Before using these rotors, please carefully read this instruction manual, "ROTORS, TUBES, BOTTLES AND CAPS", and the centrifuge instruction manual for its efficient operation and for your safety.

Keep this instruction manual for your reference and refer to it as required.

Contents

1. Specifications	_____	1
2. Applicable Tubes/Bottles	_____	2
3. How to Use	_____	3
4. Maintenance	_____	6
5. Using Cesium Chloride	_____	7
6. Packing List	_____	8
7. Decontamination	_____	9
8. Rotor Retirement	_____	9

Order no. 5720 213 001 (P100VT with RLM adapter)
Order no. 5720 213 104 (P100VT with Optical adapter)

2024.02
S99918613

Original instruction manual

Safety Reminder

- Centrifuge rotors rotating at high speed have considerable potential for damage to personal properties if used improperly.
For safe and proper use of this rotor, carefully read the centrifuge instruction manual and this rotor instruction manual before use and observe the instructions.

 **WARNING :** and  **CAUTION :** notes are used to call your attention in this manual to prevent personal injury or damage to the rotor and the centrifuge. These notes are defined as follows.

 **WARNING:** indicates a potentially hazardous situation which, if not avoided, could result in personal severe injury or possible death.

 **CAUTION:** indicates a hazardous situation which, if not avoided, could result in personal injury or severe damage to the instrument.

WARNING

- Never use any material capable of producing flammable or explosive vapors. Your centrifuge is not explosion proof centrifuge and your rotor is not explosion proof rotor.
- Your centrifuge and rotor are not designed to confine any sample particles dispersed due to leakage. Therefore, when using toxic or radioactive samples or pathogenic or infectious blood samples, make sure to prepare necessary safety measures at your own responsibility.
- Never exceed the maximum speed of the rotor(mentioned on the rotor).
Always reduce rotor speed as instructed in this manual when rotor speed is limited due to sample density or kinds of tubes.
- Check the chemical resistance chart attached the rotor, and do not use any sample inapplicable to the rotor. Using such a sample could corrode the rotor.
- Do not remove the RLM adapter or optical adapter from rotor, or replace it with the adapter for another rotor. The adapter is critical component that detects the over-speed of rotor. If an adapter that is not compatible with the rotor is attached, the rotor could break, resulting in damage to the centrifuge.
- Do not allow the rotor temperature to rise 100°C.: This would cause the material brittle.
- Never sterilize the rotor by autoclaving or boiling. Otherwise the strength of rotor may be significantly decreased.
- If the centrifuge, rotor, or an accessory is contaminated by samples that toxic or radioactive, or blood samples that pathogenic or infectious, be sure to decontaminate the item according to good laboratory procedures and methods.
- If there is a possibility that the centrifuge, rotor, or an accessory is contaminated by samples that might impair human health (for example, samples that are toxic or radioactive, or blood samples that are pathogenic or infectious), it is your responsibility to sterilize or decontaminate the centrifuge, rotor, or the accessory properly before requesting repairs from an authorized sales or service representative.
- It is your responsibility to sterilize and/or decontaminate the centrifuge, rotor, or parts properly before returning them to an authorized sales or service representative.

CAUTION

- Check the chemical resistance chart attached to the rotor, and do not use any sample inapplicable to the tubes or the caps for the tubes. Using such a sample could deteriorate them.
- The allowable speed may be lower depending on the density of sample or use of salt-such as cesium chloride-. Refer to this instruction manual.
- Always balance the tubes and samples within the allowable imbalance of rotor, including the caps. Do not exceed the allowable imbalance. Otherwise the rotor and the centrifuge may be damaged.
- Do not load only one tube or load tubes asymmetrically : Asymmetrically loading may cause imbalance operation and damage the centrifuge and rotor.
- Clean the inside of the drive hole (crown hole) of the rotor and the surface of the drive shaft (crown) of the centrifuge once a month. If the drive hole or the drive shaft is stained or any foreign matter is adhered, the rotor may be improperly installed and come off during operation.
- Do not use tubes that have exceeded their life expectancy. Failure to do so could result in damage of tubes and the rotor and the centrifuge. The life expectancy of tubes depends on factors such as the characteristics of samples, speed of the rotor used, and temperature. Always check for deterioration and damage (cracks, deformation, and so on) on tubes before using them. Do not use the tubes if you find such a problem.
- Inspect and maintain the rotor after use. If abnormality is observed, do not use it. Contact an authorized sales or service representative.

1. Specifications

1.1 Construction

⚠ CAUTION:
If the number of tubes to be set is small, do not install the space cap or the cap in the tube cavity into which no tube is inserted, but stay the cavities uninstalled for operation.

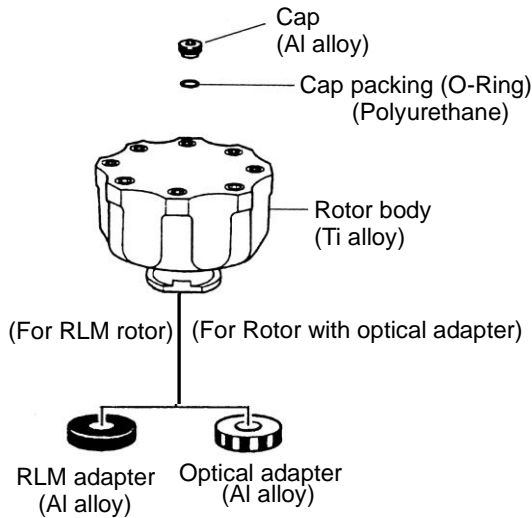
Fig.1-1. Construction of P100VT rotor

1.2 Specifications

- Max. speed: 100,000 rpm
- Max. RCF: 700,000 × g
- Allowable density at maximum speed of the rotor: 1.7g/mL
- K Factor at maximum speed of the rotor: 6
- Nominal rotor capacity: 5 mL × 8 tubes = 40 mL
- Material of rotor: Ti alloy
- Tube size: φ 1.3 × 5.3 cm
- Weight: 4.0 kg
- Max. diameter: 13.7 cm
- Max. height: 12.0 cm

⚠ CAUTION:

- Do not use a RLM rotor with the centrifuge other than a series of a CP-NX/WX/MX/α ultracentrifuge.
- Do not use rotor with optical adapter with model 55P and 40P (old type centrifuge).

1.3 Cross-sectional view of rotor

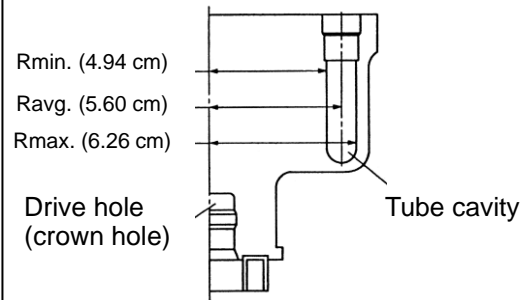
Fig.1-2. Cross-sectional view of P100VT rotor

⚠ CAUTION:
Clean the inside of the drive hole (crown hole) of the rotor and the surface of the drive shaft (crown) of the centrifuge once a month. If the drive hole or the drive shaft is stained or any foreign matter is adhered, the rotor may be improperly installed and come off during operation.

1.4 Characteristic of rotor

Table 1-1. Characteristic of P100VT

Rotor Speed (rpm)	RCF(xg)			K factor
	Rmax.	Ravg.	Rmin.	
10,000	7,000	6,260	5,520	600
20,000	28,000	25,000	22,100	150
30,000	63,000	56,300	49,700	67
40,000	112,000	100,000	88,400	37
50,000	175,000	157,000	138,000	24
60,000	252,000	225,000	199,000	17
70,000	343,000	307,000	271,000	12
80,000	448,000	401,000	353,000	9
90,000	567,000	507,000	447,000	7
100,000	700,000	626,000	552,000	6

※ For detail about K factor, refer to “ROTORS, TUBES, BOTTLES AND CAPS (Part No.S999204)” attached to the centrifuge.

Calculation formula

$$RCF = 1.118 \times 10^{-5} \times R \times N^2 (\times g)$$

R: Rotating radius (cm)
N: Speed (rpm)

2. Applicable Tubes/Bottles

2.1 Applicable Tubes/Bottles

Table 2-1. Applicable Tubes/Bottles

Nominal capacity (mL)	Tubes/Bottles				Tube adapter		Cap		Max. speed (rpm)	Max. RCF (xg)
	Order no.	Name	Size (ϕ X L cm)	Actual capacity (mL/tube)	Order no.	Name	Order no.	Name		
5	5720 411 008	5PP seal tube	1.3 X 5.3	5.1	---	---	5720 411 397	B2 space cap	100,000	700,000

1. When using 5PP seal tubes, we recommend the purchase of the tube kit, what is the set of tubes and necessary tools. Refer to Chapter 6 for the tube kit.

CAUTION :


Chemical resistance and the strength of tubes varies with speed, temperature, and so on. Before using sample, fill tubes with water, buffer solution, etc. instead of sample and run them at the intended speed, temperature, and so on to ensure that there is no abnormality.

Refer to "ROTORS, TUBES, BOTTLES AND CAPS (Part No. S999204)" for the cleaning and sterilization of the tubes/bottles.

3. How to use

How to use tubes/bottles varies in the kinds of tubes/bottles. The brief procedure of tubes/bottles handling is described as follows.

For the detail about handling tubes/bottles, refer to "ROTORS, TUBES, BOTTLES AND CAPS (Part No.S999204)" attached to the centrifuge.

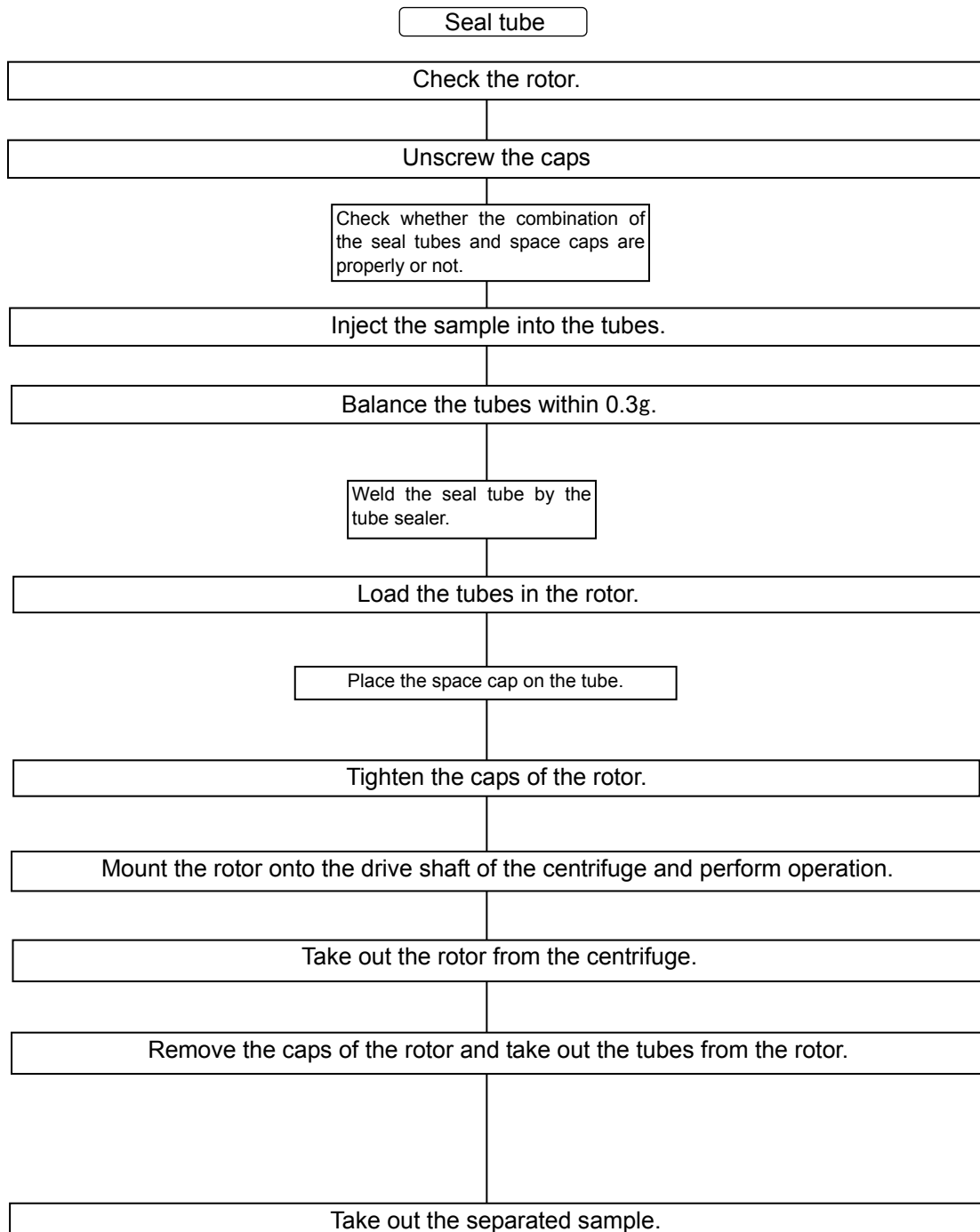=For handling=

● Allowable speed of rotors

1. To centrifuge a sample of an average density more than 1.7g/mL .
Determine the allowable speed from the following equation.

$$\text{Allowable speed(rpm)} = 100,000(\text{rpm}) \times \sqrt{\frac{1.7(\text{g/mL})}{\text{Mean density of sample(g/mL)}}}$$

2. The allowable speed may be lower depending on the density of a sample when using of salt-such as cesium chloride. Refer to "5. Using Cesium Chloride" for the detail.

● Tube

Before operation, check whether the combination of tubes, space caps, and spacers is properly or not. When using seal tubes, fill them with samples.

Weld the seal tube completely according to the instruction manual of the STF-1, the STF2, or the STF3 tube sealer.

Refer to "2. Applicable tubes/bottles" in this manual and "ROTORS, TUBES, BOTTLES AND CAPS (Part No. S999204)".

● Handling the rotor

WARNING :

- Never exceed the maximum speed of the rotor(mentioned on the rotor). Otherwise, rotor may be damaged and result in damage to the centrifuge.
- Check the chemical resistance chart attached the rotor, and do not use any sample inapplicable to the rotor. Using such a sample could corrode the rotor.


CAUTION :

- Do not load only one tube or load tubes asymmetrically : Asymmetrically loading may cause imbalance operation and damage the centrifuge and rotor.
- Balance the tubes and the samples including the caps within the allowable imbalance of the rotor. Do not exceed the allowable imbalance. Otherwise, the centrifuge or the rotor may be damaged due to excessive imbalance operation.
- If foreign matter adheres to the rotor, wipe it off promptly. Otherwise, the rotor may be corroded.
- If the number of tubes to be set is small, do not install the space cap or the cap in the tube cavity into which no tube is inserted, but stay the cavities uninstalled for operation.
- Inspect and maintain the rotor after use. If abnormality is observed, do not use it. Contact an authorized sales or service representative.

Load the balanced tubes symmetrically with the respect to the center of the rotor.
Tighten the caps of the rotor securely (see the figure below).

• Using seal tubes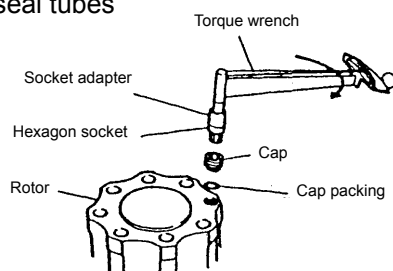

<How to use tighten the caps of the rotor>

1. Fix the holder base (one of accessories) to a desk, etc.
2. Fix the rotor to the ditch of the holder base.
3. Place the cap on the rotor and tighten to 12N·m with the torque wrench.

$$*1\text{N} \cdot \text{m} = 10.2\text{kgf} \cdot \text{cm}$$

Do the maintenance of the rotor and check the rotor after every use.

● Sample limitation

 WARNING :

- Never use any material capable of producing flammable or explosive vapors. Your centrifuge is not explosion proof centrifuge and your rotor is not explosion proof rotor.
- When using toxic or radioactive samples or pathogenic or infectious blood samples, make sure to prepare necessary safety measures at your own responsibility.

● The rotor life management of the rotor with optical adapter

 CAUTION :

Each time the rotor with optical adapter is used, be sure to enter the results in the “rotor log book”. Failure to do so will disable management of the life of rotor.
No entry is need for RLM rotor.

● RLM adapter and Optical adapter

 WARNING :

- Do not remove the RLM adapter or optical adapter from rotor, or replace it with the adapter for another rotor. The adapter is a critical component that detects the over-speed of rotor:
If an adapter that is compatible with the rotor is attached, the rotor could break, resulting in damage to the ultracentrifuge.


 CAUTION :

- Do not bring the RLM adapter near a magnet, or scratch it: Doing so will erase the memory stored in adapter, and make the rotor unusable. To prevent the adapter from being scratched, store the rotor with RLM adapter, using the stand provided with the rotor (rotor stand for protecting adapter).
If the adapter is scratched, immediately stop using rotor and contact an authorized sales or service representative.
- Take great care not to scratch the optical adapter: Any scratch on adapter will make the rotor unusable. To prevent the adapter from being scratched, store the rotor with optical adapter, using the stand provided with the rotor (rotor stand for protecting adapter). If the adapter is scratched, immediately stop using rotor and contact an authorized sales or service representative.

4. Maintenance

● Maintenance of rotor

- ⚠ WARNING :**
Do not allow the rotor temperature to rise over 100°C.
Otherwise, the rotor can be brittle.
- ⚠ CAUTION :**
Use a neutral detergent having a pH between 5 and 9.
Otherwise, the rotor can be corroded.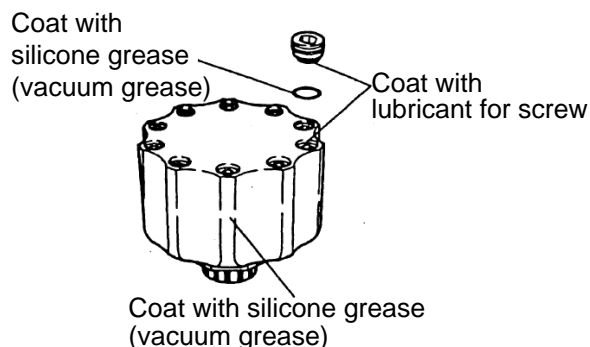

Wash the rotor with tap water or a dilute solution of neutral detergent and rinse it out with distilled water. Turn the rotor upside down with the caps removed to dry it well. Check the rotor is completely dried and then put the light coat of silicone grease (vacuum grease) on the surface of the rotor, the light coat of lubricant for screw on the thread portion, the light coat of silicone grease (vacuum grease) on the rotor packing, Store them in a dry place. If foreign matter adheres to the rotor, soak the rotor in warm water in one or two hour and remove foreign matter with a soft brush etc. Use a neutral detergent having a pH between 5 and 9. If you can not remove it, contact an authorized sales or service representative.

● Sterilizing rotor

- ⚠ WARNING :**
Never sterilize the rotor by autoclaving or boiling. Otherwise the strength of rotor may be significantly decreased.

Sterilize the rotor according to either gas sterilization method (ethylene oxide or formaldehyde) or chemical sterilization method (70% ethanol, 3% hydrogen peroxide, 3% formalin).

* Do not dip the rotor in the formalin (3%) solution more than 2 hours.

● Inspecting rotor

Check the rotor after every run as follows.

- The thread portion is worn. Contact an authorized sales or service representative.
- The tube cavity is corroded. Contact an authorized sales or service representative.
- The packing is deteriorated. Replace it with the new one.
- The adapter is stained. Wipe it out with a soft cloth.

Check the rotor once a month as follows.

- The inside of drive hole(crown hole) is stained. . . . Wipe it out with a soft cloth dampened with water.

- ⚠ CAUTION :**
Clean the inside of the drive hole (crown hole) of the rotor and the surface of the drive shaft (crown) of the centrifuge once a month. If the drive hole or the drive shaft is stained or any foreign matter is adhered, the rotor may be improperly installed and come off during operation.

● Life of rotor

While using rotor repeatedly, it's strength decreases gradually due to fatigue and creep of material, by which the rotor life is determined.

This rotor must not be used when they reached 5,000 operation times or 10,000 accumulated hours. If you use the RLM rotor at a speed lower than its maximum permissible speed, the life of the RLM rotor will be prolonged automatically by calculating from the strength of the rotor material.

5. Using Cesium Chloride

The cesium chloride (CsCl) solution is frequently used as a density gradient medium, but if the CsCl solution with a high density is used, it may saturate during the rotation of the rotor depending on the rotation condition which may cause CsCl crystals to be educed. Educated crystals have a high density (approx. 4g/mL) and apply an excessive load to the rotor which is very dangerous.

The crystallization will also greatly affect the density gradient of the CsCl solution and separation status of the sample. Therefore, always use a solution which will not crystallize the allowable speed.

Fig.5-1 shows the relationship between the maximum allowable density and the speed at which cesium chloride does not crystallize in the condition that standard tubes is filled with the cesium chloride solution. In this diagram, crystallization will occur above the curves.

Be sure to select the density of the cesium chloride solution and the rotor speed at which cesium chloride does not crystallize. The curve in this diagram vary in temperature etc. Before actual operation, centrifuge the cesium chloride solution with the same density to ensure that crystallization will not occur.

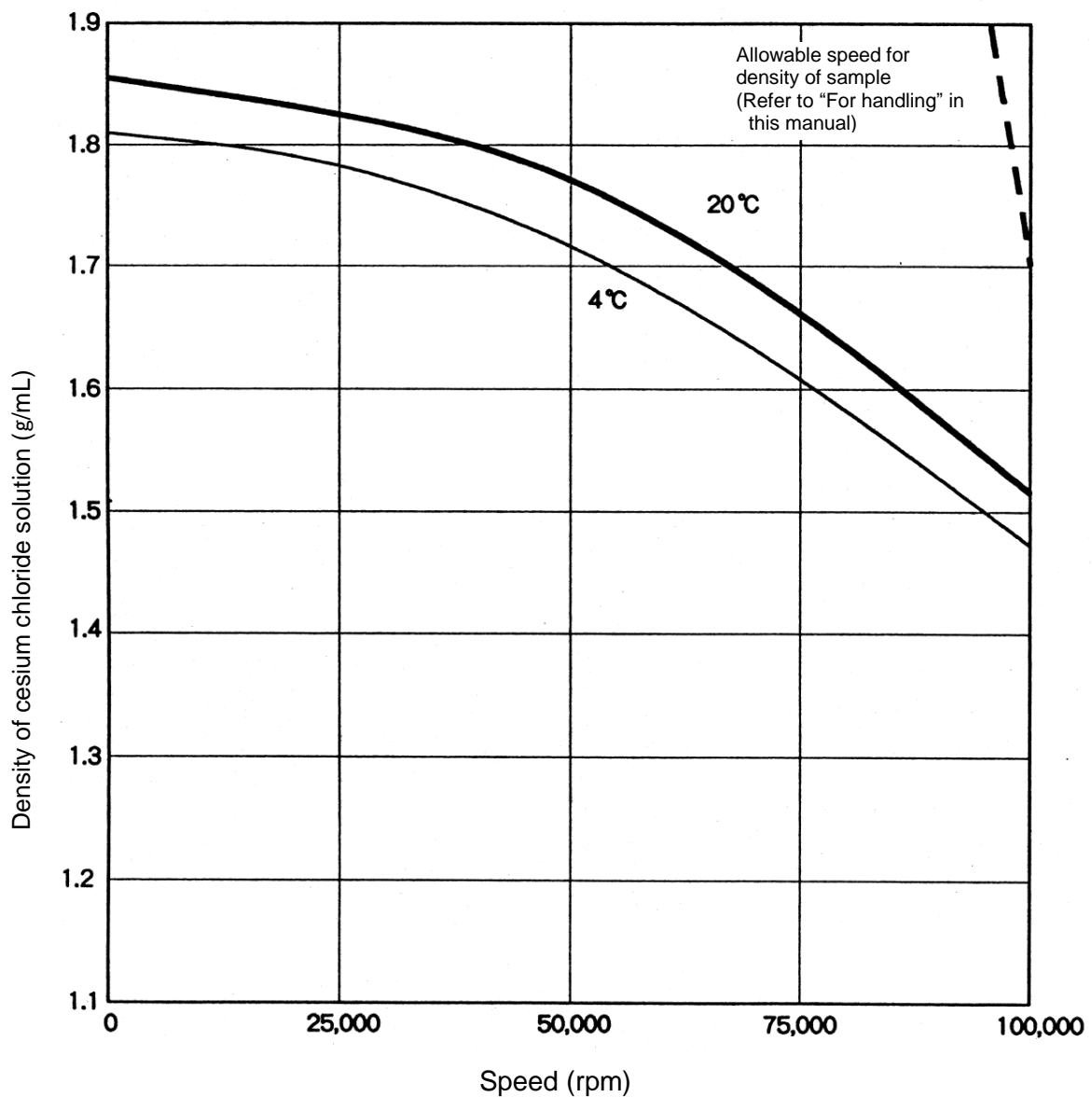
Fig.5-1. Relationship between density of cesium chloride and speed

6. Packing list

Check the parts and accessories of rotor when receiving the rotor.
Contact an authorized sales or service representative when the need arises.

No.	Name	Qty	Remark	Order no. (Qty)	No.	Name	Qty	Order no. (Qty)
①	Rotor	1			⑪	Lubricant for screw	1	5720 411 230 (1)
②	Cap (D)	8*1			⑫	Silicone grease (vacuum grease)	1	5720 850 317 (1)
③		Cap packing (D) (O-Ring)	8*1	5720 411 277 (10)	⑬	Tool box(C)	1	
④					⑭	Torque wrench	1	5720 411 282 (1)
⑤					⑮	Socket adapter		5720 411 283 (1)
⑥	Cap packing (D) (O-Ring)	16	For spare		⑯	Hexagon socket	1	
⑦					⑰	Holder base	1	5721 850 368 (1)
⑧					⑱	Rotor log book *2	1	5720 411 287 (1)
⑨					⑲	Instruction manual	1	
⑩	Rotor stand	1		5720 410 232 (1)	⑳			

*1: the quantities of the parts that are mounted to the rotor

*2: The "rotor log book" is not included in RLM rotor.

Tubes are not included in this rotor. Please purchase the applicable tubes.

When using 5PP seal tubes, we recommend the purchase of under-mentioned tube kit, what is the set of tubes and necessary tools.

※ When using 5PP seal tubes, Tube sealer (STF3, STF2, STF-1) and Tube rack (B) (Order no. 5720 410 004) are needed.


● 5PP seal tube kit (A) (Order no. 5720 411 070)

(component)

5720 411 008 5PP seal tube : 2(50pcs/set)

5720 411 397 B2-space cap : 8

5720 411 232 Tube setter : 1

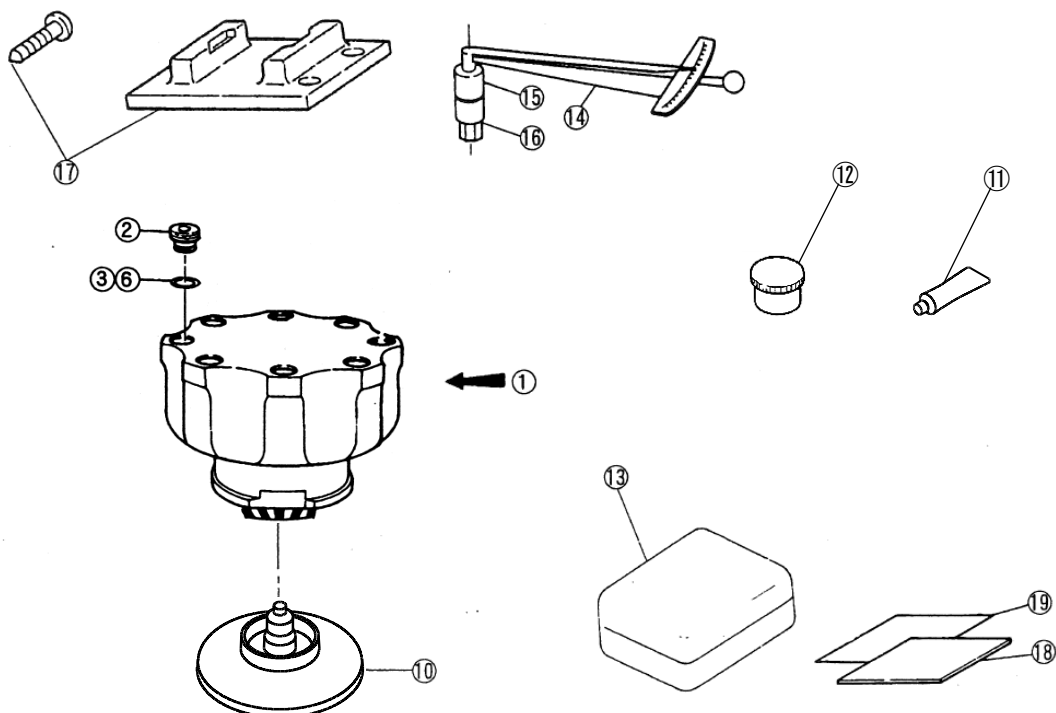7. Decontamination

⚠ WARNING:

- If the centrifuge, rotor or an accessory is contaminated by samples that are toxic or radioactive, or blood samples that are pathogenic or infectious, be sure to decontaminate the item according to good laboratory procedures and methods.
- If there is a possibility that the rotor or an accessory is contaminated by samples that might impair human health (for example, samples that are toxic or radioactive, or blood samples that are pathogenic or infectious), it is your responsibility to sterilize or decontaminate the rotor or the accessory properly before requesting repairs from an authorized sales or service representative. Note that we cannot repair the centrifuge, rotor or the accessory unless sterilization or decontamination is completed.
- It is your responsibility to sterilize and/or decontaminate the rotor or parts properly before returning them to an authorized sales or service representative. In such cases, copy the attached decontamination sheet and fill out the copied sheet, then attach it to the item to be returned. We may ask you about the treatment for the rotor or parts if the decontamination is checked and judged as insufficient by us. It is your responsibility to bear the cost of sterilization or decontamination. Note that we cannot repair or inspect the rotor or the accessory unless sterilization or decontamination is completed.

8. Rotor Retirement

After many years of use, there will be inevitably some corrosion or stress corrosion. At some points, the combination of such damage and metal fatigue could make the rotor vulnerable to a failure. Although a rotor may appear to be in a good condition, you should follow the rotor retirement recommendation shown below.

Rotor	Material	Retire After Years
Angle rotor	Titanium alloy	12
Vertical rotor	Aluminum alloy	10
Swing rotor	Titanium alloy Aluminum alloy	10
Zonal rotor	Titanium alloy	10
Continuous flow rotor	Titanium alloy	10

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Copyright © 2024 Eppendorf Himac Technologies Co., Ltd. All rights reserved.

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