

## Installation guide

epBlue™ GxP

English (EN)

This information is a supplement to the operating manuals of epBlue and epMotion and it does not replace these operating manuals. The epBlue GxP modification kit is an extension to epBlue version 40.x and higher.

## 1 Activate licenses

For the epBlue GxP modification kit it is necessary to activate a computer-specific license to run the software functionality.


### 1.1 Request an activation license

Prerequisites

- epMotion and MultiCon were installed by a service technician.
- A USB-storage medium is connected to the MultiCon.
- A user belonging to the user group *Administrator Lab* is logged in.

1. Click on the symbol *Settings* on the start screen of epBlue.
2. Select the tab *License Management*.
3. Press the button *Show computer information*.

The window *Computer Identity Information* opens.

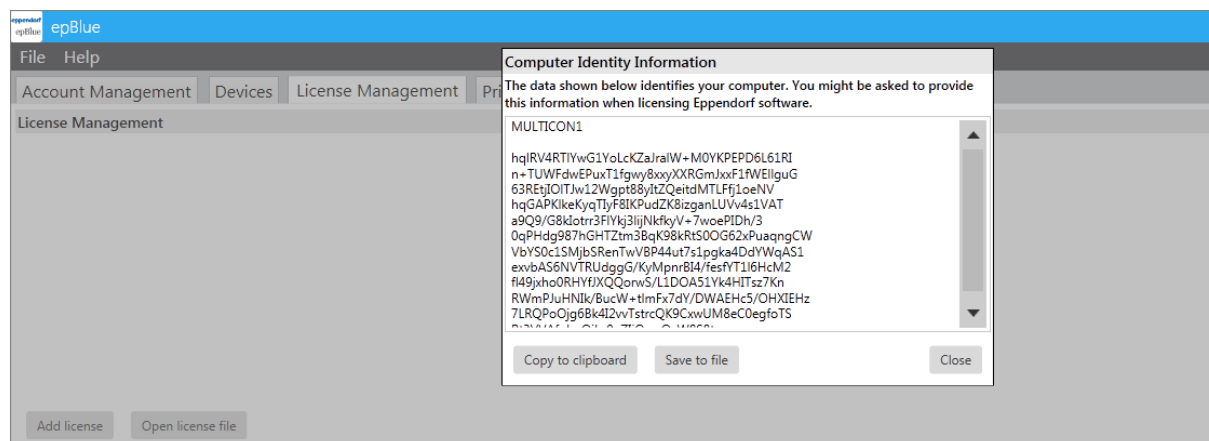

4. Press the button *Save to file*.
5. Select a directory on the USB storage medium.
6. Press the button *Save*.  
A text file containing the computer identity information is saved.
7. To request your epBlue GxP license, register on webpage [www.eppendorf.com/software-licensing](http://www.eppendorf.com/software-licensing).
8. Copy the content of the text file from the USB storage medium to the appropriate field on the webpage.




You will receive an epBlue GxP license file per e-mail within 48 hours.

## Installation guide

epBlue™ GxP

English (EN)

### 1.2 Import license file

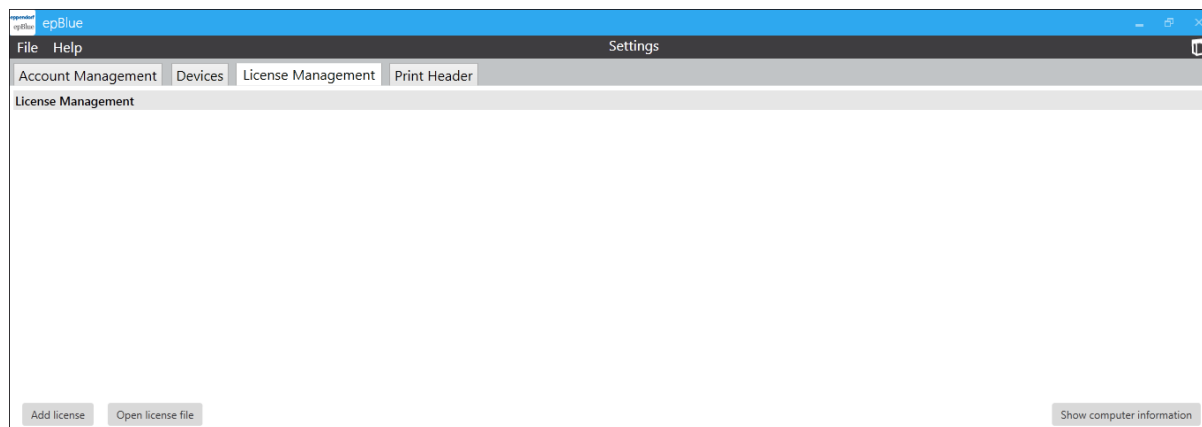

#### Prerequisites

- A user belonging to the user group *Administrator Lab* is logged in.
- The tab *Settings > License Management* is open.
- The USB storage medium containing the epBlue GxP license is connected to the MultiCon.

1. Press the button *Open license file*.


The window *Open Eppendorf License* opens.

2. Select the license to be imported.

3. Press the button *Open*.

License is inserted.

4. Restart epBlue.

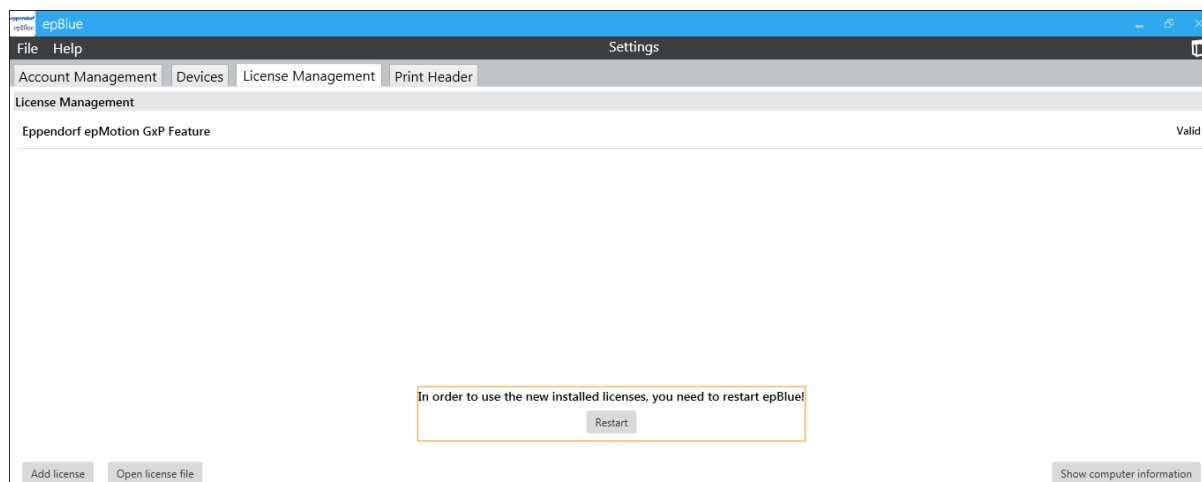

**Your local distributor: [www.eppendorf.com/contact](http://www.eppendorf.com/contact)**

Eppendorf AG · 22331 Hamburg · Germany

[eppendorf@eppendorf.com](mailto:eppendorf@eppendorf.com) · [www.eppendorf.com](http://www.eppendorf.com)