New Brunswick™ TCA-3 Temperature Monitoring System

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Copyright

Copyright © 2014 Eppendorf AG, Germany. No part of this publication may be reproduced without the prior permission of the copyright owner.

The company reserves the right to change information in this document without notice. Updates to information in this document reflect our commitment to continuing product development and improvement.

Trademarks

Eppendorf® and the Eppendorf logo are registered trademarks of Eppendorf AG, Germany.

New Brunswick™ is a trademark of Eppendorf AG, Germany.

VeriSign Secured® is a registered trademark of VeriSign, Inc., USA

Microsoft® and Excel® are registered trademarks of Microsoft Corporation in the United States and other countries.

Trademarks are not marked in all cases with ™ or ® in this manual.

Eppendorf has attempted to identify the ownership of all trademarks from public records. Any omissions or errors are unintentional.

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Operação

TCA-3 Temperature Monitoring System
Brazilian Portuguese (BR)

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The image shows a login interface for New Brunswick. On the left, there is a form with the following fields: "Gateway Mac:" with a text input field, "Login Name:" with a text input field, "Password:" with a text input field, and a checkbox labeled "Remember Gateway Mac". Below these fields is a "Login" button. A red oval highlights the "Gateway Mac", "Login Name", "Password", and "Remember Gateway Mac" fields. To the right of the form, the text "Monitoring made easy." is displayed. At the bottom left of the form area, there is a link: "Click [HERE](#) to try the demo site."

Fig. 4-2: Visor de entrada

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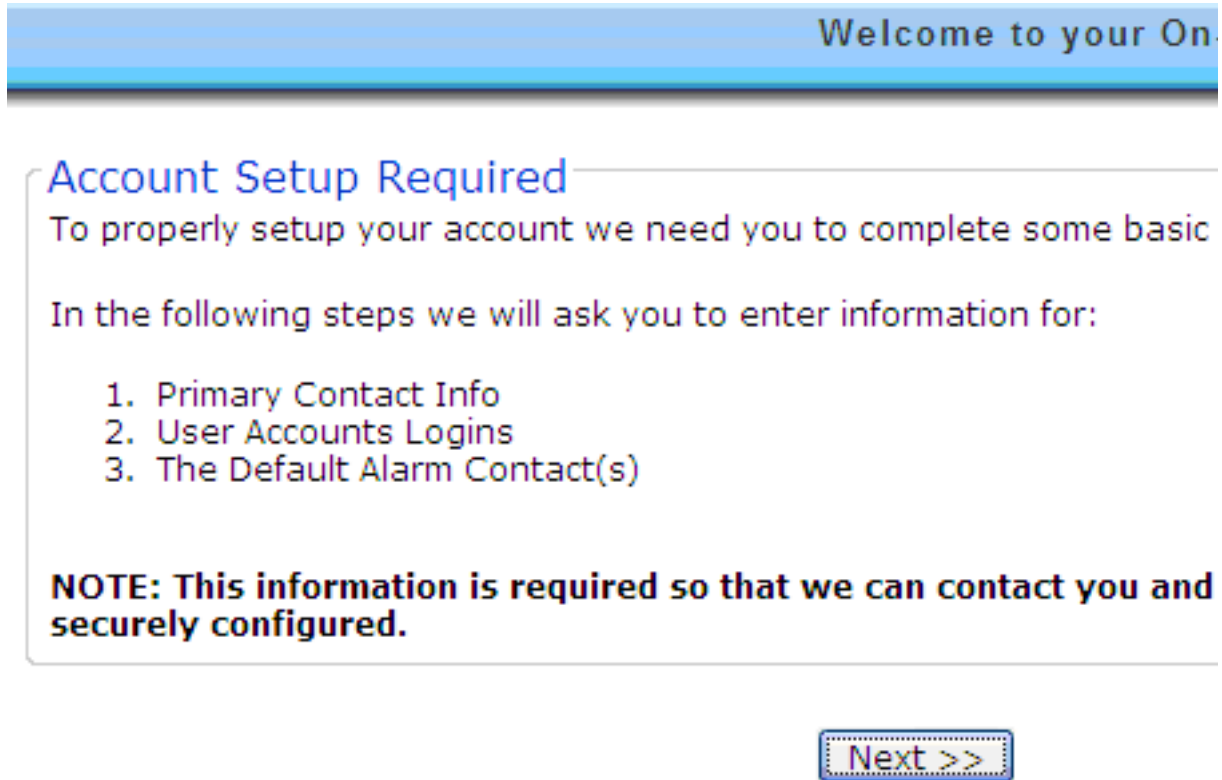
The screenshot shows a web page with a blue header bar containing the text "Welcome to your On...". Below the header, there is a section titled "Account Setup Required" in blue text. The text below the title reads: "To properly setup your account we need you to complete some basic". This is followed by the sentence "In the following steps we will ask you to enter information for:". Below this, there is a numbered list with three items: "1. Primary Contact Info", "2. User Accounts Logins", and "3. The Default Alarm Contact(s)". At the bottom of the section, there is a bold note: "NOTE: This information is required so that we can contact you and securely configured.". At the bottom right of the page, there is a button with a blue border and a dotted background, containing the text "Next >>".

Welcome to your On...

Account Setup Required

To properly setup your account we need you to complete some basic

In the following steps we will ask you to enter information for:

1. Primary Contact Info
2. User Accounts Logins
3. The Default Alarm Contact(s)

NOTE: This information is required so that we can contact you and securely configured.

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Account Se

Primary Contact

From time to time we may need to notify you of system updates, schedules, and other important information. Who is the primary contact for this monitoring system account?

Who is the primary contact for this monitoring system account?

Full Name*:	<input type="text"/>
Company*:	<input type="text"/>
Address Line1*:	<input type="text"/> <small>Street address, P.O. box, company name, c/o</small>
Address Line2:	<input type="text"/> <small>Apartment, suite, unit, building, floor, etc.</small>
City*:	<input type="text"/>
State/Province/Region*:	<input type="text"/>
ZIP/Postal Code*:	<input type="text"/>
Country*:	<input type="text" value="United States"/>
Timezone*:	<input type="text" value="(GMT-05:00) Eastern Standard Time (DST)"/>
Phone Number*:	<input type="text"/>
Email Address*:	<input type="text"/>
Repeat Email Address*:	<input type="text"/>

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Account Se

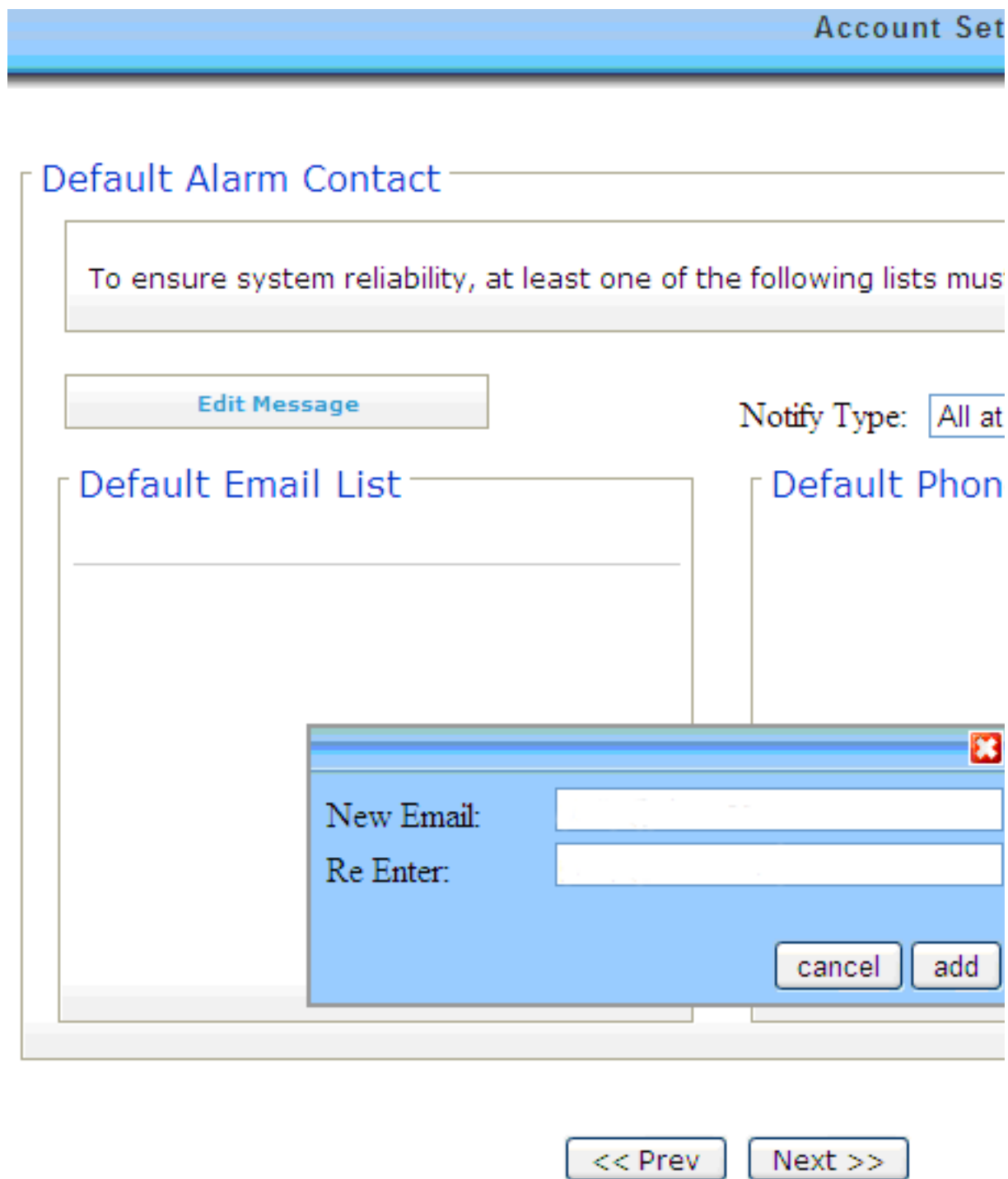
User Accounts

Your Admin and Reader accounts have been setup with default passwords, you must change these passwords, if you haven't already.

<h4 style="margin: 0;">Administrator</h4> <p style="margin: 5px 0;">Login Name*: <input style="width: 80%;" type="text" value="Admin"/></p> <p style="margin: 5px 0;">Password: <input style="width: 80%;" type="text"/></p> <p style="margin: 5px 0;">Repeat Password: <input style="width: 80%;" type="text"/></p> <p style="margin: 5px 0;">First Name: <input style="width: 80%;" type="text"/></p> <p style="margin: 5px 0;">Last Name: <input style="width: 80%;" type="text"/></p>	<h4 style="margin: 0;">Read-Only Account</h4> <p style="margin: 5px 0;">Login Name*: <input style="width: 80%;" type="text"/></p> <p style="margin: 5px 0;">Password: <input style="width: 80%;" type="text"/></p> <p style="margin: 5px 0;">Repeat Password: <input style="width: 80%;" type="text"/></p> <p style="margin: 5px 0;">First Name: <input style="width: 80%;" type="text"/></p> <p style="margin: 5px 0;">Last Name: <input style="width: 80%;" type="text"/></p>
---	---

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Account Set

Default Alarm Contact

To ensure system reliability, at least one of the following lists must

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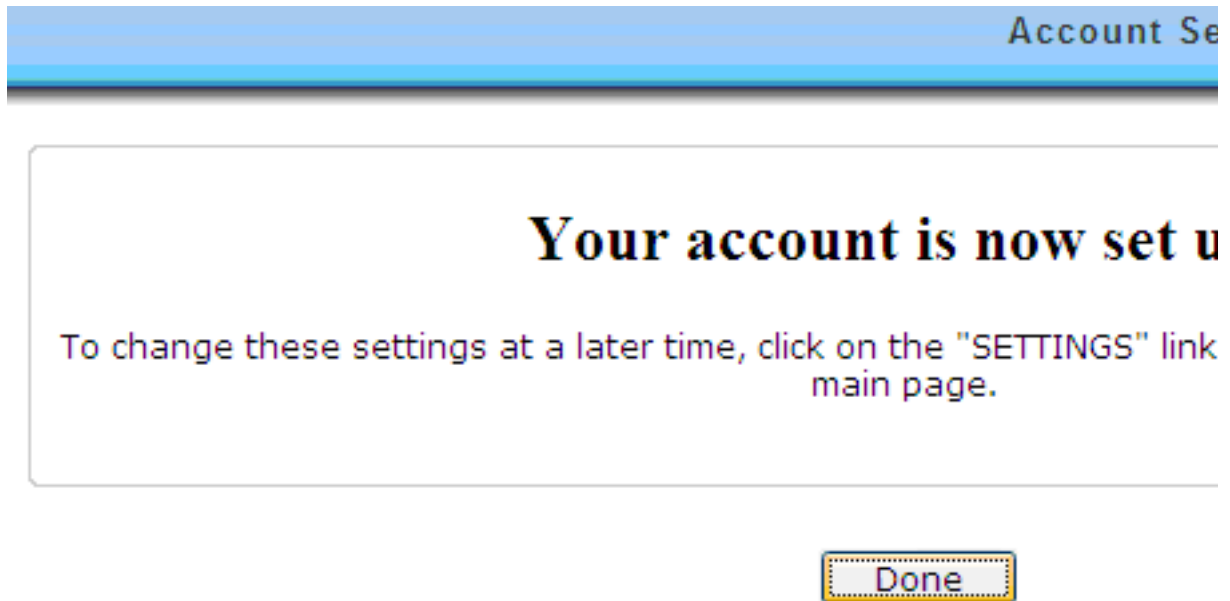
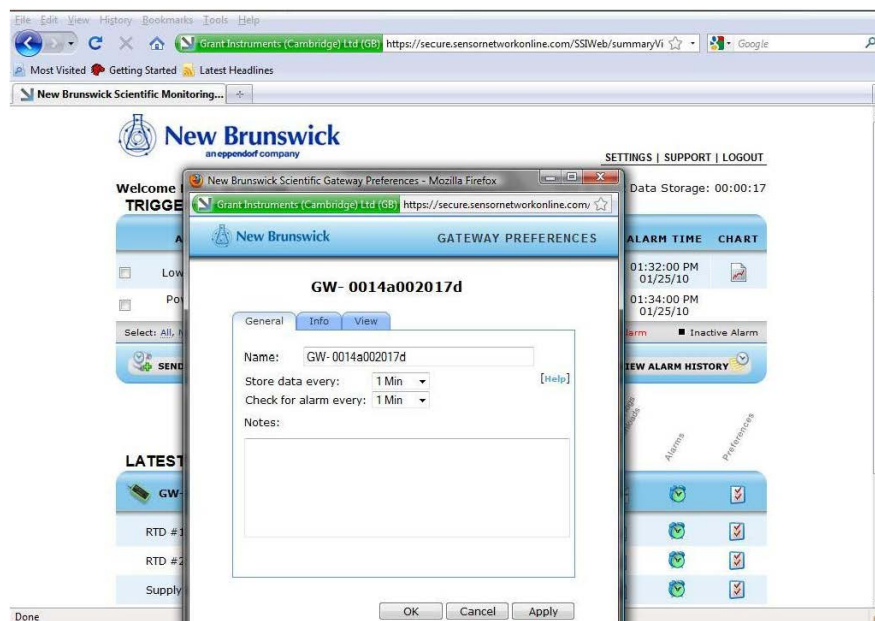
Se clicar nos ícones do visor principal abre as janelas pop up.

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The screenshot shows a software window titled 'New Brunswick SYSTEM SETTINGS'. The 'User Accounts' tab is selected. It contains two main sections: 'Administrator' and 'Read-Only Account'. Each section has a 'Login Name*' field with a value, a 'Password:' field, a 'Repeat Password:' field, a 'First Name:' field with 'Accsense', and a 'Last Name:' field. For the Administrator, the Login Name is 'Admin' and the Last Name is 'Administrator'. For the Read-Only Account, the Login Name is 'Reader' and the Last Name is 'Reader'. At the bottom of the window are 'OK', 'Cancel', and 'Apply' buttons. A 'Done' button is visible at the very bottom left of the window frame.

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The image shows two screenshots from the New Brunswick TCA-3 monitoring system. The top screenshot displays the 'TRIGGERED ALARMS' section with a table of active alarms. The bottom screenshot shows the configuration window for a 'GATEWAY COMMUNICATION ALARM'.

ALARM NAME	POD	MEASUREMENT	VALUE	ALARM TIME	CHART
<input type="checkbox"/> Low supply voltage	GW- 0014a002017d	Supply Voltage (Battery)	2.28	01:32:00 PM 01/25/10	
<input type="checkbox"/> Power or Internet Outage	N/A	N/A		01:34:00 PM 01/25/10	

GATEWAY COMMUNICATION ALARM
GW- 0014a002017d

General

Alarm Enabled

Alarm Description:
Name:
[Add Notes](#)

Notify List:
This will notify everyone in the group.
The list is located in the page.

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The screenshot shows the New Brunswick TCA-3 Temperature Monitoring System interface. At the top, there is a logo for New Brunswick (an eppendorf company) and navigation links for SETTINGS, SUPPORT, and LOGOUT. Below the logo, it says 'Welcome' and 'Time to Next Data Storage: 00:00:25'. The main section is titled 'TRIGGERED ALARMS' and contains a table with columns: ALARM NAME, POD, MEASUREMENT, VALUE, ALARM TIME, and CHART. The table shows one alarm: 'low' for POD 'GW- 0014a002016a', measurement 'RTD #1', value '-245.69', and alarm time '08:58:00 AM 01/05/12'. Below the table are buttons for 'SEND TO HISTORY' and 'VIEW ALARM HISTORY'. Below this is the 'LATEST MEASUREMENTS' section, which shows a table of measurements for device 'GW- 0014a002016a' on '01/05/12 08:58:00 AM'. The measurements are: RTD #1 (-245.69 C), RTD #2 (-237.56 C), Supply Voltage (Battery) (4.05 Volts), Supply Voltage (External) (4.73 Volts), and Thermocouple (1,177.49 C). Each measurement row has three icons: a chart icon, a status icon, and a preference icon. The preference icon for 'RTD #2' is circled in red. Above the preference icons, there are labels: 'Configure Data Log and Downloadable', 'Alarms', and 'Preferences'.

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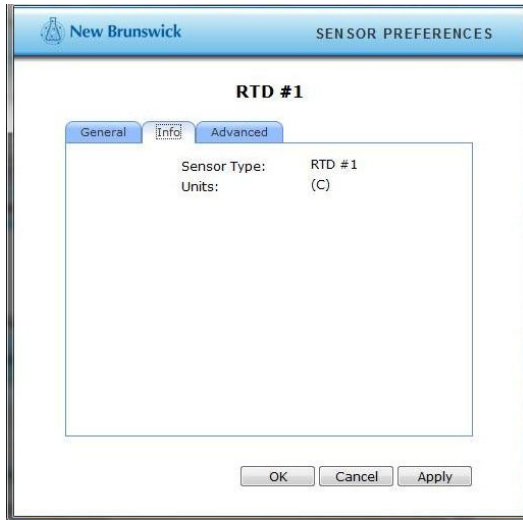
The screenshot shows the 'SENSOR PREFERENCES' dialog box for 'RTD #1'. The dialog has a title bar with the New Brunswick logo and the text 'SENSOR PREFERENCES'. Below the title bar, there are three tabs: 'General', 'Info', and 'Advanced'. The 'General' tab is selected. The 'Name' field contains 'RTD #1'. The 'Units' dropdown menu is set to 'C'. There is a 'Notes' text area below the units. At the bottom of the dialog, there are three buttons: 'OK', 'Cancel', and 'Apply'.

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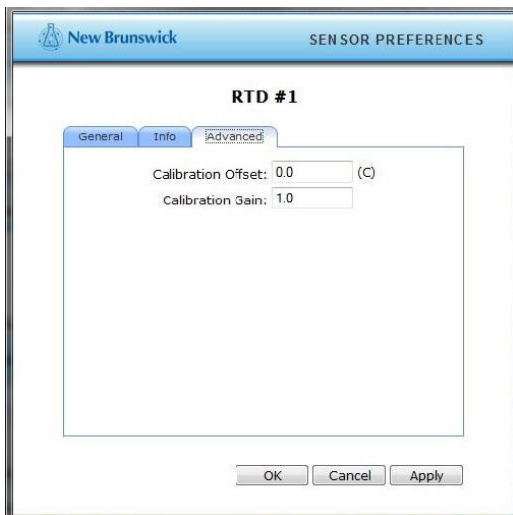
4.11 Alarmes de sensores

The screenshot displays the New Brunswick TCA-3 Temperature Monitoring System interface. At the top, there is a navigation bar with 'SETTINGS | SUPPORT | LOGOUT' and a 'Welcome' message. Below this, the 'TRIGGERED ALARMS' section shows a table with columns: ALARM NAME, POD, MEASUREMENT, VALUE, ALARM TIME, and CHART. A message states 'There are no triggered alarms'. Below this, there are filters for 'Active Alarm' and 'Inactive Alarm', and buttons for 'SEND TO HISTORY' and 'VIEW ALARM HISTORY'.

The 'LATEST MEASUREMENTS' section shows a table with columns: Sensor Name, Measurement Value, and three status icons. The table includes the following data:

Sensor Name	Measurement Value	Status 1	Status 2	Status 3
GW- 0014a002016a	01/05/12 09:19:00 AM	[Green Check]	[Green Check]	[Green Check]
RTD #1	-245.76 C	[Green Check]	[Green Check]	[Green Check]
RTD #2	-237.61 C	[Green Check]	[Alarm Icon]	[Green Check]
Supply Voltage (Battery)	4.05 Volts	[Green Check]	[Green Check]	[Green Check]
Supply Voltage (External)	4.73 Volts	[Green Check]	[Green Check]	[Green Check]
Thermocouple	1,177.51 C	[Green Check]	[Green Check]	[Green Check]

The 'Alarm Icon' in the 'RTD #2' row is circled in red. The interface also includes a 'Generate Data Logs and Downloads' button and a 'Preferences' link.

Fig. 4-4: Ícone sensor de alarme

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Alarm Description

Name: Fridge [Add Notes](#)

Notify List: Default List
This will notify everyone on the "Default Notify List".
The list is located in the "Settings" menu on the main page.

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New Brunswick ALARM DEFINITION

GW- 0014a002016a - RTD #1

General Trigger Email List Phone List

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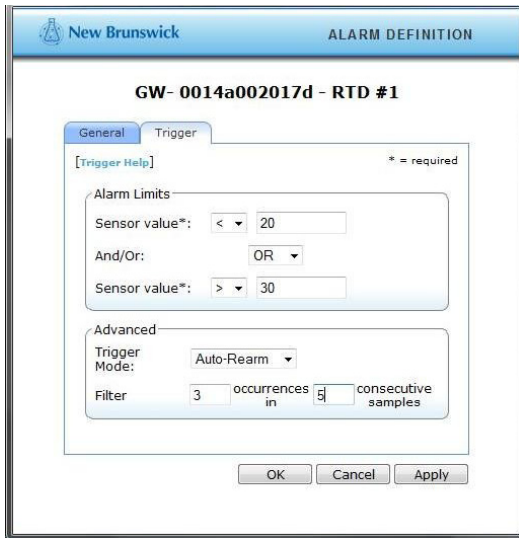
4.12 Gráfico sensorial

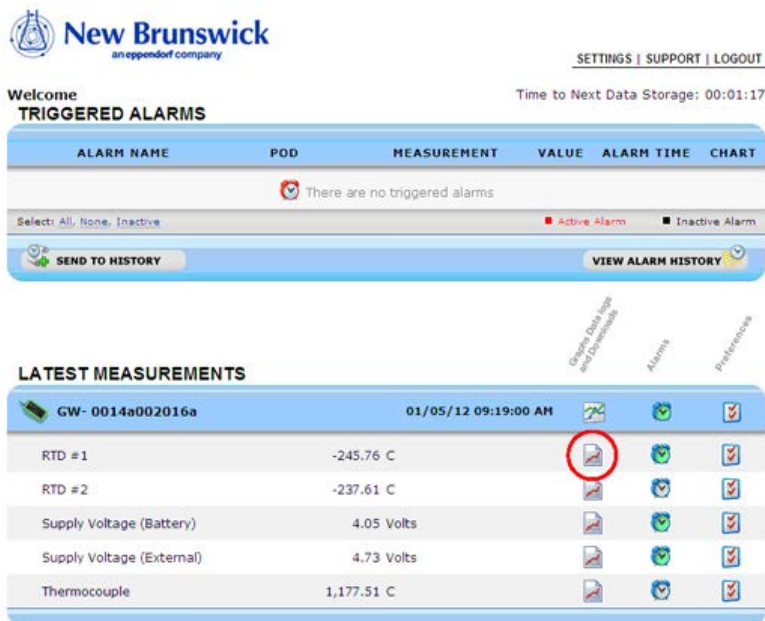
Fig. 4-5: Ícone de gráfico sensorial

OperaçãoTCA-3 Temperature Monitoring System
Brazilian Portuguese (BR)

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7 Certificados

Declaration of Conformity

The product named below fulfills the requirements of directives and standards listed. In the case of unauthorized modifications to the product or an unintended use this declaration becomes invalid.

Product name:

TCA-3
including accessories

Product type:

Remote logging temperature monitor

Relevant directives / standards:

2004/108/EC EN 61326-1
2006/95/EC EN 61010-1
2011/65/EU

Management Board

Portfolio Management

Date: January 06, 2014

Your local distributor: www.eppendorf.com/contact
Eppendorf AG · 22331 Hamburg · Germany
eppendorf@eppendorf.com

ISO 9001
Certified

ISO 13485
Certified

ISO 14001
Certified

MI288-9999-00

Eppendorf® and the Eppendorf Logo are registered trademarks of Eppendorf AG, Hamburg/Germany. All rights reserved incl. graphics and pictures. Copyright 2013 © by Eppendorf AG.

Certificados

TCA-3 Temperature Monitoring System
Brazilian Português (BR)

Evaluate your manual

Give us your feedback.

www.eppendorf.com/manualfeedback